LABCONCO CORPORATION

8811 Prospect Avenue
Kansas City, MO 64132
(800) 821-5525, (816) 333-8811
(816) 363-0130 fax
labconco@labconco.com

User's Manual

Protector[®] Pass-Through Laboratory Fume Hoods

Models

1134 Series
1135 Series
1136 Series

To receive important product updates,
complete your product registration card
online at register.labconco.com

Please read the User's Manual before operating the equipment.

Copyright © 2021 Labconco Corporation. The information contained in this manual and the accompanying products are copyrighted and all rights reserved by Labconco Corporation. Labconco Corporation reserves the right to make periodic design changes without obligation to notify any person or entity of such change.

Warranty

Labconco Corporation provides a warranty to the original buyer for the repair or replacement of parts and reasonable labor as a result of normal and proper use of the equipment with compatible chemicals. Broken glassware and maintenance items, such as filters, gaskets, light bulbs, finishes and lubrication are not warranted. Excluded from warranty are products with improper installation, erratic electrical or utility supply, unauthorized repair and products used with incompatible chemicals.

The warranty for Protector® Pass-Through Laboratory Fume Hoods will expire one year from date of installation or two years from date of shipment from Labconco, whichever is sooner. Warranty is non-transferable and only applies to the owner (organization) of record.

Buyer is exclusively responsible for the set-up, installation, verification, decontamination or calibration of equipment. This limited warranty covers parts and labor, but not transportation and insurance charges. If the failure is determined to be covered under this warranty, the dealer or Labconco Corporation will authorize repair or replacement of all defective parts to restore the unit to operation. Repairs may be completed by 3rd party service agents approved by Labconco Corporation. Labconco Corporation reserves the rights to limit this warranty based on a service agent's travel, working hours, the site's entry restrictions and unobstructed access to serviceable components of the product.

Under no circumstances shall Labconco Corporation be liable for indirect, consequential, or special damages of any kind. This warranty is exclusive and in lieu of all other warranties whether oral, or implied.

Returned or Damaged Goods

Do not return goods without the prior authorization from Labconco. Unauthorized returns will not be accepted. If your shipment was damaged in transit, you must file a claim directly with the freight carrier. Labconco Corporation and its dealers are not responsible for shipping damages.

The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

Limitation of Liability

The disposal and/or emission of substances used in connection with this equipment may be governed by various federal, state, or local regulations. All users of this equipment are required to become familiar with any regulations that apply in the user's area concerning the dumping of waste materials in or upon water, land, or air and to comply with such regulations. Labconco Corporation is held harmless with respect to user's compliance with such regulations.

Contacting Labconco Corporation

If you have questions that are not addressed in this manual, or if you need technical assistance, contact Labconco's Customer Service Department or Labconco's Product Service Department at 1-800-821-5525 or 1-816-333-8811, between the hours of 7:30 a.m. and 5:30 p.m., Central Standard Time.

TABLE OF CONTENTS

CHAPTER 1: INTRODUCTION	1
About This Manual	2
Typographical Conventions (Conventions Typographiques)	3
Your Next Step	4
CHAPTER 2: PREREQUISITES	5
Location Requirements	6
Support Requirements	6
Exhaust Requirements (Exigences électriques)	6
Airflows & Static Pressure	7
Electrical Requirements	7
Service Line Requirements	8
Space Requirements	8
CHAPTER 3: GETTING STARTED	9
Unpacking Your Fume Hood	10
Removing the Shipping Skid	10
Sash Weight Release	11
Installing the Hood on a Supporting Structure & Work Surface	12
Connecting to the Hood Exhaust System	13
Connecting the Electrical Supply Source	13
Connecting the Service Lines to the Protector Hood	16
Sealing the Protector Hood to the Work Surface	17
Certifying the Protector Fume Hood	17
CHAPTER 4: PERFORMANCE FEATURES AND SAFETY PRECAUTIONS	18
Performance Features	18
Safety Precautions (Consignes de sécurité)	21

CHAPTER 5: USING YOUR PROTECTOR PASS-THROUGH	25
Operating the Vertical Rising Sash	25
Operating the Blower	25
Operating the Lights	25
Working in your Protector Fume Hood	26
CHAPTER 6: MAINTAINING YOUR PROTECTOR PASS-THROUGH	28
Service Safety Precautions (Mesures de sécurité des services)	28
Routine Maintenance Schedule	29
Routine Service Operations	30
CHAPTER 7: MODIFYING YOUR PROTECTOR PASS-THROUGH	31
Installing Work Surfaces	31
Installing Ceiling Enclosures Above the Fume Hood	32
Installing Additional Service Fixtures	32
Installing Guardian Airflow Monitors	33
Sash Stop Kit – Field Installation	33
Installing an Electrical Duplex Outlet (Installation d'une sortie Duplex électrique)	34
CHAPTER 8: TROUBLESHOOTING	35
APPENDIX A: PROTECTOR PASS-THROUGH COMPONENTS	38
APPENDIX B: PROTECTOR PASS-THROUGH DIMENSIONS	41
APPENDIX C: PROTECTOR PASS-THROUGH SPECIFICATIONS	42
(ANNEXE C: SPÉCIFICATIONS PROTECTEUR DU PASS-THROUGH MINISTRE)	
APPENDIX D: SERIAL NUMBER TAG DESCRIPTION	44
(ANNEXE D: DESCRIPTION DE SÉRIE TAG NUMBER)	
APPENDIX E: REFERENCES	50
APPENDIX F: DISASSEMBLY AND REASSEMBLY INSTRUCTIONS	53

CHAPTER 1

INTRODUCTION

Congratulations on your purchase of a Labconco Protector® Pass-Through Laboratory Fume Hood. Your Protector Pass-Through Laboratory Fume Hood is designed to protect you. It is the result of Labconco's more than 60 years experience in manufacturing fume hoods, and users like you suggested many of its features to us.

The Labconco Protector Pass-Through Fume Hood has been engineered to provide maximum visibility in a laboratory, and effectively contain toxic, noxious, or other harmful materials when properly installed. The Protector offers many unique features to enhance safety, performance, and visibility. To take full advantage of them, please acquaint yourself with this manual and keep it handy for future reference. If you are unfamiliar with how fume hoods operate, please review *Chapter 4: Performance Features and Safety Precautions* before you begin working in the fume hood. Even if you are an experienced fume hood user, please review *Chapter 5: Using Your Protector Pass-Through*, which describes your Protector Hood's features so that you can use the hood efficiently.

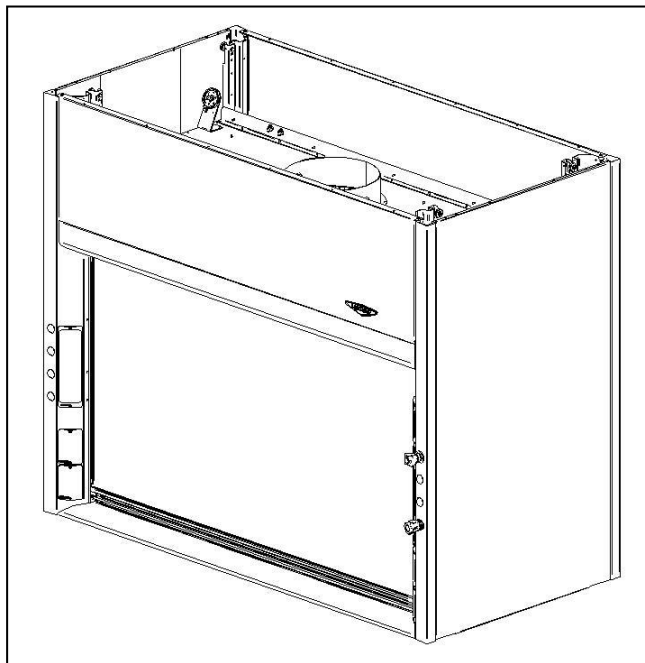
Figure 1-1

About This Manual

This manual is designed to help you learn how to install, use, and maintain your laboratory fume hood. Instructions for installing optional equipment on your hood are also included.

Chapter 1: Introduction provides a brief overview of the laboratory fume hood, explains the organization of the manual, and defines the typographical conventions used in the manual.

Chapter 2: Prerequisites explains what you need to do to prepare your site before you install your laboratory fume hood. Electrical and service requirements are discussed.

Chapter 3: Getting Started contains the information you need to properly unpack, inspect, install, and certify your laboratory fume hood.

Chapter 4: Performance Features and Safety Precautions explains how the Protector operates and the appropriate precautions you should take when using the fume hood.

Chapter 5: Using Your Protector Pass-Through discusses the basic operation of your fume hood. Information on how to prepare, use and shut down your Protector Hood are included.

Chapter 6: Maintaining Your Protector Pass-Through explains how to perform routine maintenance on your fume hood.

Chapter 7: Modifying Your Protector Pass-Through explains how to modify the fume hood or add accessories.

Chapter 8: Troubleshooting contains a table of problems you may encounter while using your laboratory fume hood including the probable causes of the problems and suggested corrective actions.

Appendix A: Protector Pass-Through Components contains labeled diagrams of all of the components of the fume hoods.

Appendix B: Protector Pass-Through Dimensions contains comprehensive diagrams showing all of the dimensions for the laboratory fume hoods.

Appendix C: Protector Pass-Through Specifications contains the electrical requirements for laboratory fume hood. Wiring diagrams are also included.

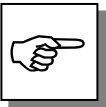
Appendix D: Serial Number Tag Description provides current rating code used on serial number tag.

Appendix E: References lists the various resources available that deal with laboratory fume hoods.

Appendix F: Disassembly and Reassembly Instructions lists the required tools as well as step by step instructions for disassembling and reassembling the Protector XL Hood.

Typographical Conventions


Recognizing the following typographical conventions will help you understand and use this manual:

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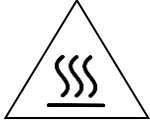- CAUTION – Hot Surface.
- AVERTIR – Surface Chaude

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Your Next Step

If your Fume Hood needs to be installed, proceed to *Chapter 2: Prerequisites* to ensure your installation site meets all of the requirements. Then, go to *Chapter 3: Getting Started* for instructions on how to install your laboratory fume hood and make all of the necessary connections.

If you would like to review how laboratory fume hoods operate, go to *Chapter 4: Performance Features and Safety Precautions*.

For information on the operational characteristics of your laboratory fume hood, go to *Chapter 5: Using Your Protector Pass-Through*.

If your laboratory fume hood is installed and you need to perform routine maintenance on the cabinet, proceed to *Chapter 6: Maintaining Your Protector Pass-Through*.

For information on making modifications to the configuration of your fume hood, go to *Chapter 7: Modifying Your Protector Pass-Through*.

Refer to *Chapter 8: Troubleshooting* if you are experiencing problems with your fume hood.

CHAPTER 2

PREREQUISITES

Before you install your laboratory fume hood, you need to prepare your site for installation. Carefully examine the location where you intend to install your hood. You must be certain that the area is level and of solid construction. In addition, a dedicated source of electrical power must be located near the installation site.

Carefully read this chapter to learn the requirements for your installation site:

- The location requirements.
- The support requirements.
- The exhaust requirements.
- The electrical power requirements.
- The service line requirements.
- The space requirements.

Refer to *Appendix B: Protector Pass-Through Dimensions* for complete fume hood dimensions.

Refer to *Appendix C: Protector Pass-Through Specifications* for complete laboratory fume hood electrical and environmental conditions, specifications and requirements.

Location Requirements

The fume hood should be located away from traffic patterns, doors, windows, fans, ventilation registers, and any other air-handling device that could disrupt its airflow patterns. All windows in the room should be closed.

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Support Requirements

DO NOT install the fume hood on a cart, dolly, or mobile bench. **ALL** Protector Hood installations must be permanent and stationary. The supporting structure usually consists of a base cabinet and chemically resistant work surface.

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Exhaust Requirements

The exhaust duct connection has been designed for 12" nominal duct (12.75" OD) to allow for minimum static pressure loss while operating at 60 to 100 fpm face velocities. The 12" diameter exhaust duct also allows for proper transport velocities away from the hood in the 1000 fpm to 2500 fpm range. The proper exhaust volume and static pressure loss are listed next for each hood model:

Airflows and Static Pressure

Table 2-1 lists the Airflow and Static Pressure for Protector Pass-Through fume hoods based upon the desired sash opening height.

Table 2-1

Face Velocity (fpm)		Airflow Volumetric Rate (CFM) @ Static Pressure (inches H ₂ O)		
One Sash Full Open at 25"	One Sash 62.5% Open at 18"	4' Hood	5' Hood	6' Hood
100	130	775 @ 0.33	1020 @ 0.42	1250 @ 0.59
NA	100	590 @ 0.19	775 @ 0.24	950 @ 0.34

Proper blower selection can be determined from these exhaust requirements and the total system static pressure loss. Contact Labconco Customer Service for assistance in sizing a blower system.

Electrical Requirements

The Protector Hood models feature internal wiring for the fluorescent light assembly and light switch. All internal wiring is terminated at the single point wiring junction box for hook-up by a qualified electrician. The blower switch, and light switch wires are also terminated at the single point wiring junction box for hook-up by a qualified electrician. Refer to *Chapter 3: Getting Started* and *Appendix C: Protector Pass-Through Specifications* for the wiring diagram for proper electrical installation.

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Service Line Requirements

All service lines to the laboratory fume hood should be ¼ inch outside diameter, copper (brass for natural gas), and equipped with an easily accessible shut-off valve, should disconnection be required. Recommended operating pressure is 40 PSI, with a maximum allowable pressure of 200 PSI. Consider a pressure regulator to reduce line pressure to 40 PSI. Please check with local codes for other requirements.

Space Requirements

The dimensions for the different models are shown in *Appendix B: Protector Pass-Through Dimensions*.

CHAPTER 3

GETTING STARTED

Now that the site for your laboratory fume hood is properly prepared, you are ready to unpack, inspect, install, and certify your unit. Read this chapter to learn how to:

- Unpack and move your Protector Hood.
- Set up the fume hood with the supporting structure and work surface.
- Connect to an exhaust system.
- Connect the electrical supply source.
- Connect the service lines.
- Sealing the Protector Hood to the work surface.
- Arrange certification of your Protector Hood.

Depending upon which model you are installing, you may need common plumbing and electrical installation tools in addition to 5/16", 3/8", 7/16", and 1/2" wrenches, ratchets, sockets, a nut driver set, a flat-blade screwdriver, a Phillips screwdriver, and a carpenter level to complete the instructions in the chapter.

The Protector Pass-Through Hood models weigh between 400 to 800 lbs. (182-363 kg). The shipping skid allows for lifting with a mechanical lift truck or floor jack. If you must lift the fume hood manually, follow safe-lifting guidelines. Normally, the fume hood can be slid off a hydraulic lift table and be placed into position on top of the work surface. Do not lift by the front air foil.

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
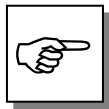
Unpacking Your Fume Hood

Carefully remove the shrink-wrap or carton on your fume hood and inspect it for damage that may have occurred in transit. If your unit is damaged, notify the delivery carrier immediately and retain the entire shipment intact for inspection by the carrier.

The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

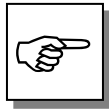DO NOT RETURN GOODS WITHOUT THE PRIOR AUTHORIZATION OF LABCONCO. UNAUTHORIZED RETURNS WILL NOT BE ACCEPTED.

IF YOUR HOOD WAS DAMAGED IN TRANSIT, YOU MUST FILE A CLAIM DIRECTLY WITH THE FREIGHT CARRIER. LABCONCO CORPORATION AND ITS DEALERS ARE NOT RESPONSIBLE FOR SHIPPING DAMAGES.

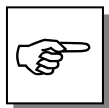
Do not discard the shipping skid or packing material for your fume hood until you have checked all of the components and installed and tested the unit. Do not remove the fume hood from its shipping skid until it is ready to be placed into its final location. Move the unit by placing a flat, low dolly under the shipping skid, or by using a floor jack.

Do not move the hood by tilting it onto a hand truck.

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Removing the Shipping Skid

LEAVE THE FUME HOOD ATTACHED TO ITS SHIPPING SKID UNTIL IT IS AS CLOSE TO ITS FINAL LOCATION AS POSSIBLE. MOVE THE HOOD BY USING A SUITABLE FLOOR JACK, OR BY PLACING A FURNITURE DOLLY UNDERNEATH THE SKID. DO NOT MOVE THE HOOD BY TILTING IT ONTO A HAND TRUCK.

After you verify the fume hood components, move your hood to the location where you want to install it. Should you require disassembly to move the hood, then follow the instructions in *Appendix F*. Then, follow the steps listed next to remove the shipping skid from your unit.

1. Remove side panels by removing the two (2) Phillips screws atop each side panel.

2. Find the hardware (bolts, washers, nuts) that attach the fume hood to the skid and remove the hardware. The hardware is located on the sides of the fume hood now exposed with the side panels removed.

Sash Weight Release

To protect the fume hood from damage in shipment, the sash weights have been disconnected from the sash cables and secured to the side frame with a bolt. As shown in Fig. 3-1 remove the two bolts and ensure sash cables are on pulleys before attaching the weights. To access sash weights, remove Weight Cover.


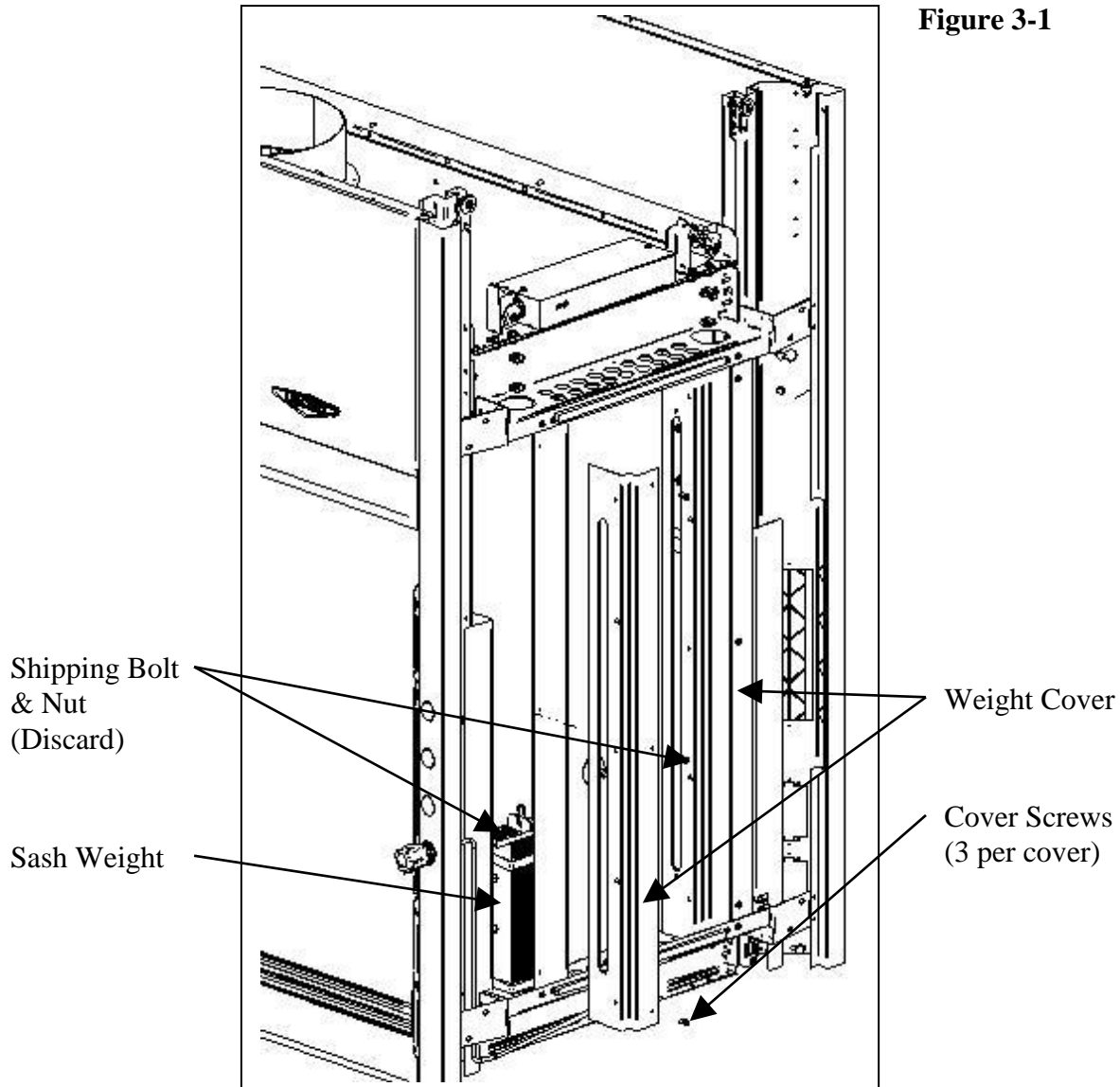
Figure 3-1

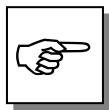NOTE: THE SASH WEIGHTS ARE INDIVIDUALLY MATCHED FOR THIS SPECIFIC HOOD AND SHOULD NOT BE EXCHANGED ON ANY OTHER UNIT.

Installing the Hood on a Supporting Structure and Work Surface

The Protector Hood is heavy! Use caution when lifting or moving the unit.

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When installing the Protector Fume Hood onto a chemically-resistant work surface or benchtop, ensure that the structure can safely support the combined weight of the fume hood and any related equipment. The work surface should be at least as wide as the hood to properly support it. **Position front edge of fume hood 1.7" past front edge of worksurface. Lower base cabinets or casework is placed flush with front edge of worksurface as shown in Figure 3-2.**

The work surface should be smooth and durable, such as a chemically-resistant epoxy resin. The surface should be nonporous and resistant to the acids, solvents, and chemicals used in conjunction with the Protector Fume Hood. The work surface should also contain a dished recessed area for containing primary spills.

NOTE: Labconco recommends supports by others for underneath the work surface at time of installation.

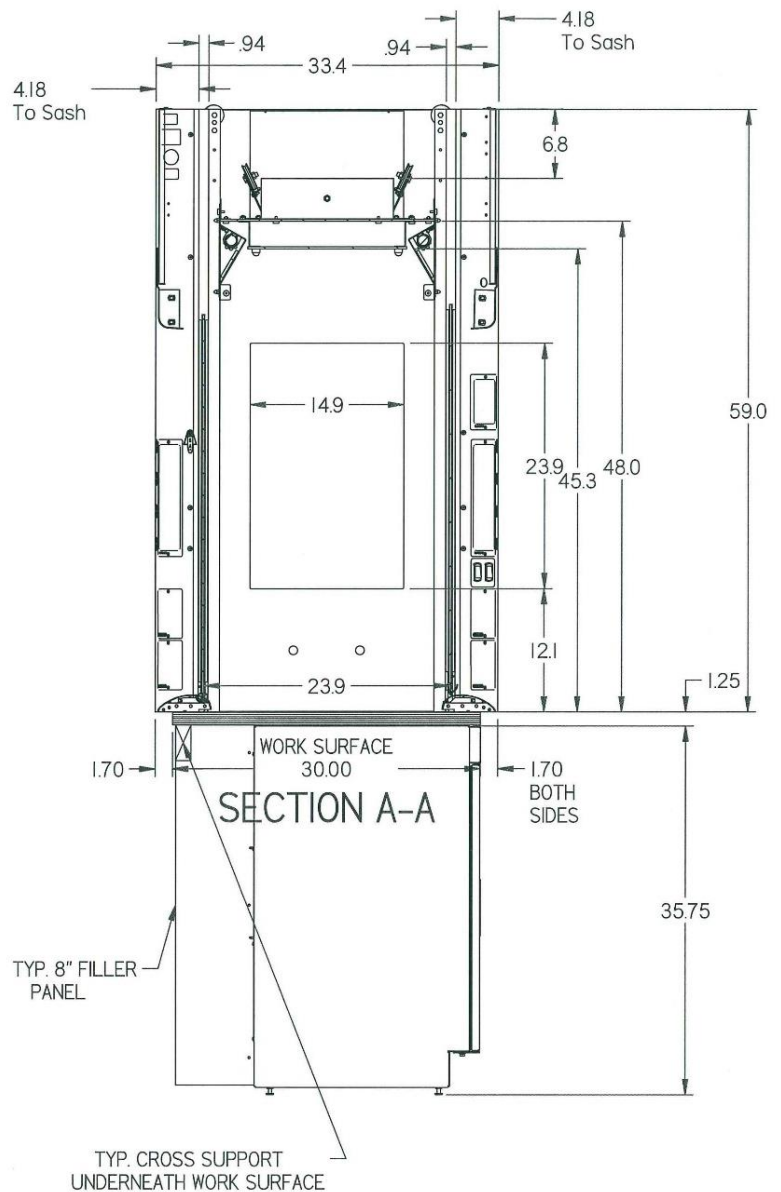
Figure 3-2

Connecting to the Hood Exhaust System

WARNING: The weight of the exhaust ductwork system must be supported independently of the hood superstructure. Do not allow this weight to be supported by the hood structure as damage to the hood may occur.

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The exhaust connection should be installed by a qualified HVAC contractor. The exhaust connection on your hood has been designed for 12" nominal pipe (12.75" OD) to allow for minimum static pressure loss with proper transport velocities away from the hood. Consult Labconco Customer Service should you require help sizing your blower for the exhaust volume and total system static pressure loss.

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The selected exhaust duct material should match the hood procedures and chemicals used to ensure compatibility.

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Connecting the Electrical Supply Source to the Protector Fume Hood

Prior to connecting any electrical wiring to the fume hood structure, refer to the hood identification plate for the proper electrical requirements of your specific model.

Chapter 3: Getting Started

WARNING: The building electrical supply system for Protector Hoods should include overload protection. A switch or circuit breaker should be in close proximity to the equipment and within easy reach of the operator. The switch or circuit breaker is to be marked as the disconnecting device for the equipment.

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The identification plate, model number, serial number, and electrical connection boxes are accessible from the front of the fume hood by removing the front panel.

The Protector Pass-Through Hood is normally wired for 115 Volt, 50/60 Hz, 20 Amp or 230Volt, 50/60 Hz, 10 Amp electrical service. Check the I.D. plate behind the front panel for voltage verification. The number of circuits varies depending on the model. All of the electrical connections are terminated at the field wiring terminal box for hook-up by a qualified electrician. We recommend each circuit be a dedicated branch circuit. However, if wired together the maximum load allowed is the sum of individual outlets plus the rating of the unit (i.e. 2 Amps). The single point internal junction box is used for the connection of the lights, blower, and duplex outlets. Refer to the wiring diagram for your Protector XL in *Appendix C: Protector Pass-Through Specifications*.

The fume hood is required to be grounded to the MAINS protective earthing ground for safe operation. Using a ring terminal sized for a 10-24 machine screw, connect the MAINS ground conductor to the grounding lug marked with the protective earthing symbol, . Only MAINS ground conductors should be connected to the protective earthing ground lug, no other conductors should be connected to this grounding lug. Using wire nuts connect the MAINS supply conductors to the fume hood supply wires. Insure that the wires are connected as per the appropriate wire color codes for the input voltage. For 115V Phase (Hot) is black and Neutral is white, for 230V Phase1 is brown and Phase2 is blue. Refer to the wiring diagram for your Protector XL in *Appendix C: Protector Pass-Through Specifications*.

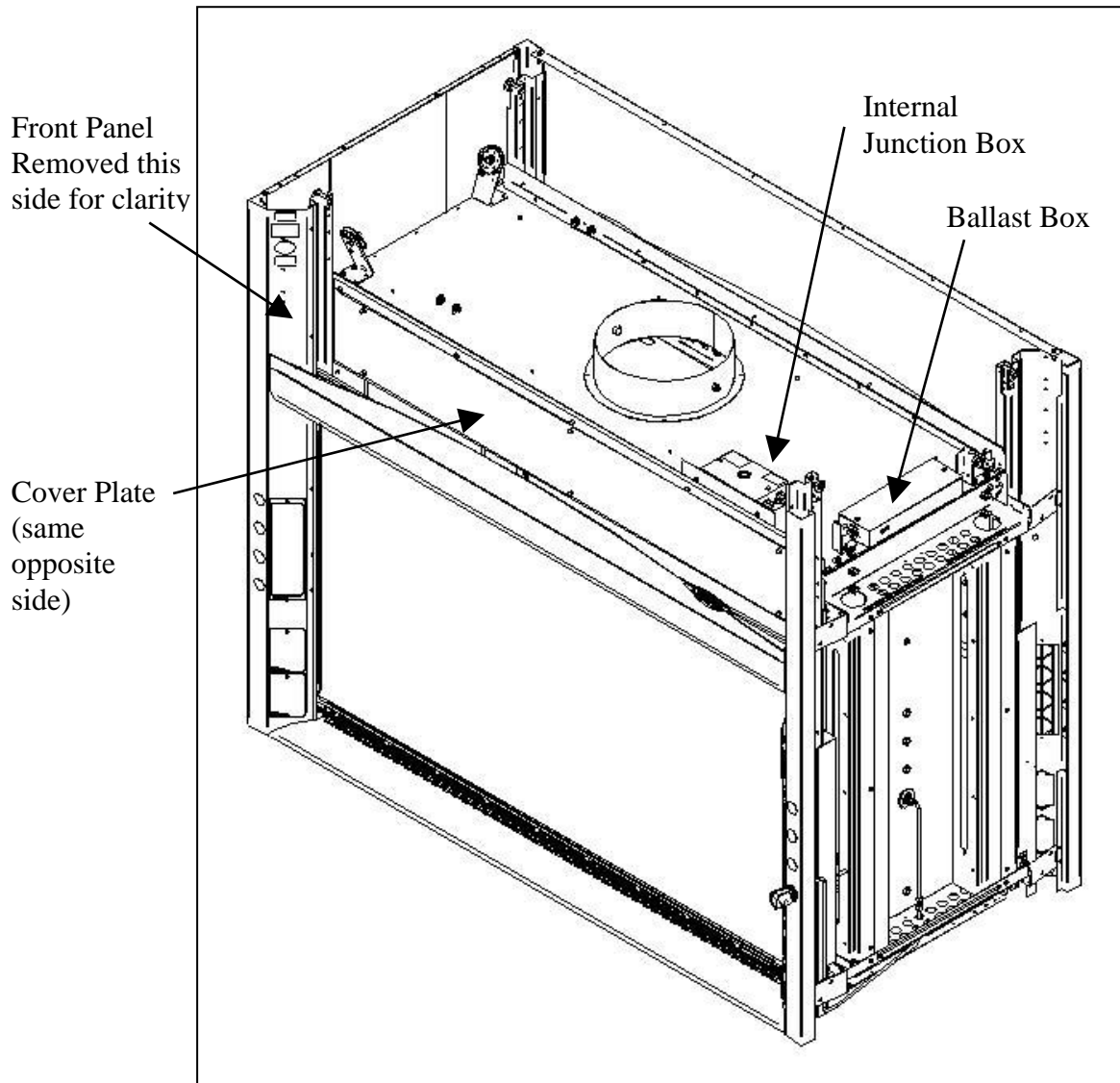
Figure 3-3

All wiring for the fume hood SHOULD be performed by a licensed electrician and conform to all local codes. In most cases, the hood will require the use of shielded conduit to protect the wiring into the hood. The grounding connection shall not be made to the terminal box cover.

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Chapter 3: Getting Started

The fluorescent light has been mounted behind the front panel and is sealed from vapors inside the hood structure. The fluorescent ballast is located in a separate box on top of the fume hood adjacent to the main junction box. To change the fluorescent light bulbs in your hood, you must first remove the front panel(s) and then the cover plate(s), which is secured with six (6) acorn nuts (See Fig. 3-3). Twist and remove any defective bulbs. Reverse order to reassemble.

Connecting the Service Lines to the Protector Fume Hood

The hoods with service fixtures have been plumbed from the valve to the hose connector or gooseneck for your installation convenience. Supply tubing shall be provided by the qualified installer. Tubing can enter the hood from above, through the back, or through the work surface to make these connections to the service fixtures.

NOTE: Inspect all fittings for leakage. Tighten the fittings slightly if needed.

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CAUTION: Do not use oxygen with any standard service fixture. Contact Labconco Customer Service for oxygen fixture information.

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Should access to the hood plumbing fixture bodies be required, remove the service access plate on the hood front corner posts by loosening their individual screws (see item 11, Figure A-1 in *Appendix A*). The valve body will now be fully exposed for any service work that may be necessary. The service fixtures supplied on your laboratory hood are designed for use with the following services:

- Air
- Cold Water
- Hot Water
- Natural Gas – See caution below
- Vacuum

WARNING: Contact Labconco Customer Service directly before using any service other than those listed above in these valves to assure full compatibility.

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CAUTION: Natural gas should be used only in the service fixture that has been pre-plumbed with brass tubing. Sulfur content of the gas could cause deterioration of standard copper supply lines.

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Sealing the Protector Hood to the Work Surface

When the hood has been set in place, ducted, wired, and plumbed, it should be sealed at the work surface to prevent spilled materials from collecting under the walls of the hood. Materials such as silicone sealants are recommended to seal the hood structure.

Certifying the Protector Fume Hood

The combination of your laboratory hood, exhaust ductwork, and exhaust blower gives you the flexibility to change the airflow at the sash opening of your hood. To determine the actual face velocity at the sash opening, airflow velocity readings will need to be taken. This should be done across the sash opening of the hood in accordance with the *Industrial Ventilation Manual* section on laboratory hoods (see *Appendix E: References*). For best containment on Pass-Through Hoods, Labconco recommends an average face velocity at the sash opening of 100 feet per minute. Consult Chapter 2 for proper airflow volumes for your particular model.

Your Protector Fume Hood has been tested at the factory per ASHRAE 110-1995. All hoods achieve an “as manufactured rating” of less than 0.05 part per million (ppm) at 4 liters per minute (lpm); AM<0.05 (consult Labconco for individual fume hood ratings). For “field use” ASHRAE testing contact Labconco Sales Engineering Team or Customer Service for a certified on-site contractor.

NOTE: Face velocity profiles and smoke testing should be done periodically to ensure safe performance.

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CHAPTER 4

PERFORMANCE FEATURES AND SAFETY PRECAUTIONS

Performance Features

The Protector Pass-Through Laboratory Hood is designed to meet the needs of the laboratory scientist at OSHA approved velocities at 100 feet per minute. The laboratory fume hood has been designed to effectively contain toxic, noxious, or other harmful materials when properly installed. A fully featured by-pass hood with baffle and air foil, this hood maintains safe airflow. Optional energy saving A-Style Combination Sash models are also available. The hood features by-pass airflow design that promotes full containment as the sash is moved. Airflow is diverted behind the front panel and under the air foil to help control fluctuations in face velocity, which occur as the sash is closed.

1. **Unique sash provides maximum visibility of 37.5" high while conserving energy by limiting sash travel to 25"**. Vertical-rising sash may be raised from a closed to 28" operating height. The fume hood features a containment-enhancing sash handle. Optional sash stops are available to limit sash height and reduce energy usage.
2. **By-pass airflow design** ensures containment as the sash moves.
3. **Large usable interior work depth and interior height of 48"** provides ample working space.
4. **Exterior access cover plates** are removable for easy access to plumbing valves and sash adjustment hardware when access through the sides is not available.
5. **Lift-Away™ front panel** provides easy access to electrical wiring, and lighting fixtures.

6. **Energy efficient fluorescent lighting** is located behind a laminated safety glass shield mounted to the top of the hood. The factory-wired instant start T8 lighting is serviceable from outside the hood cavity. If needed, more energy efficient, long lasting 50,000 hour LED direct replacement T8 bulbs can be ordered. See Appendix A.
7. **Low mounted, factory-wired light and blower switches** are ADA compliant.
8. **All Air Foils** allow air to sweep the work surface for maximum containment. The Clean-Sweep™ openings create a constant barrier from contaminants. In addition, should the operator inadvertently block the airflow entering, the air continues to enter from under the air foil and through the Clean-Sweep openings. See Appendix A for alternate PVC Eco-Air Foil as PVC material resists corrosion from mineral acids.
9. **Streamlined corner posts** provide maximum visibility and the flexibility to add services after installation.
10. **All hoods are factory prepared for up to 8 service fixtures** (4 service fixtures if Side View Window option is selected).
11. **Duplex electrical receptacles** are mounted on the right and left corner posts as requested. Receptacles are factory-wired to hood single point junction box.
12. **Shipped fully assembled** and eliminates the need for costly onsite assembly.
13. **Accessory Guardian™ Digital Airflow Monitor** or **Guardian Airflow Monitor** continuously monitors face velocity. An audio/visual alarm alerts the user to low airflow conditions. The right corner post is factory prepared to accommodate the Guardian Monitor (sold separately).
14. **Sash Alarm.** Audible alarm sounds if both sashes are raised to alert user.
15. **Frame of epoxy-coated steel and aluminum** is durable and corrosion resistant.
16. **Exhaust Connection.** The hood features 12" (12.75" OD pipe) exhaust connections sized to allow for a minimum static pressure loss through the hood structure while providing a good transport velocity through the exhaust system.
17. **Spillstopper™ Solid Epoxy Work Surface** is dished to contain spills. (Work surface is sold separately).
18. **Optional Ceiling Enclosure Kits** are available for a decorative facade between the hood and the ceiling.
19. **Optional Sash Stops** provide a means of controlling the operating height of the sash.

Chapter 4: Performance Features & Safety Precautions

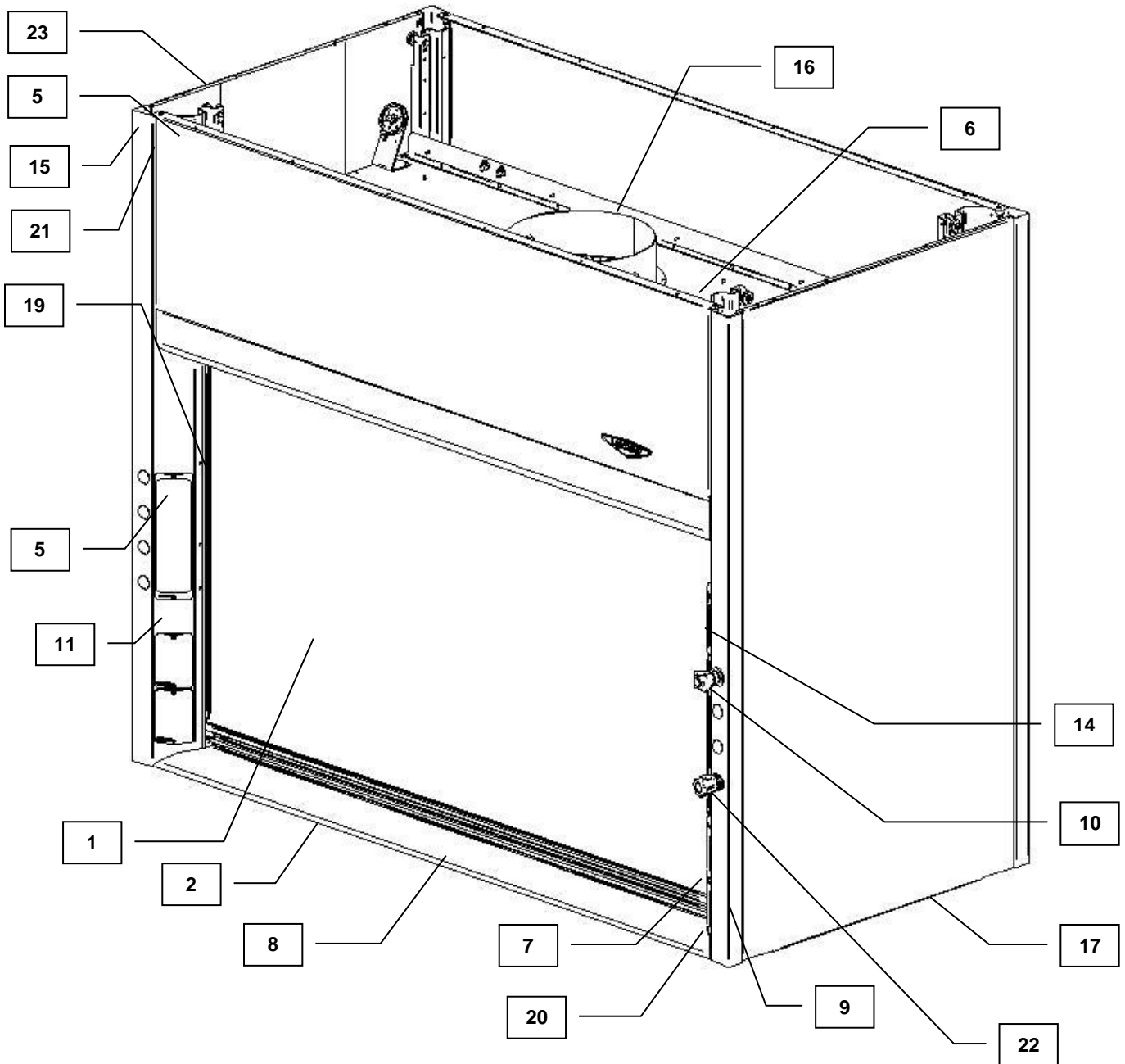
Figure 4-1

See Diagram on Page 20.

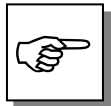
20.  **CAUTION** – Current rating of receptacle is specified in *Appendix D*.
AVERTIR – Classification des prises de courant est spécifié dans l'Annexe D.
21.  **CAUTION** – See Appendix C and D for complete current rating.
AVERTIR – Voir Annexe C et D pour la classification complète du courant.
22.  **CAUTION** – Flammable Gas.
AVERTIR – Gaz Inflammable.
23. **Optional Side Viewing Window** is available on some models and provides an additional observation vantage point.

Safety Precautions

Although the laboratory hood has been engineered to maintain optimum operator safety, caution should always be used while working in the hood. Prior to using the hood, check to make sure that the exhaust blower is operating and that air is entering the hood at its specified face velocity.

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USE GOOD HOUSEKEEPING IN THE HOOD AT ALL TIMES. CLEAN UP SPILLS IMMEDIATELY WITH A MILD DETERGENT. PERIODICALLY CLEAN HOOD INTERIOR, INCLUDING FLUORESCENT LIGHT GLASS PANEL. REPLACE BURNED OUT LIGHT BULBS TO MAINTAIN MAXIMUM ILLUMINATION.

THE LAMP(S) IN THIS PRODUCT CONTAIN MERCURY

Manage in accordance with local disposal laws. DO NOT place lamps in trash. Dispose as a hazardous waste. For information regarding safe handling, recycling and disposal, consult www.lamprecycle.org

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Chapter 4: Performance Features & Safety Precautions

DO NOT OVERLOAD THE WORK SURFACE WITH APPARATUS OR WORK MATERIAL. THE SAFE OPERATION OF THE LABORATORY HOOD IS BASED UPON HAVING PROPER AIRFLOW THROUGH THE STRUCTURE. DO NOT PLACE LARGE, BULKY OBJECTS SUCH AS BLOCK HEATERS, DIRECTLY ON THE HOOD WORK SURFACE. INSTEAD, ELEVATE THE OBJECT 2" TO 3" ON BLOCKS TO ALLOW A FLOW OF AIR UNDER THE OBJECT AND INTO THE LOWER REAR BAFFLE EXHAUST SLOT. ENSURE BLOCKS ARE LEVEL AND SECURED IN PLACE.

Blocking the bottom of the baffle at rear of hood will change the airflow pattern in the hood causing turbulence and possible leakage at the face of the hood. (Don't store containers or supplies against baffles, as this will affect airflow through the hood).

Avoid placing your head inside hood. Keep hands out of hood as much as possible.

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Always work as far back in hood as possible. It is best to keep all chemicals and apparatus 6" inside the front of the hood.

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This hood does not feature explosion-proof electrical components, unless ordered separately. Therefore, use of flammable or explosive materials in quantities above the explosive limit are not recommended.

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Do not work with chemicals in this hood without the exhaust system running. Do not store chemicals in a fume hood.

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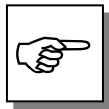
Perchloric acid use in this hood is prohibited.

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High level radioisotope materials are prohibited for usage in this hood.

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AVOID CROSS DRAFTS AND LIMIT TRAFFIC IN FRONT OF THE HOOD. AIR DISTURBANCES CREATED MAY DRAW FUMES OUT OF THE HOOD.

The use of heat-generating equipment in this hood without the exhaust system operating properly can cause damage to the hood.

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The Protector Laboratory Hood should be certified by a qualified certification technician before it is initially used. The hood should be re-certified whenever it is relocated, serviced or at least annually thereafter.

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Ensure that the unit is connected to electrical service in accordance with local and national electrical codes. Failure to do so may create a fire or electrical hazard. Do not remove or service any electrical components without first disconnecting the hood from electrical service.

Chapter 4: Performance Features & Safety Precautions

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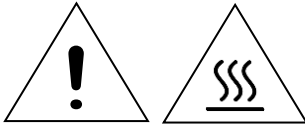
Proper operation of the fume hood depends largely upon the hood's location and the operator's work habits. Consult the Reference Manual in *Appendix E*.

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If the unit is not operated as specified in this manual it may impair the protection provided by the unit.

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Do not touch the blower motor. The surfaces of the motor can become hot and could cause burns.

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Do not position the fume hood so that it is difficult to operate the main disconnect device.

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To prevent the possibility of minor injury keep hands and fingers clear of sprockets at the top of the four corners.

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USING YOUR PROTECTOR PASS-THROUGH

S

Operating the Vertical-Rising Sash

Because of the Protector Hood counterbalanced sash mechanism, it will take only a few pounds of force to move the sash up or down, and you can operate the sash smoothly with one or two hands positioned any where along the handle. The vertical-rising sashes may be raised to a maximum 28" operating height. The airflow requirements should be sized for the 28" operating height; if using sash stops then the airflow requirements can be reduced by approximately 40% at 18". See *Chapter 2* for airflows.

S

Operating the Blower

Your Protector Fume Hood utilizes a remote style blower, which can be activated by turning the blower switch to "ON." You can validate the hood performance by watching smoke drawn into the hood face opening.

S

Operating the Lights

Your Protector Fume Hood utilizes a factory-wired fluorescent light to illuminate the hood interior. Optional direct replacement LED T8 bulbs are available per Appendix A. Simply turn the light switch to "ON" to operate.

Working in your Protector Fume Hood

Planning

- Thoroughly understand procedures and equipment required before beginning work.
- Arrange for minimal disruptions, such as room traffic or entry into the room while the hood is in use.

Start-up

Labconco follows the recommendations of the American National Standard Institute with regards to fume hood exhaust systems. A full list of requirements can be found in the ANSI Standard No. Z9.5 *Laboratory Ventilation*. If you are unsure of the proper use of the exhaust blower for your fume hood, contact your safety officer or industrial hygienist before making any modifications to the blower's operational status (on/off/low, etc.). Keeping the exhaust system under constant negative pressure is an excellent way to ensure hazardous fumes in the fume hood do not enter the laboratory space inadvertently.

- Turn on the light.
- Slowly raise the sash and verify the fume hood has proper airflow.
- Check the baffle air slots for obstructions.
- Allow the hood to operate unobstructed for two minutes.
- Wear a long sleeved lab coat and rubber gloves. Use protective eyewear. Wear a protective mask if appropriate.

Loading Materials and Equipment

- Only load the materials required for the procedure. Do not overload the hood.
- Do not obstruct the front air foil (sill), or rear baffle slots.
- Large objects should not be placed close together and spaced above the work surface to permit airflow to sweep under the equipment.
- After loading the hood, wait one minute to purge airborne contaminants from the work area.

Work Techniques

- Keep all materials at least 6 inches inside of the sash and at least two inches from the side walls and perform all contaminated operations as far to the rear of the work area as possible.
- Keep all clean and contaminated materials in the work area separate for clear identification.
- Avoid using techniques or procedures that disrupt the airflow patterns of the hood.

Final Purging

- Upon completion of the work, the hood should be allowed to operate for two minutes undisturbed, to purge airborne contaminants from the work area before closing the sash.

Unloading Materials and Equipment

- Objects in contact with contaminated material should be surface decontaminated before removal from the hood.
- All open trays or containers should be covered before being removed from the hood.

Shutdown

- Turn off the light and close the sash.

CHAPTER 6

MAINTAINING YOUR PROTECTOR PASS-THROUGH

Now that you have an understanding of how to work in the fume hood, we will review the suggested maintenance schedule and the common service operations necessary to maintain your fume hood for peak performance.

Service Safety Precautions

- If performing any electrical maintenance, always disconnect the power at the main disconnect.
- If performing decontamination inside the fume hood, consult your safety officer for proper personal protective equipment and procedure.
- Since some service operations require a step ladder, always use proper safety and consult your safety officer
- If performing maintenance on any service lines, always shut off the supply first.
- Some removable components may be heavy, follow safe-lifting guidelines.
- Verify all components are installed correctly with performance verified before conducting normal operations.

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Routine Maintenance Schedule

Weekly

- Using ordinary dish soap to clean the surface inside of the fume hood, and the work surface.
- Using an appropriate glass cleaner, clean the sash and all glass surfaces.
- Operate the fume hood blower, noting the airflow velocity through the hood using a source of visible smoke.

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Chapter 6: Maintaining Your Protector Pass-Through

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Biannually

- The sash assembly should be checked to ensure proper operation and to make sure there are no signs of abnormal wear on the sash pulleys, cables and clamps.

Routine Service Operations

Front Panel Removal:

1. Simply lift the front panel up and then away from the hood to provide access to the top.


Changing the Fluorescent Lamp or LED Direct Replacement Lamp:

1. Turn light switch to “OFF”.
2. Remove the front panel as noted earlier.
3. Remove the cover plate enclosing the fluorescent lamp (see Fig. 3-3). It is secured with six (6) acorn nuts, remove these nuts to remove the cover plate.
4. Remove the lamp by rotating it out of the two sockets.
5. Install the new lamp by reversing the removal procedure.

CHAPTER 7

MODIFYING YOUR PROTECTOR PASS-THROUGH

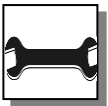There are several ways to modify the fume hood for your individual requirements. These include the addition of work surfaces, service fixtures, air monitor, electrical duplex outlets, and ceiling enclosures.

Installing Work Surfaces

Your Protector Fume Hood requires a work surface to work properly. Contact Labconco Customer Service for ordering information as listed.

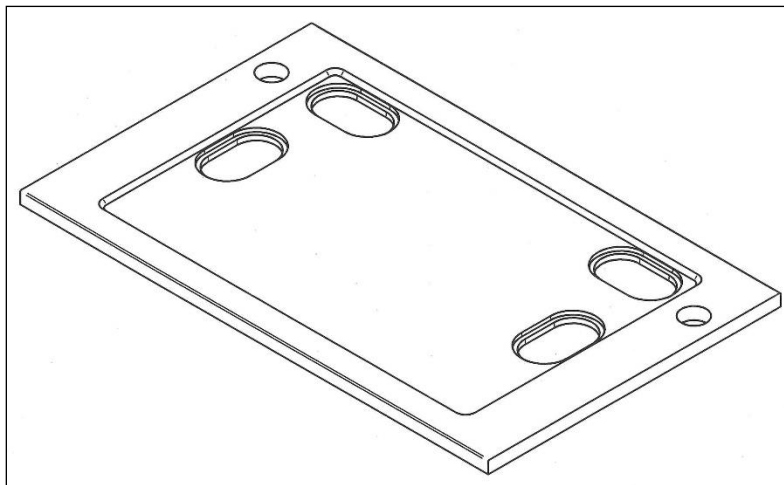
Figure 7-1

Ordering Information:

9570

4 – 4'	0 – None	0 – None
5 – 5'	1 – Rear	1 – Rear
6 – 6'	2 – Side	2 – Side
Width	Left	Right
	Cupsink	Cupsink
	Cutout	Cutout

Installing Ceiling Enclosures Above the Fume Hood

Your Protector Fume Hood has mounting holes to accept a ceiling enclosure to close off the area between the top of the hood and the ceiling. Contact Labconco Customer Service for ordering information. Ceiling Enclosures for Pass-Through Hoods are handled as a specials request.

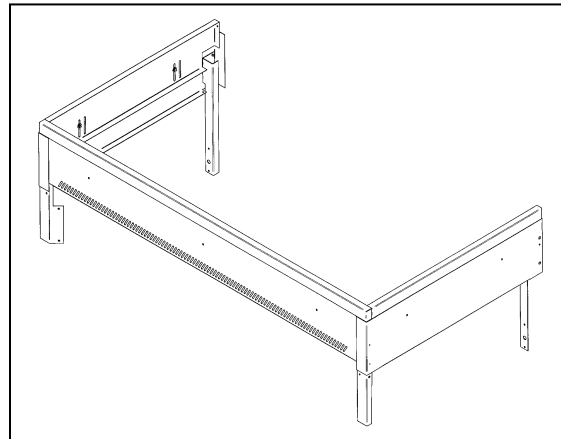
Figure 7-2

Installing Additional Service Fixtures

Additional service fixtures can be installed in the available service fixture holes in both sidewalls and corner posts. The fume hood is factory set to accept up to four valves per side. Contact Labconco Customer Service for information.

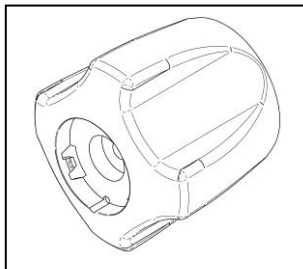
Figure 7-3 - Knob

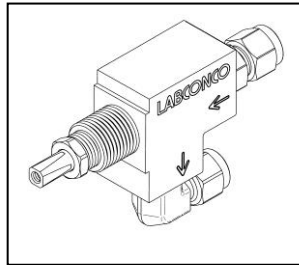
Figure 7-4 - Valve

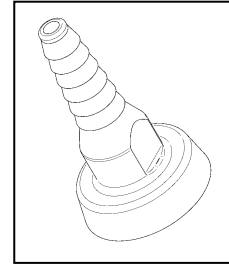
Figure 7-5 – Hose Connector

9827900	Cold Water Gooseneck, any position
9571000	Service Fixture CW, lower right, Pos. 1
9571001	Service Fixture CW, mid-lower right, Pos. 2
9571002	Service Fixture CW, mid-upper right, Pos. 3
9571003	Service Fixture CW, upper right, Pos. 4
9571004	Service Fixture CW, lower left, Pos. 1
9571005	Service Fixture CW, mid-lower left, Pos. 2
9571006	Service Fixture CW, mid-upper left, Pos. 3
9571007	Service Fixture CW, upper left, Pos. 4
9571100	Service Fixture GAS, lower right, Pos. 1

9571101	Service Fixture GAS, mid-lower right, Pos. 2
9571102	Service Fixture GAS, mid-upper right, Pos. 3
9571103	Service Fixture GAS, upper right, Pos. 4
9571104	Service Fixture GAS, lower left, Pos. 1
9571105	Service Fixture GAS, mid-lower left, Pos. 2
9571106	Service Fixture GAS, mid-upper left, Pos. 3
9571107	Service Fixture GAS, upper left, Pos. 4

Note: Any fixture can be converted to Air, Vacuum, etc. See *Appendix A* of this manual.

Installing Guardian™ Digital Airflow Monitor or Guardian™ Airflow Monitor

The Guardian Airflow Monitors continuously monitor face velocity through the fume hood opening. The fume hood right corner post is factory prepared to mount either monitor. Contact Labconco Customer Service to order.

Guardian Digital Airflow Monitors

9413400	100-115V, 50/60 Hz Digital
9413401	208-230V, 50/60 Hz Digital

Guardian Airflow Monitors

9413300	100-115V, 50/60 Hz Analog
9413301	208-230V, 50/60 Hz Analog

Note: Order the following for proper installation with Pass-Through Hoods

9571300	Kit, Guardian AFM Extension, Pass-Through
---------	---

Sash Stop Kit – Field Installation P/N 9410300

The sash stop kit restricts how far a vertical-rising sash may be opened. This small plastic device may be easily field installed on the fixture corner post of any fume hood.

Installing an Electrical Duplex Outlet

Your Protector Fume Hood can be ordered with duplex outlets, however, if you ordered a model without an electrical duplex outlet you can have one installed in the field by a qualified electrician. Contact Labconco Customer Service for ordering information. (Not acceptable on explosion-proof hoods). Order 9571200 for Pass-Through Hoods.

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CHAPTER 8

TROUBLESHOOTING

Refer to the following table if your fume hood fails to operate properly. If the suggested corrective actions do not solve your problem, contact Labconco for additional assistance.

PROBLEM	CAUSE	CORRECTIVE ACTION
Remote blower and lights won't operate	Wires not connected at junction boxes or switches.	Check connection of switches. Check connection to control box on top of unit.
	Circuit breakers tripped in building electrical supply.	Reset circuit breakers.
Remote blower won't operate, but lights work	Blower wiring is disconnected.	Inspect blower wiring and switch.
	Belt broken	Replace belt.
	Blower motor is defective.	Replace blower motor.
Fume hood blower operates but lights will not operate	Lamp not installed correctly.	Inspect lamp installation.
	Lamp is defective.	Replace lamp.
	Lamp circuit breaker in building is tripped.	Reset the lamp circuit breaker.

Chapter 8: Troubleshooting

PROBLEM	CAUSE	CORRECTIVE ACTION
Fume hood blower operates but lights will not operate	Lamp wiring is disconnected.	Inspect lamp wiring.
	Defective lamp ballast.	Replace lamp ballast.
Contaminants outside of fume hood	Improper user techniques for the fume hood.	See “Certifying the Hood” Chapter 3 and “Safety Precautions” Chapter 4 sections in the manual. (Ref. Appendix E)
	Restriction of the baffle air slots or – blockage of the exhaust outlet.	Remove baffles to ensure that all air slots, and the exhaust outlet are unobstructed.
	External factors are disrupting the fume hood airflow patterns or acting as a source of contamination.	See “Location Requirements” Chapter 2, “Certifying the Hood” Chapter 3, and “Safety Precautions” Chapter 4 sections of this manual. (Ref. Appendix E)
	Fume hood has improper face velocity.	Have fume hood re-certified and check remote blower exhaust system. Hood should have average face velocity of 60-100 fpm.
Vertical sash no longer operates smoothly	Cable is frayed or plastic protection is damaged.	Inspect cable and replace cable if worn or damaged immediately; otherwise injury could result.
	Pulley bearing is damaged.	Replace pulley, bearing or add grease.
	Cable has slipped off the pulleys.	Re-install, cable must be replaced immediately if damaged.
	Weight has broken pulleys.	Replace pulley(s).
	Vertical sash frame is distorted.	Place horizontal glass symmetrically and pull sash down to air foil. Straighten damaged frame.

PROBLEM	CAUSE	CORRECTIVE ACTION
	Cable is frayed or has slipped off the pulleys.	Re-install, cable must be replaced immediately if damaged.
Electrical duplex outlets no longer have power	Wires not connected or faulty duplex.	Check wire connection or replace duplex.
	Circuit breakers tripped in building electrical supply.	Reset circuit breakers.
Service valves no longer operate	Faulty building supply.	Inspect building supply shut off valves and appropriate pressures below the recommended 40 PSI.
	Valve no longer operates.	Replace valve and check for leaks.
	Supply line or outlet line has leaks.	Inspect line for leaks and fix any leaking plumbing connections.
Audible Alarm Sounding	Both sashes are too high.	Close one sash completely.

APPENDIX A

PROTECTOR PASS-THROUGH COMPONENTS

Illustration A-1 indicate the location of the following service parts:

Protector Pass-Through Replacement Parts

Item	Quantity	Part No.	Description
1A	1	9823700	Valve, Labconco (Water) 1/4" Compression Fitting
1B	1	9823701	Valve, Labconco (Water) 3/8" Compression Fitting
1C	1	9817000	Valve, Labconco 1/4" Compression Fitting (AIR, GAS, VAC, NIT, etc.)
1D	1	9817001	Valve, Labconco 3/8" Compression Fitting (AIR, GAS, VAC, NIT, etc.)
1E	1	9823702	Valve, Labconco Deionized 1/4" Compression Fitting
1F	1	9823703	Valve, Labconco Deionized 3/8" Compression Fitting
1G	1	9818000	Nut, Valve Mounting (Labconco)
2A	1	9826800	WaterSaver Valve/Gooseneck – GRN
2B	1	9826801	WaterSaver Valve/Connector (VAC) – YEL
2C	1	9826802	WaterSaver Valve/Connector (AIR) – ORG
2D	1	9826803	WaterSaver Valve/Connector (GAS) – BLU
2E	1	9826805	WaterSaver Valve/Connector (HOT WATER) – RED
2F	1	9826806	WaterSaver Valve/Connector (CW) – GRN
2G	1	9826807	WaterSaver Valve/Connector (STEAM) – BLK
2H	1	9826808	WaterSaver Valve/Connector (NITROGEN) – BRN
2I	1	9826809	WaterSaver Valve/Connector (OXYGEN) – LIGHT GREEN
2J	1	9826810	Swivel Gooseneck only – GRN
2K	1	9826812	Swivel Gooseneck only – WHITE
3	1	9818700 thru 08	Knobs (GRAY, GRN, BLU, ORG, YEL, RED, WHT, BLK, BRN)
4A	1	9818800	Hose Barb, GRAY (NEUTRAL OR ARGON) – NOT SHOWN
4B	1	9818801	Hose Barb, GREEN (COLD WATER) – NOT SHOWN
4C	1	9818802	Hose Barb, BLUE (GAS) – NOT SHOWN
4D	1	9818803	Hose Barb, ORANGE (AIR) – NOT SHOWN
4E	1	9818804	Hose Barb, YELLOW (VACUUM) – NOT SHOWN
4F	1	9818805	Hose Barb, RED (HOT WATER) – NOT SHOWN
4G	1	9818806	Hose Barb, WHITE (DEIONIZED WATER) – NOT SHOWN
4H	1	9818807	Hose Barb, BLACK (NEUTRAL OR STEAM) – NOT SHOWN
4I	1	9818808	Hose Barb, BROWN (NITROGEN) – NOT SHOWN
4J	1	9819000	Nut, Hose Barb – NOT SHOWN
5	1	9825500	Label, Knob (contains all the labels)
6	1	9818900	Lens, Knob
7A	1	9947100, 01	115V Duplex Receptacle (GRAY) Right (00), Left (01) w/ wires
7B	1	9818200	Cover Plate 115V Duplex
7C	1	9947103, 04	115V GFCI Duplex Receptacle (GRAY) Right (03), Left (04) w/ wires
7D	1	9818100	Cover Plate, 115V GFCI
7E	1	9818300	Cover Plate, Blank

Appendix A: Protector Pass-Through Components

Item	Quantity	Part No.	Description
8A	2	9721902	Lamp, Fluorescent (T8 x 2') – use on 4' Hoods – NOT SHOWN
8B	2	9721901	Lamp, Fluorescent (T8 x 3') – use on 5' Hoods – NOT SHOWN
8C	2	9721900	Lamp, Fluorescent (T8 x 4') – use on 6' Hoods – NOT SHOWN
8D	1	9721911	LED, Direct Replacement T8 x 3' – use on 5' Hoods
8E	1	9721910	LED, Direct Replacement T8 x 4' – use on 6' Hoods
8F	1	9721912	LED, Direct Replacement T8 x 2' – use on 4' Hoods
9	1	3738300	Limit Switch (Sash Position alarm)
10A	1	1302301	Switch, Rocker
10B	1	1327500	Switch, Plug (Fills cutout when switch is not used)
11	1	9818400	Access Cover
12A	1	9568300	Side Panel, no window
12B	1	9568301	Side Panel, 30", side view window
12D	4	1885308	Screw, Machine #10-24 x .50 Phillips
13A	1	9436501	Eco-Foil 4'
13B	1	9436502	Eco-Foil 5'
13C	1	9436503	Eco-Foil 6'
13D	1	9436521	PVC Eco-Foil, 4'
13E	1	9436522	PVC Eco-Foil, 5'
13F	1	9436523	PVC Eco-Foil, 6'
14	4	1861400 or 1850000	Pulley, Front, 1-3/16 Dia. or Nylon 2"
15A	2	4949901	Cable, Sash 80"
15B	2	4949904	Cable, Sash 190"
16	4	9742000	Pulley, Rear, 2-3/4" Dia.
17	2	9713300	Bumper, Rubber (lower sash bumper) – NOT SHOWN
18	8	1920100	Clamp, Cable (Replacement)
19	2	9935800	Threaded Connector (to attach weight to cable) – NOT SHOWN
20	1	9569500 or 01	Wiring Harness, Main (115v or 230v) - NOT SHOWN
21	1	9572500	Wiring Harness, Ballast (same all voltages) - NOT SHOWN
22	1	9409601, 02, 03	Front Panel, 4', 5', 6'
23	1	1294000	Ballast (115-277v, 50/60 Hz)
24A	1	9413300, 01	Guardian Airflow Monitor (115v or 230v) - NOT SHOWN
24B	1	9571300	Guardian Airflow Monitor Extension Kit (Required on Pass-Through)
25	1	9413400, 01	Guardian Digital Airflow Monitor (115v or 230v) - NOT SHOWN
26A	1	1328102	Sound, Alarm 115V
26B	1	1328101	Sound, Alarm 230V
27	2	3738300	Switch, Trip Alarm

Appendix A: Protector Pass-Through Components

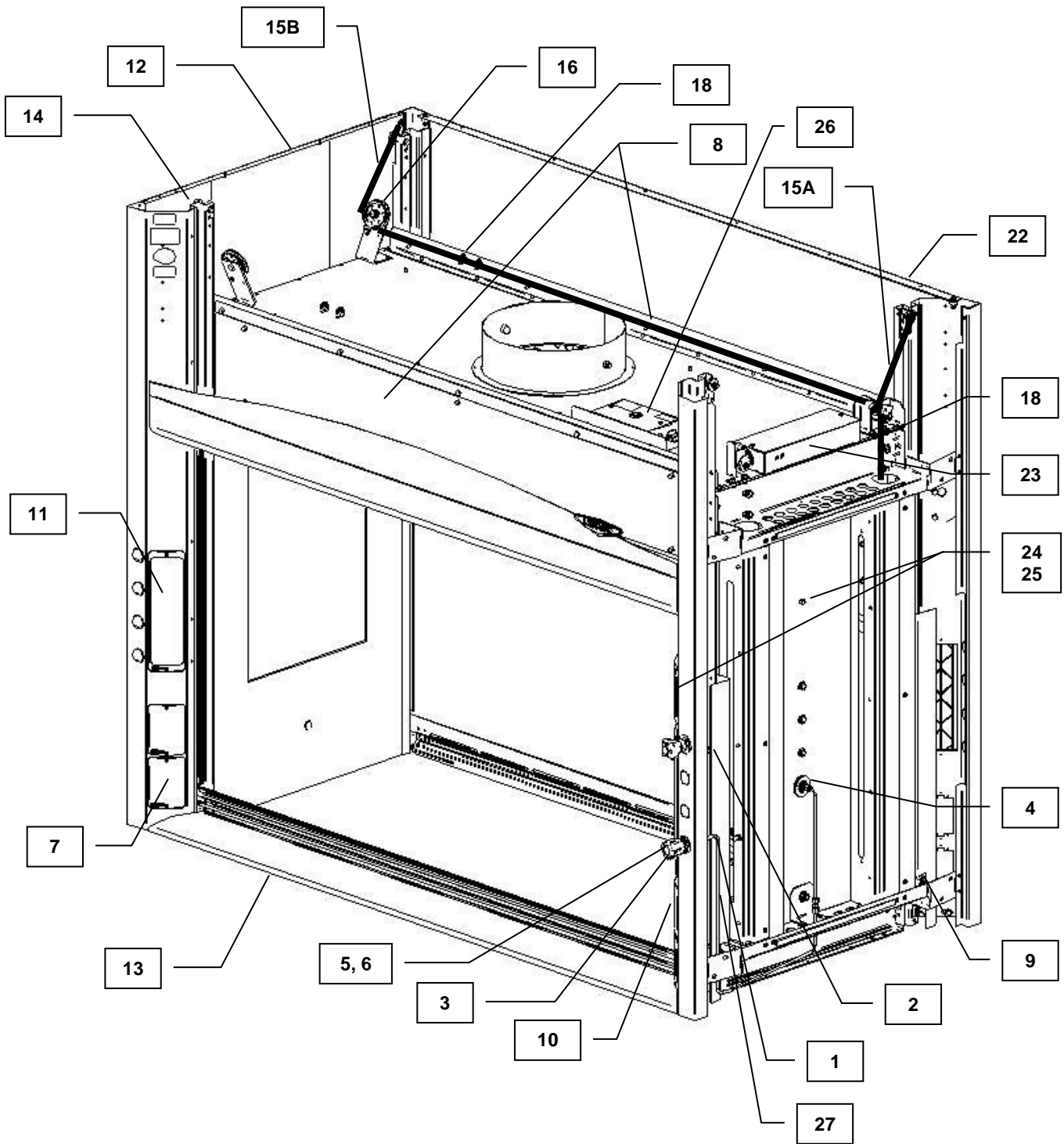
Figure A-1

APPENDIX C

PROTECTOR PASS-THROUGH

SPECIFICATIONS


Environmental Conditions

- Indoor use only.
- Maximum altitude: 10,000 feet (3,048 meters).
- Ambient temperature range: 41° to 104°F (5° to 40°C).
- Maximum relative humidity: 80% for temperatures up to 88°F (31°C), decreasing linearly to 50% relative humidity at 104°F (40°C).
- Main supply voltage fluctuations not to exceed $\pm 10\%$ of the nominal voltage.
- Transient over-voltages according to Installation Categories II (Over-voltage Categories per IEC 1010). Temporary voltage spikes on the AC input line that may be as high as 1500V for 115V models and 2500V for 230V models are allowed.
- Used in an environment of Pollution degrees 2 (i.e., where normally only non-conductive atmospheres are present). Occasionally, however, a temporary conductivity caused by condensation must be expected, in accordance with IEC 664.
- Electrical Ratings

<u>Volts AC</u>	<u>Phase</u>	<u>Cycle</u>	<u>AMP*</u>
115	1	50/60	10
115	1	50	10
115	1	60	10
230	1	50/60	5
230	1	50	5
230	1	60	5

*Does not include current rating of receptacles.

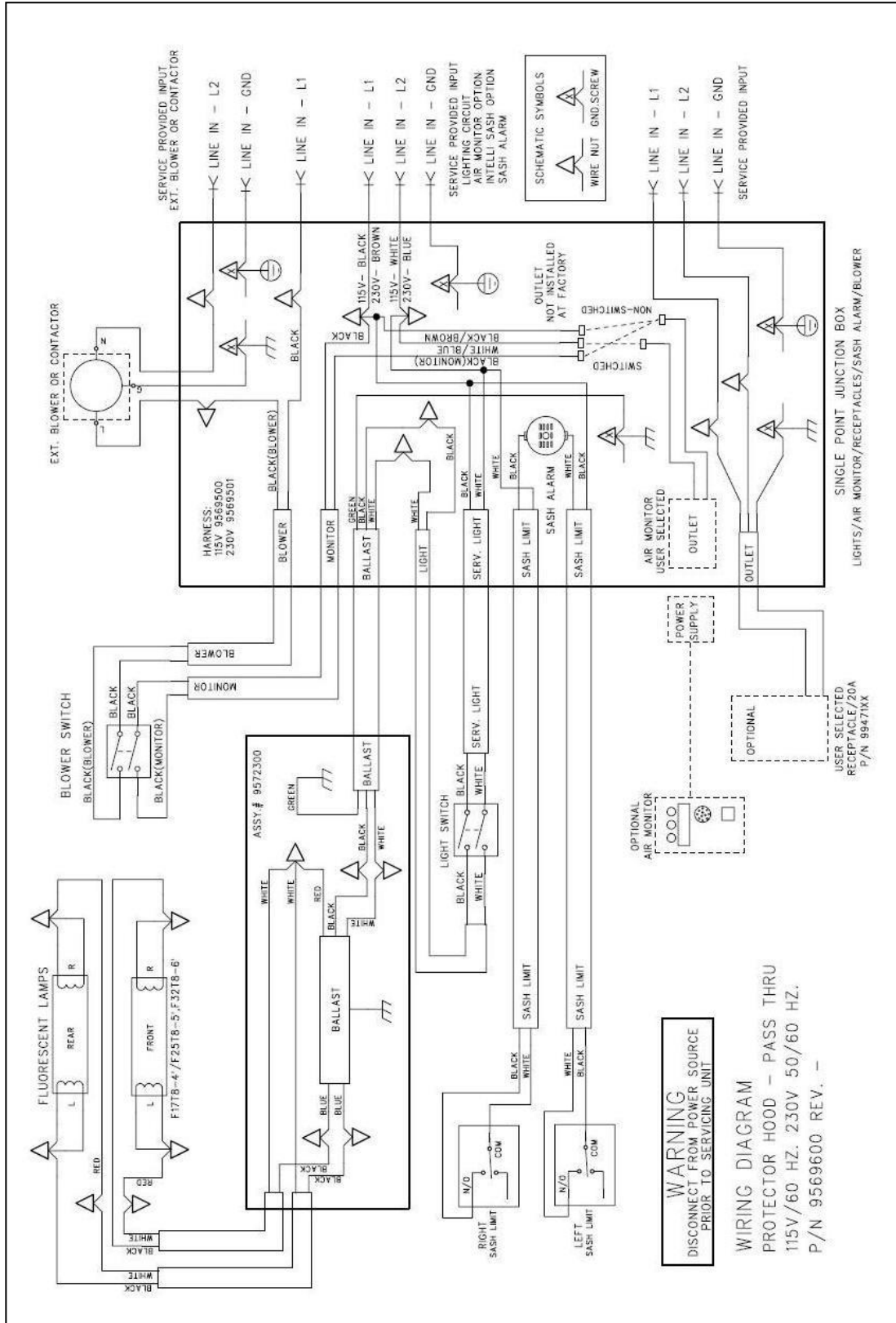Appendix C: Protector Pass-Through Specifications

APPENDIX D

SERIAL NUMBER TAG

DESCRIPTION

Serial tag includes standard information with the following changes:

The portion of the tag for electrical information is labeled with the following units: Volts AC, Phase, Cycle, AMP.

Each hood will include one of the following:

<u>Volts AC</u>	<u>Phase</u>	<u>Cycle</u>	<u>AM!</u> _x
115	1	50/60	10X
115	1	50	10X
115	1	60	10X
230	1	50/60	5X
230	1	50	5X
230	1	60	5X

Where  may be a separate label indicating: Caution, See Manual.

The X above will be one of several possible alpha characters as defined in the manual as follows:

- X: No additional duplexes on hood, noted amperage applies.
- B: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- C: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.

- D: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- E: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- F: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- G: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- H: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- J: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- K: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- L: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.

Appendix D: Serial Number Tag Description

- M: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- N: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.

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APPENDIX E

REFERENCES

Many excellent reference texts and booklets are currently available. The following is a brief listing:

Laboratory Ventilation Standards

Federal Register 29 CFR Part 1910

Non-mandatory recommendations from "Prudent Practices".

- Fume hoods should have a continuous monitoring device
- Face velocities should be between 60-100 linear feet per minute (lfpm)
- Average 2.5 linear feet of hood space per person

Occupational Health and Safety

U.S. Department of Labor

200 Constitution Avenue N.W.

Washington, DC 20210

(202) 523-1452

Industrial Ventilation-ACGIH

- Fume hood face velocities between 60-100 lfpm
- Maximum of 125 lfpm for radioisotope hoods
- Duct velocities of 1000-2000 fpm for vapors, gasses and smoke
- Stack discharge height 1.3-2.0 x building height
- Well designed fume hood containment loss, <0.10 ppm

Industrial Ventilation, A Manual of Recommended Practice.

24th Edition, 2001

American Conference of Governmental Industrial Hygienists

1330 Kemper Meadow drive

Cincinnati, OH 45240-1634

(513) 742-2020

ASHRAE 110-1995 Method of Testing Performance of Fume Hoods

Evaluates fume hood's containment characteristics

- Three part test: Smoke generation, Face velocity profile, Tracer gas release @ 4 liters per minute
- Rated As Manufactured (AM), As Installed (AI) and As Used (AU)

American Society of Heating, Refrigerating, and Air Conditioning Engineers

1791 Tullie Circle N.E.

Atlanta, GA 30329

(404) 636-8400

ANSI Z9.5-2011 Laboratory Standard

Covers entire laboratory ventilation system.

- New and remodeled hoods shall have a monitoring device
- Ductless hoods should only be used with non-hazardous materials

American Industrial Hygiene Association

2700 Prosperity Avenue, Suite 250

Fairfax, VA 22031

(703) 849-8888

SEFA 1-2002

- Fume hood face velocities based on toxicity levels of chemicals
 - Class A – 125 to 150 fpm
 - Class B – 80 to 100 fpm
 - Class C – 75 to 80 fpm
- Test method – face velocity profile and smoke generation

Scientific Equipment & Furniture Association

1028 Duchess Drive

McLean, VA 22102

(703) 538-6007

NFPA 45 – 2002 Fire Protection for Laboratories Using Chemicals

- Laboratory hoods should not be relied on for explosion protection
- Exhaust air from fume hoods should not be recirculated
- Services should be external to the hood
- Canopy hoods only for non-hazardous applications
- Materials of construction should have flame spread of 25 or less
- 80 to 120 fpm to prevent escape

NFPA 30 – 2000 Flammable and Combustible Liquids Code

- Approved cabinets may be metal or wood
- Vent location on cabinets are required
- Venting of cabinets not a requirement

National Fire Protection Association

1 Batterymarch Park

P.O. Box 9101

Quincy, MA 02269-9101

(800) 344-3555

General References

American Conference of Governmental Industrial Hygienists. *Industrial Ventilation, A Manual of Recommended Practice*, Cincinnati, OH

ASHRAE Standard Committee. *ASHRAE Standard* Atlanta: ASHRAE Publications Sales Department, 1995

British Standards Institution, *Laboratory Fume Cupboards*. Parts 1, 2 and 3, London: 1990

Department of Labor, Occupational Safety and Health Administration, *29 CFR Part 1910, Occupational Exposures to Hazardous Chemicals in Laboratories, Final Rule*. Vol. 55, No. 21. Washington D.C.:1990

Appendix E: References

DiBerardinis, L. et al. *Guides for Laboratory Design, Health and Safety Considerations*. Wiley & Sons, 1987

McDermott, Henry, *Handbook of Ventilation for Contaminant Control, 2nd Edition*. Butterworth Publishers, 1985.

Miller, Brinton M. et al. *Laboratory Safety: Principles and Practices*. American Society for Microbiology, Washington, D.C.: 1986

NIH Guidelines for the Laboratory Use of Chemical Carcinogens. NIH Publication No. 81-2385.

Rayburn, Stephen R. *The Foundation of Laboratory Safety, A Guide for the Biomedical Laboratory*. Springer-Verlag, New York: 1990

Sax, N. Irving and Lewis, JR., Richard J. *Rapid Guide to Hazardous Chemicals in the Workplace*. Van Nostrand Reinhold, 1987.

Schilt, Alfred A. *Perchloric Acid and Perchlorates*. The G. Frederick Smith Chemical Company, Columbus, OH: 1979.

Steere, Norman. *CRC Handbook of Laboratory Safety, 2nd Edition*. CRC Press, 1971.

APPENDIX F

DISASSEMBLY AND REASSEMBLY INSTRUCTIONS

Required Tools

The following tools will be required for disassembly and reassembly:

Cordless Drills and Phillips bits.

1/4", 5/16", 3/8", 7/16" Sockets	Small Screwdriver
Phillips Screwdriver	Right Angle Screwdriver
1/4", 5/16" Box Ratchet	1/4" Extension
5/16", 3/8", 7/16" Combination Wrench	1/4" Ratchet
Small Right Angle Needle Nose Pliers	Nut Driver Set

Decorative Panel Removal

1. Remove side panels by removing #10-24 screws and lifting away.
2. Remove front panels by lifting up and away.

Blower Wiring

1. All the removable wires are part of the wiring harness to the corner post switches. Be sure to label all wires and refer to the wiring diagrams for proper reassembly.

Header Removal

1. The header panels are fastened to the corner posts by four #12 screws each. Simply reach behind the corner posts from the side and remove the screws while another person supports the header.

Appendix F: Disassembly & Reassembly Instructions

Sash Removal

There are two (2) sashes to be removed on the Pass-Through hoods. Please repeat the procedure for each sash.

1. Remove the sash end stop brackets by removing the two (2) #10-24 screws on each end stop bracket. The sash cables are attached through these end stop brackets.
2. Now the glass sash assembly can be removed by sliding it up and out.

Note - Sash weights are mounted in retention tracks on one side of the Pass-Through hood. They do not need to be removed to remove the sash(es). If sash weight removal is desired, see *Chapter 3: Sash Weight Release*.

Airfoil Removal

1. Remove lower corner post screw located under the airfoil.
2. Remove the airfoil by removing the rear screws located on top of the airfoil on each side.

Transport and Reassembly

The various subassemblies are now ready to be transported and reassembled in the opposite way the hood was disassembled. All hoods will require a minimum of five people to lift the hood, two people on each side and one in front (for balance only). A hydraulic lift table should be used as an aid in lifting. If in doubt about lifting the hood, always reassemble the hood where it will be placed.

Sash Weight Reassembly

After the hood is assembled, the sash weights are easily installed before the hood is installed in a wall, because the wall may reduce access to the side panel and the sash weights behind the side panel. **Please keep this in mind and install the sash weights now.**

Hood Replacement and Service Installation

The hood is now ready for placement and the HVAC, plumbing and electrical service can be installed.