

Instruction Manual

HI 5221 & HI 5222 pH/mV/ISE/Temperature Bench Meters



www.hannainst.com

Dear Customer,

Thank you for choosing a Hanna Instruments product.

Please read this instruction manual carefully before using these instruments. This manual will provide you with the necessary information for correct use of these instruments, as well as a precise idea of their versatility.

If you need additional technical information, do not hesitate to e-mail us at tech@hannainst.com or view our worldwide contact list at www.hannainst.com.

WARRANTY

HI 5221 and **HI 5222** are guaranteed for two years against defects in workmanship and materials when used for their intended purpose and maintained according to instructions. Electrodes and probes are guaranteed for six months. This warranty is limited to repair or replacement free of charge.

Damage due to accidents, misuse, tampering or lack of prescribed maintenance is not covered.

If service is required, contact the dealer from whom you purchased the instrument. If under warranty, report the model number, date of purchase, serial number and the nature of the problem. If the repair is not covered by the warranty, you will be notified of the charges incurred. If the instrument is to be returned to Hanna Instruments, first obtain a Returned Goods Authorization number from the Technical Service department and then send it with shipping costs prepaid. When shipping any instrument, make sure it is properly packed for complete protection.

To validate your warranty, fill out and return the enclosed warranty card within 14 days from the date of purchase.

All rights are reserved. Reproduction in whole or in part is prohibited without the written consent of the copyright owner.

TABLE OF CONTENTS

WARRANTY	2
PRELIMINARY EXAMINATION	3
GENERAL DESCRIPTION	4
FUNCTIONAL DESCRIPTION	5
SPECIFICATIONS	8
OPERATIONAL GUIDE	9
DISPLAYING MODES	11
SYSTEM SETUP	14
pH SETUP	21
mV SETUP	34
ISE SETUP (HI 5222 only)	36
pH CALIBRATION	44
pH MEASUREMENT	47
mV & Relative mV MEASUREMENTS	49
ISE CALIBRATION (HI 5222 only)	51
ISE MEASUREMENT (HI 5222 only)	54
TEMPERATURE CALIBRATION	60
LOGGING	61
PC INTERFACE	64
ADDITIONAL INFORMATION	65
pH BUFFER TEMPERATURE DEPENDENCE	68
ELECTRODE CONDITIONING AND MAINTENANCE	69
TROUBLESHOOTING GUIDE	71
TEMPERATURE CORRELATION FOR pH SENSITIVE GLASS	72
ACCESSORIES	73

PRELIMINARY EXAMINATION

Remove the instrument from the packing material and examine it carefully to make sure that no damage has occurred during shipping. If there is any damage, notify your Dealer or the nearest Hanna Customer Service Center.

The meters are supplied complete with:

- **HI 1131B** Glass body Combination pH Electrode
- **HI 7662-W** Temperature probe
- **HI 7082S** Electrolyte Solution
- **HI 76404W** Electrode Holder
- pH Calibration Solutions Kit
- One capillary dropper pipette
- 12 Vdc Power Adapter
- Instruction Manual

HI 5221-01 and **HI 5222-01** are supplied with 12 Vdc/115 Vac adapter.

HI 5221-02 and **HI 5222-02** are supplied with 12 Vdc/230 Vac adapter.

Note: Save all packing material until you are sure that the instrument works properly. Any defective item must be returned in the original packing with the supplied accessories.

GENERAL DESCRIPTION

HI 5221 and **HI 5222** are professional bench meters with a color graphic LCD for pH (with Calibration Check), ORP (Oxidation Reduction Potential), ISE (**HI 5222** only) and Temperature measurements.

The display can be configured as a single channel or dual channel display (**HI 5222** only). Each channel can be configured as pH, mV, Relative mV or ISE (**HI 5222** only).

The main features of the instruments are:

- Single (**HI 5221**) or two (**HI 5222**) input channels;
- Capacitive touch keypad;
- pH calibration using up to five pH buffers: choose from the list of Hanna buffers (1.68, 3.00, 4.01, 6.86, 7.01, 9.18, 10.01 and 12.45) or custom buffers (up to 5 custom buffers);
- ISE calibration using up to five standards: choose from the list of nominal standard values (e.g. for ppm: 0.010, 0.100, 1.00, 10.0, 100, 1000, 10000 ppm) and/or from 5 user-supplied custom standard values (up to 5 custom solutions);
- AutoHold feature to freeze first stable reading on the LCD;
- Two selectable alarm limits;
- User Selectable logging modes: Automatic logging, Manual logging with or without AutoHold;
- Continuous Lot logging directly on meter, with selectable log interval: Store up to 100,000 total data points;
- Up to 100 logging lots; a single lot can have up to 50,000 data points; **HI 5222** (same for each measurement channel). Log files include GLP information;
- Basic Measurement can be viewed with detailed GLP information, or with a Graph or a Log History;
- User-friendly interface on large color graphic LCD (240 x 320 pixels) with user selectable color palette;
- PC interface via USB; download logged data to PC or use for Real time logging (**HI 92000** PC application required);

- Dedicated Help key with contextual message;
- Profile feature: store up to ten different user setup (HI 5222 - five on each channel).


FUNCTIONAL DESCRIPTION

HI 5221 DESCRIPTION

FRONT PANEL



REAR PANEL

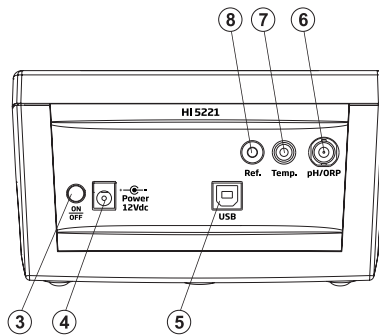


- 1) Liquid Crystal Display (LCD)
- 2) Capacitive touch keypad
- 3) ON/OFF switch
- 4) Power adapter socket
- 5) USB connector
- 6) BNC electrode connector for pH/ORP measurements
- 7) Temperature probe socket
- 8) Reference input socket

HI 5222 DESCRIPTION

FRONT PANEL



REAR PANEL

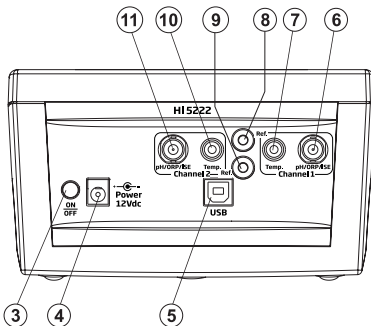

- 1) Liquid Crystal Display (LCD)
- 2) Capacitive touch keypad
- 3) ON/OFF switch
- 4) Power adapter socket
- 5) USB connector
- 6) BNC electrode connector for pH/ORP/ISE measurements (Channel 1)
- 7) Temperature probe socket (Channel 1)
- 8) Reference input socket (Channel 1)
- 9) Reference input socket (Channel 2)
- 10) Temperature probe socket (Channel 2)
- 11) BNC electrode connector for pH/ORP/ISE measurements (Channel 2)

KEYBOARD DESCRIPTION


FUNCTION KEYS

- CAL** To enter/exit calibration mode.
- MODE** To select the desired measurement mode, pH, mV, Rel mV (or ISE – HI 5222 only).
- SETUP** To enter Setup (System, pH, mV or ISE) and to access Log Recall function.
- HELP** To obtain general information about the selected option/operation.

VIRTUAL KEYS

The upper row keys are assigned to the **virtual keys** placed on the bottom of the LCD, which allow you to perform the displayed function, depending on the current menu (e.g. **Display**, **Start Log 1** and **Channel** in *Measure* mode).

Note: All the virtual keys are assigned to the highlighted channel (HI 5222 only).

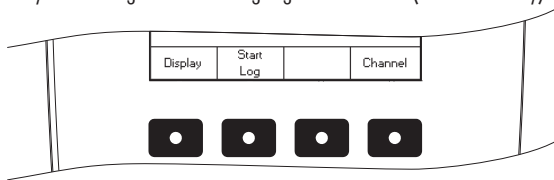

LCD GENERAL DESCRIPTION

The diagram shows a detailed view of the LCD screen with various elements labeled:

- Time & Date:** 01:37:49 PM, Dec 15, 2014
- Selected mode:** Measure
- Stability indicator:** Stable
- Measured value:** 10.015
- Measurement unit:** pH
- Short GLP information:** Last Cal.: Dec 15, 2014 01:37 PM, Electrode Cond.: 100%
- Temperature probe measurement status:** ATC2
- Temperature value & unit:** 24.5°C
- Reminder messages area:** (Empty area)
- Virtual keys:** Display, Start Log, Channel

SPECIFICATIONS

	HI 5221	HI 5222
Range	-2.0 to 20.0 pH / -2.00 to 20.00 pH / -2.000 to 20.000 pH	
	± 2000.0 mV	
	-	e.g. 10^{-7} to 10 M, 0.005 to 10^5 ppm, $5 \cdot 10^{-7}$ to $5 \cdot 10^7$ conc.
	-20.0 to 120.0 °C / -4.0 to 248.0 °F / 253.15 to 393.15 K	
Resolution	0.1 pH / 0.01 pH / 0.001 pH	
	0.1 mV	
	-	1, 2, 3 significant digits
	0.1 °C / 0.1 °F / 0.1 K	
Accuracy @25 °C / 77 °F	± 0.1 pH / ± 0.01 pH / ± 0.002 pH ± 1 last significant digit	
	± 0.2 mV ± 1 last significant digit	
	-	$\pm 0.5\%$ (monovalent ions) $\pm 1\%$ (divalent ions)
	± 0.2 °C / ± 0.4 °F / ± 0.2 K	
Relative mV offset range	± 2000.0 mV	
pH Calibration	Up to five points calibration, 8 standard buffers available (1.68, 3.00, 4.01, 6.86, 7.01, 9.18, 10.01, 12.45), and 5 custom buffers	
ISE Calibration	-	Up to five points calibration, seven fixed standard solutions available, for each measurement unit, and five custom solutions
Temperature User Calibration	3 points (0, 50, 100 °C)	
Input Channels	1 pH/mV	2 pH/mV/ISE
Temperature Compensation	Manual or Automatic	
GLP	Electrode offset / slope, calibration points, calibration time stamp	
Logging Record	Up to 100 lots, 50,000 records max/lot / maximum 100,000 data points/channel	
Logging Interval	14 selectable between 1 second and 180 minutes	
Logging Type	Automatic, Log on demand and Auto Hold	
pH Electrode	HI 1131B	
Temperature Probe	HI 7662-W	
PC Interface	Opto-isolated USB	
Input Impedance	10^{12} ohms	
Display	Color Graphic LCD (240 x 320 pixels)	
Power Supply	12 Vdc adapter	
Environment	0 - 50 °C / 32 - 122 °F / 273 - 323 K max. 95% RH non-condensing	
Dimensions	160 x 231 x 94 mm (6.3 x 9.1 x 3.7")	
Weight	1.2 Kg (2.6 lb.)	

OPERATIONAL GUIDE

POWER CONNECTION

Plug the 12 Vdc adapter into the power supply socket.

- Notes:**
- These instruments use non-volatile memory to retain the pH, ISE calibrations and all other settings, even when unplugged.
 - Make sure a fuse protects the main line.

ELECTRODE AND PROBE CONNECTIONS

For pH or ORP measurements, connect a pH/ORP electrode with internal reference to the BNC connector located on the rear panel of the instrument (to the desired channel - **HI 5222** only).

For ISE measurements, (**HI 5222** only) connect, to the desired channel, an ISE electrode with internal reference to the BNC connector located on the rear panel of the instrument.

For electrodes with a separate reference, connect the electrode's BNC to the BNC connector and the electrode's reference to the reference input socket.

For temperature measurement and automatic temperature compensation, connect the temperature probe to the appropriate socket (to desired channel - **HI 5222** only).

INSTRUMENT START UP

- Please ensure that the capacitive keypad is not covered by hand or other objects at the meter power on.
- Turn the instrument on from the power button located on the rear panel of the instrument.
- Please wait until the instrument finishes the initialization process.

Note: It is normal for the loading process to take a few seconds. If the instrument doesn't display the next screen, restart the meter using the power button. If the problem persists, contact your dealer.



CHANNEL SELECTION (HI 5222 only)

- Press **Channel** while in *Measure* mode to access channel selection menu. Four available options will be displayed: Channel 1, Channel 2, or multi-channel with the first or the second channel focused. The “**Choose Channel Configuration**” message is displayed in the Reminder messages area.
- Select the desired option by pressing the appropriate key: **Channel 1**, **Channel 2**, **Channel 1 Channel 2** or **Channel 1 Channel 2**. The instrument will display in the selected option *Measure* mode.

11:23:44 AM Dec 15, 2014		Measure	
Channel 1			
9.66		Stable E-3 ppm	
Last Cal.: Dec 15, 2014 11:12 AM ISE: Fluoride		TEMP1 24.8°C	
Channel 2			
7.090		Stable pH	
Last Cal.: Dec 15, 2014 11:22 AM Electrode Cond.: 100%		ATC2 25.2°C	
Choose Channel Configuration			
Channel 1	Channel 2	Channel 1 Channel 2	Channel 1 Channel 2

DISPLAYING MODES



For each measurement mode (pH, mV, Rel mV or ISE) the following display configurations are available: Basic, Good Laboratory Practice (GLP) (pH, ISE only), Graph and Log History.

Basic

The main measured value and its units are displayed on the LCD, along with the temperature value, temperature probe status and basic calibration information when this option is accessed.

To choose the Basic display mode:

- Press Display while in *Measure* mode. The “Choose Display Configuration” message will be displayed in the Reminder messages area.
- Press Basic. The instrument will display the basic information for the selected measurement mode.

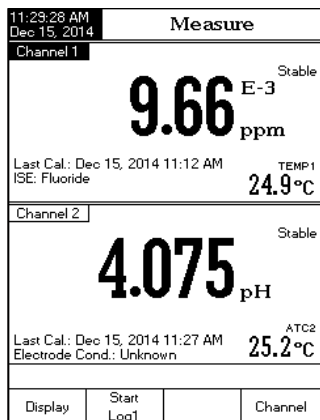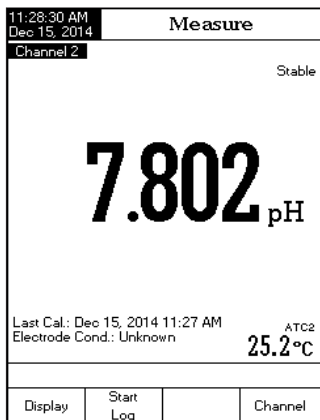



GLP

Detailed GLP data will be displayed on the LCD for pH Measure and ISE Measure modes only when this option is selected: Last Calibration date and time, Offset and Slope values, Calibration Buffers/Standards and general information regarding the buffers/standards: the calibration temperature, temperature compensation mode, date and time. For *pH Measure*, the Electrode Condition is also displayed on the LCD in percent.

Note: If a one point pH calibration is performed or the current calibration does not include at least two consecutive standard buffers of pH 4.01, 7.01 (6.86) and 10.01 (9.18) buffers, the Electrode Condition will be unknown. Electrode condition remains active for 24 hours after a calibration.

To access the GLP display option:

- Press Display while in *Measure* mode. The “Choose Display Configuration” message will be displayed in the Reminder messages area.
- Press GLP. The instrument will display the detailed GLP data.

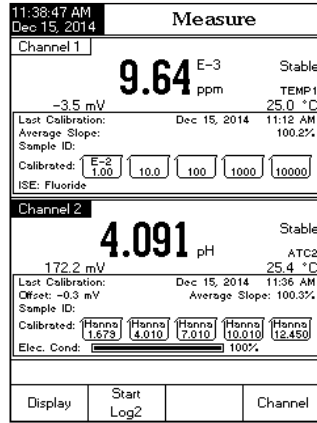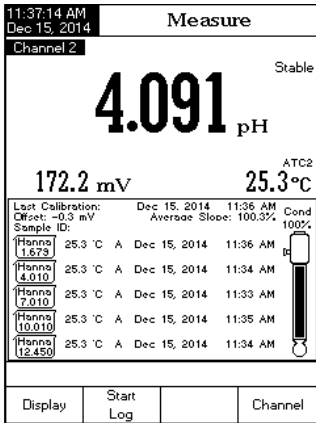



Graph

The on-line graph with currently logged values (pH, mV, Rel mV, or ISE vs. Seconds) will be displayed when this option is selected.

If there is no active log, the previously logged data for the selected parameter will be shown.

To access the off-line / on-line graph:

- Press **Display** while in *Measure / Logging* mode. The “Choose Display Configuration” message will be displayed in the Reminder messages area.
- Press **Graph**. Press **Start Log** to begin an online graph.

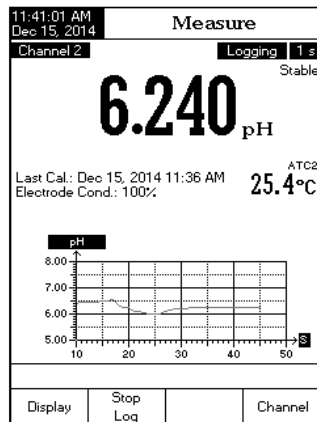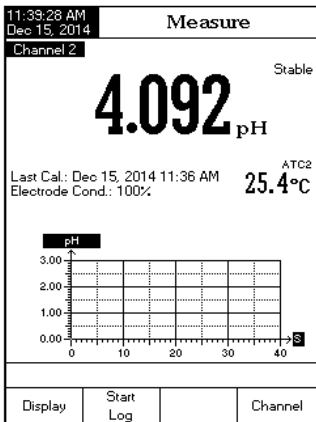

To Zoom Graph

- Press **Display** then **Graph**, **<** and **<** will appear in virtual keys.
- Use **<** and **>** to move along the X (Time) axis of the graph.
- Press **SETUP** to access the zoom menu for Y axis. Use **Zoom IN** or **Zoom OUT** for zooming Y (parameter) axis.
- Press **Escape** to return to the main menu.

When the **off-line graph** is displayed:

- Use the arrow keys to move along the X (Time) and Y (parameter) axes of the graph.
- Press **SETUP** to access the zoom menu for X and Y axes. Use **Zoom Time** or **Zoom pH** / **Zoom mV** / **Zoom Rel mV** / **Zoom ISE** to switch between the active zooming axes. Press **Zoom IN** or **Zoom OUT** to zoom the selected axis.

Note: While in zoom graph menu the **MODE** key is not accessible.

- Press **Escape** to return to the main menu.

Log History

The measurement, along with **LOG History**, will be visible when this option is selected:

- 1) The last stored logged data (Not actively logging) or
- 2) The last data logged from an active logging lot or
- 3) An empty display - NO LOTS saved, NOT currently logging

The log history list also contains the appropriate mV values, the logged temperature, the temperature probe status source, as well as the records time stamp.

To access the Log History display option:

- Press **Display** while in *Measure* mode. The “Choose Display Configuration” message will be displayed in the Reminder messages area.
- Press **Log History**. The instrument will display the log history regarding the selected *Measure* mode.

- Notes:**
- When an alarm condition is active, all logged records will have an exclamation mark “!”.
 - If logged in Auto Hold, logged records will have an “H”.

12:12:53 PM Dec 15, 2014		Measure	
Channel 2		Stable	
5.782		pH	
Last Cal: Dec 15, 2014 11:36 AM		ATC2 25.7 °C	
Electrode Cond.: 100%			
pH	mV	Temp(°C)	Time
Display	Start Log	Channel	

12:13:57 PM Dec 15, 2014		Measure	
Channel 2		Logging 1 s	
5.622		pH	
Last Cal: Dec 15, 2014 11:36 AM		ATC2 25.7 °C	
Electrode Cond.: 100%			
pH	mV	Temp(°C)	Time
5.622	81.5	25.7 A	12:13:57PM
5.622	81.5	25.7 A	12:13:56PM
5.622	81.5	25.7 A	12:13:55PM
5.622	81.5	25.7 A	12:13:54PM
5.622	81.5	25.7 A	12:13:53PM
5.622	81.5	25.7 A	12:13:52PM
5.622	81.5	25.7 A	12:13:51PM
5.629	81.1	25.7 A	12:13:50PM
5.632	81.0	25.7 A	12:13:49PM
Display	Stop Log	Channel	

- If another *Measure* mode is selected, the Log History will reset.
- If the temperature unit is changed, all logged temperature values will be automatically displayed in the new temperature unit.

SYSTEM SETUP

The System Setup menu allows the user to customize the user interface, view meter information, set the external serial communication interface and to restore the manufacturer settings.

Accessing System Setup

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**. The system setup options will be displayed on the LCD.

To access a System Setup option:


- Use **Δ** or **▽** to highlight the desired option.
- Press **Select** to access the selected option.



The following is a detailed description of the System Setup option screen:

Beeper

This option allows the user to turn an acoustic warning signal on or off. This function can be used to signal 4 different events: a stable signal, an alarm state, when every key is pressed or when an incorrect key is pressed. Enable (or disable) the Beeper for these events. Disabling the Beeper will stop audible signals.

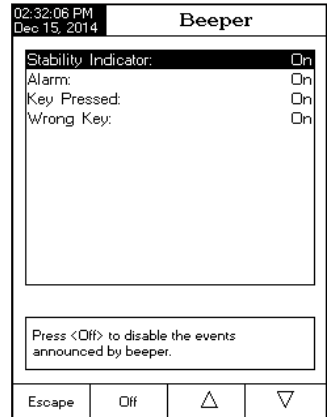

Saving Confirmation

Enable this option to force confirmation of a change made to a setting in GLP data option field or a Sample ID name. If Saving Confirmation is enabled, the user will have to accept the change with a key stroke. If Saving Confirmation is disabled, the changes made to these fields change automatically without asking for confirmation.

02:32:49 PM Dec 15, 2014		System Setup	
Beeper			
Saving Confirmation:		Enabled	
GLP Data			
Date & Time			
LCD Setup			
Color Palette:		Color 3	
Language:		English	
Serial Communication:		38400 bps	
Meter Information			
Restore Factory Settings			
Software Update			
Press <Disable> to disable the saving confirmation option.			
Escape	Disable	△	▽

GLP Data

Use this option to customize logging GLP information with specific identification data. When enabled, these ID tags will be included in the GLP section of all data logs for all modes of operation and both channels (**HI 5222**). Each data field can use up to 10 characters.

The five available fields are:

Operator ID : used to add the name of the operator

Instrument ID : used to name an instrument with a discrete name, location or number

Company Name : used to include the Company ID to the GLP data field.

Additional Info : Two data fields are available for general notes or notations.

To add the GLP Data:

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**
- Use **Δ** or **∇** to select the GLP Data option.
- Press **Select** and use **Δ** or **∇** to highlight the desired option.
- Press **Select** to edit the desired information. The Text Editor menu will be displayed on the LCD.
- Enter the desired information by accepting the highlighted character which is added to the text bar, using **Select**. The **▶** and **∇** keys help the user to select the desired character. It is also possible to delete the last character by positioning the cursor on the Backspace character (**←**) and pressing **Select**.
- Press **Escape** to return to the GLP Data options. If the Saving Confirmation is enabled, press **Yes** to accept the modified option, **No** to escape without saving or **Cancel** to return to the editing mode. Otherwise, the modified options are saved automatically.



Date & Time

Set the current date & time and the format in which they appear.

Set Date and Time

This option allows you to set the current date (year/month/day) and time (hour/minute/second).

- Notes:**
- Only years starting with 2000 are accepted.
 - The time is set using the selected time format. For 12 Hour time format only, the AM/PM can also be selected with **Δ** or **∇**.

Set Time Format












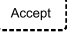
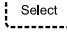



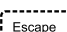
Choose between 12-Hour (AM/PM) time format or 24-Hour time format.


Set Date Format

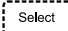Choose the desired date format from 7 available options: DD/MM/YYYY, MM/DD/YYYY, YYYY/MM/DD, YYYY-MM-DD, Mon DD, YYYY, DD-Mon-YYYY or YYYY-Mon-DD.

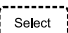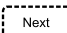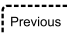To set the Date & Time:

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**

- Use  or  to select the Date & Time option.
- Press  and use  or  to highlight the desired option you want to modify.
- Press  to confirm your selection. Use  /  to select next/previous entry to be edited. Press  and use  or  to set the desired value, then press  to save the modified value (for Set Date and Time option). For the Set Date Format option press  to confirm your selection and select one of the displayed options with  or .
- Press  to confirm your selection and return to the Date & Time options or press  to cancel operation.

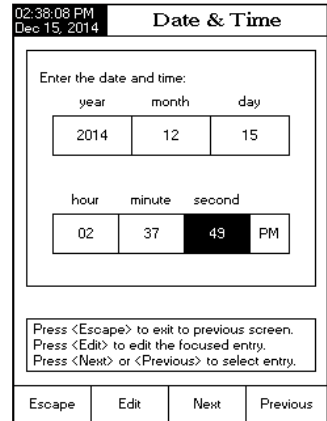


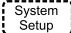







Note: If the time is changed with more than one hour before last pH/ISE calibration, a pop-up warning will appear on the LCD, notifying the user that a date/time conflict has occurred and some time-dependent modes could work improperly (e.g. *Measure*, GLP, Log).


LCD Setup

This option allows the user to set the *Contrast*, the *Backlight* of the LCD and the *Backlight Saver*. The Contrast parameter can be adjusted within 7 steps, while the Backlight parameter within 8 steps. The Backlight Saver can be set from 1 to 60 minutes or it can be OFF (disabled). All the changes are visible on the LCD for each parameter.

Note: If the instrument backlight turns off after the set period of time, press any key to turn it back on.

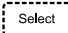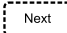To set the LCD Setup:

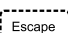- Press  while in *Measure* mode.
- Press .
- Use  or  to select the LCD Setup option.
- Press  and use  key to highlight the desired parameter.
- Use  or  to adjust the contrast / backlight or to set the desired backlight saver time.
- Press  to confirm the modified options and return to the System Setup menu.

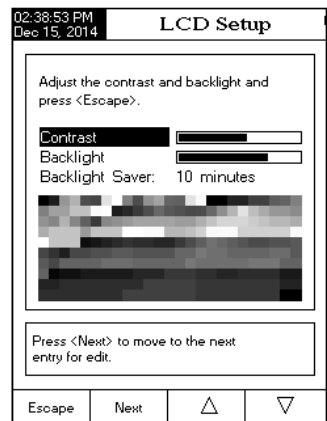


Color Palette

This option allow the user to choose desired color palette.

To select the Color Palette:

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**
- Use **△** or **▽** to select the Color Palette option.
- Press **Select** and use **△** or **▽** to highlight the desired color.
- Press **Select** to confirm your selection and return to the System Setup menu or press **Escape** to return to the System Setup menu without changing.

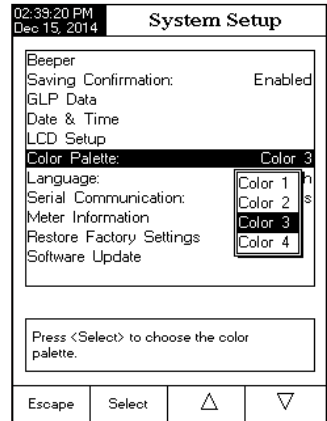

Language

This option allows the user to choose the desired language in which all information will be displayed.

To select the Language:

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**
- Use **△** or **▽** to select the Language option.
- Press **Select** and use **△** or **▽** to highlight the desired language.
- Press **Select** to confirm your selection and return to the System Setup menu or press **Escape** to return to the System Setup menu without changing.




Serial Communication

This option allows the user to set the desired speed for the serial communication (baud rate) in bps. The meter and the PC program must have the same baud rate.

To set the Serial Communication:

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**.
- Use **Δ** or **∇** to select the Serial Communication option.
- Press **Select** and use **Δ** or **∇** to highlight the desired baud rate.
- Press **Select** to confirm your selection and return to the System Setup menu or press **Escape** to return to the System Setup menu without changing.

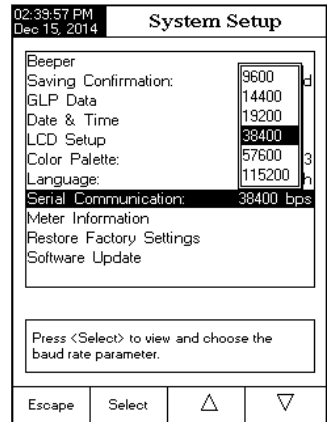


Meter Information

This option provides general information about the instrument serial number (each instrument has a unique identification serial number), the software version and the factory calibration date and time (for mV and temperature).

Note: All instruments are factory calibrated for mV and temperature. One year after factory calibration, a warning message "**Factory Calibration Expired**" will be displayed when powering up the instrument. The instrument will still function, however, it should be taken to the nearest Hanna Customer Service for factory calibration.

To view the Meter Information:

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**.
- Use **Δ** or **∇** to select the Meter Information option.
- Press **Select** to access the Meter Information menu.
- Press **Escape** to return to the System Setup menu.

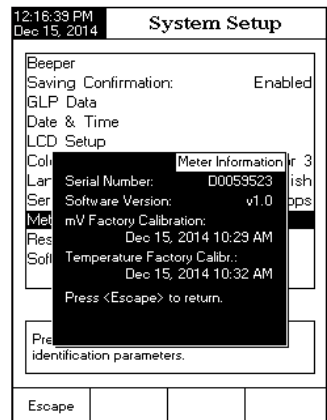


Restore Factory Settings

This option allows the user to erase all user settings and reset the instrument to the default factory settings in pH mode.


To restore the Factory Settings:

- Press **SETUP** while in *Measure* mode.
- Press **System Setup**
- Use **Δ** or **∇** to select the Restore Factory Settings option.
- Press **Select** to confirm your selection. A pop-up menu will be displayed, asking for confirmation.
- Press **Yes** to confirm your selection and return to the System Setup or press **No** to return to the System Setup menu without restoring defaults.
- Press **Escape** to return to *Measure* mode.

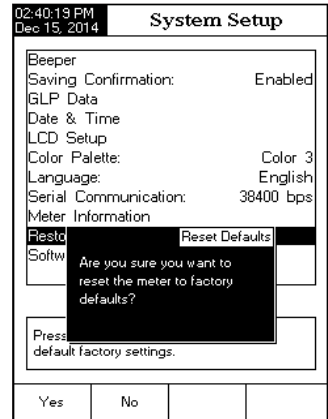

Software update

This function allows the user to update instrument software. In order to start the PC upgrade application, you need to select the proper baud rate, the software update package and start the update.

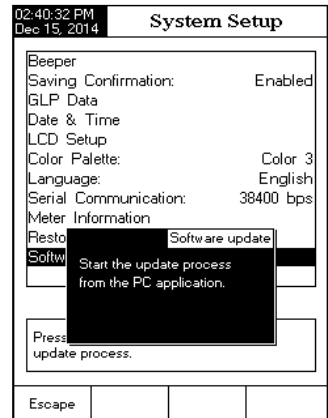

pH SETUP

The pH Setup menu allows the user to set the parameters associated with pH measurement and calibration. These parameters can be set specifically for each channel (HI 5222 only). The settings will be applied only to the active channel.

Accessing pH Setup

- Press **MODE** while in *Measure* mode and then **pH** to select pH range for the desired channel.
- Press **SETUP** and then **pH Setup** to access pH Setup menu.

To access a pH Setup option:

- Use **Δ** or **▽** to highlight the desired option.
- Press **Select** to access the selected option.

The following is a detailed description of the pH Setup option screens.

12:40:11 PM Dec 15, 2014		pH Setup	
Channel 2			
Profile			
Temperature			
Calibration			
Sample ID			
Stability Criteria:		Medium	
Reading Mode:		Direct	
Log			
Alarm			
Isopotential Point:		7.000 pH	
pH Resolution:		X.XXX	
Press <Select> to access the profiles manager.			
Escape	Select	Δ	▽

Profile

This option opens the Profile manager. Enabling Profile allows the user to Save, Load or Delete an application Profile. The Profile option allows the user to store up to ten separate profile applications (HI 5222 - five profiles for each channel). Each Profile can be named and recalled at a moment's notice. A profile is a sensor setup complete with measurement units, logging and display preferences, calibration standards (Buffer or Standards including custom), setup of the Display screen for measurement (i.e. single, dual, graphing, GLP) and any other sensor configuration. Once saved, the exact same profile can be used at another time. This is a handy feature if the meter is used occasionally for additional applications because it saves time in the setup of the meter and ensures that the same procedure will be used.

To save the measurement configuration for pH mode:

- Press **pH Setup** and use **Δ** or **▽** to highlight Profile.
- Press **Enable** / **Disable** to enable / disable this feature.

12:40:58 PM Dec 15, 2014		pH Setup	
Channel 2			
Profile Feature:		Enabled	
Save Profile			
Save Profile As ...			
Load Profile			
Delete Profile			
Press <Disable> to disable the Profile feature.			
Escape	Disable	Δ	▽

The available options are:

Save Profile: save the current profile.








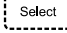
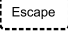
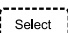
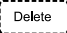
Save Profile As...: save current profile using a specific name.

Load Profile: load from available profiles.

Delete Profile: delete a profile.

Save Current Profile

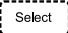To save the current profile:

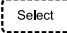- Use  or  to select Save Profile As...
- Press . The *Text Editor* box will be displayed on the LCD.
- Enter the desired profile name by using  and  to highlight the desired character and then press  to add it to the text bar. It is also possible to delete the last character by positioning the cursor on the Backspace character () and pressing .
- Press  to return to the Profile options.
- Use Save Profile to save changes made to a presently used Profile. Changes will overwrite existing configurations.
- Select Load Profile to select a profile to use from the list of saved profiles. Highlight the desired profile and press .
- Select Delete Profile to remove a selected profile from the saved list.
Highlight the profile and press .

Temperature

The temperature has a direct influence on pH. This option allows the user to choose the temperature source and units, as well as the desired manual temperature for manual temperature compensation mode.

Temperature Source (HI 5222 only)

If using a temperature probe, Automatic Temperature Compensation will be performed relative to the displayed temperature, with the “ATC” indicator displayed on the LCD. For the **HI 5222**, one temperature probe can be used for both measurement channels if desired. Select the source by selecting Manual, Channel 1 or Channel 2 . If no temperature probe is detected, Manual Temperature Compensation will be performed, with the “MTC” indicator on the LCD.

Temperature Unit


The desired temperature unit can be chosen (Celsius, Fahrenheit or Kelvin degrees) and the meter will automatically make the conversion for the selected unit.

Manual Temperature

If no temperature probe is connected, the desired temperature can be set manually. The default setting is 25 °C. If the measured temperature is different, the value can be manually adjusted to obtain an accurate pH measurement.

To set one of the Temperature options:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**.
- Use **Δ** or **∇** to select the Temperature option.
- Press **Select** and use **Δ** or **∇** to highlight the desired Temperature option you wish to modify.
- Press **Select** and use **Δ** or **∇** to highlight the desired option (for Temperature Source & Unit options) or use **Δ** or **∇** to adjust the temperature value between the displayed limits (for Manual Temperature option).
- Press **Select** to confirm your selection (for Temperature Source & Unit options) or press **Accept** to save the current value (for Manual Temperature option). Otherwise, press **Escape** to cancel operation.

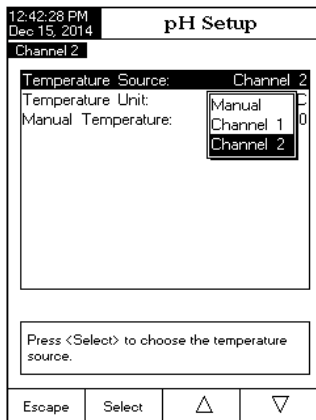

Calibration

This option allows the user to setup desired buffers and calibration reminder.

Buffer Entry Type

Three settings are available for the pH buffers used for electrode calibration:


Automatic - the instrument automatically selects the closest buffer to the measured pH value from the predefined buffers chosen in the option *Edit Buffer Group*.

Semiautomatic - the instrument automatically selects the closest buffers to the measured pH value from all available buffers and you can choose the one used to be, from standard and custom buffers.

Manual Selection - the desired pH buffer is manually selected from all available buffers (standard and custom).

To set the Buffer Entry Type:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**.
- Use **Δ** or **∇** to select the Calibration option.
- Press **Select** and use **Δ** or **∇** to highlight the Buffer Entry Type option.
- Press **Select** and use **Δ** or **∇** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

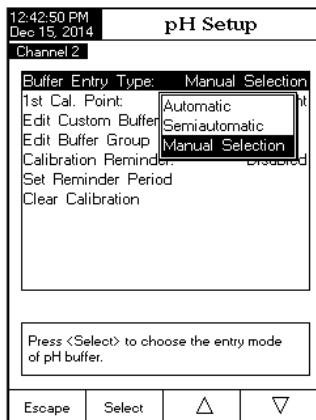

1st Cal. Point

Two options are available for the 1st Cal. Point parameter: Point and Offset.

Point: A new buffer can be added to an existing calibration. The electrode slope will be reevaluated with the addition of this buffer (normal operation).

Offset: The new buffer calibration point can create a constant offset to all existing pH calibration data (existing calibration must have a minimum of two pH buffers).

To set the 1st Cal. Point:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **△** or **▽** to select the Calibration option.
- Press **Select** and use **△** or **▽** to highlight the 1st Cal. Point option.
- Press **Point** / **Offset** as desired.
- Press **Escape** to return to previous menu.

12:43:15 PM Dec 15, 2014	pH Setup		
Channel 2			
Buffer Entry Type:	Manual Selection		
1st Cal. Point:	Point		
Edit Custom Buffers			
Edit Buffer Group			
Calibration Reminder:	Disabled		
Set Reminder Period			
Clear Calibration			
Press <Offset> to set the function of the next first pH calibration buffer.			
Escape	Offset	△	▽

Edit Custom Buffers

If special custom pH buffers are required during calibration, the Edit Custom Buffers option is available. Up to five pH custom buffers can be added. If a custom buffer is used, the user must verify it's value at temperature of calibration.

To edit/set the Custom Buffers:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **△** or **▽** to select the Calibration option.
- Press **Select** and use **△** or **▽** to highlight the Edit Custom Buffers option.
- For a previous set value, press **Invalidate Buffer** to set the custom buffer value to "----" if desired and confirm the setting by pressing **Yes**, otherwise press **Edit Buffer** to edit the selected custom buffer.
- While in edit custom buffer menu press **Reset Buffer** to set the custom buffer value to 7.000 pH and then use **△** or **▽** to set the desired custom buffer value.

12:43:33 PM Dec 15, 2014	Custom Buffers		
Channel 2			
[Cust ----]			
[Cust ----]			
[Cust ----]			
[Cust ----]			
[Cust ----]			
Press <Next Buffer> to choose the next custom buffer for edit.			
Escape	Next Buffer	Edit Buffer	


- Press **Escape** to exit custom buffer edit menu. If the Saving Confirmation is enabled, press **Yes** to accept the modified option, **No** to escape without saving or **Cancel** to return to the editing mode. Otherwise, the modified option is saved automatically.
- Use **Next Buffer** key to select the next custom buffer to be set or press **Escape** to return to Calibration options.

Edit Buffer Group

Accessing this option, the user can edit the desired group of five pH buffers for automatic buffer recognition (Automatic Buffer Entry Type). If the Buffer Group already contains five pH buffers, at least one pH buffer has to be removed in order to add another buffer.

To edit/set the Buffer Group:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Calibration option.
- Press **Select** and use **Δ** or **∇** to highlight the Edit Buffer Group option.
- Press **Select** and use **▶** and **∇** to choose the pH buffer to be included in the buffer group.
- Press **Add** or **Remove** to add/remove the selected pH buffer to/from the buffer group.
- Press **Escape** to return to Calibration options and to save the changes.

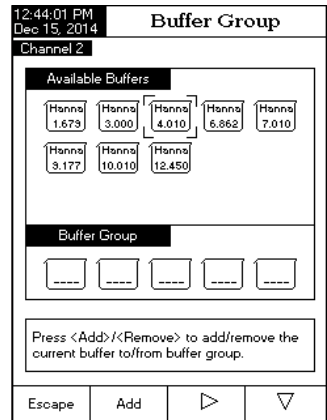


Calibration Reminder

This option allows the user to select a calibration reminder schedule if desired. Three options are available for the calibration reminder: Daily, Periodic or Disabled.

To set the Calibration Reminder:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Calibration option.
- Press **Select** and use **Δ** or **∇** to highlight the Calibration Reminder option.
- Press **Select** and use **Δ** or **∇** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

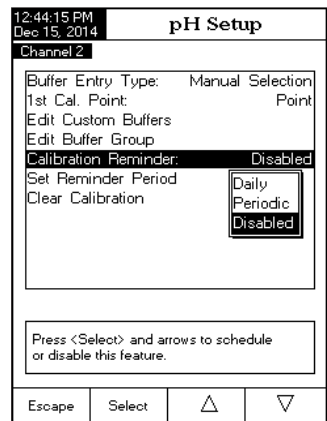

Set Reminder Period

Schedule the calibration reminder timing with this option (verify Daily or Periodic is set for Calibration Reminder).

If a Daily reminder is desired, set the time of day you wish the reminder to occur.

If a Periodic reminder is desired, schedule time in days, hours and/or minutes after the last calibration for the reminder to occur.

To set the Reminder Period:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**.
- Use **Δ** or **∇** to select the Calibration option.
- Press **Select** and use **Δ** or **∇** to highlight the Set Reminder Period option.
- Press **Select** and use **Next** / **Previous** to select next/previous entry to be edited.
- Press **Edit** and use **Δ** or **∇** to set the desired value, then press **Accept** to save the modified value.
- Press **Escape** to return to the Calibration options. If the Saving Confirmation is enabled, press **Yes** to accept the modified option, **No** to escape without saving or **Cancel** to return to the editing mode. Otherwise, the modified option is saved automatically.

12:44:41 PM Dec 15, 2014		Periodic Reminder	
Channel 2			
Enter the time period that must be passed since the last calibration before the time reminder will appear.			
days	hours	minutes	
00	01	00	
Press <Escape> to exit to previous screen. Press <Edit> to edit the focused entry. Press <Next> or <Previous> to select entry.			
Escape	Edit	Next	Previous

Clear Calibration

This feature deletes the pH electrode calibration for the attached electrode and channel (**HI 5222**). A default meter calibration will replace the actual electrode calibration until a new electrode calibration is made.

To clear Calibration:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**.
- Use **Δ** or **∇** to select the Calibration option.
- Press **Select** and use **Δ** or **∇** to highlight the Clear Calibration option.
- Press **Select** to clear calibration. A pop-up menu will be displayed asking for confirmation (when a calibration is available).
- Press **Yes** to confirm or press **No** to escape without saving and return to the Calibration options.

12:44:53 PM Dec 15, 2014		pH Setup	
Channel 2			
Buffer Entry Type:	Manual Selection		
1st Cal. Point:	Point		
Edit Custom Buffers			
Edit Buffer Group			
Calibration Reminder:	Periodic		
Set Reminder Period			
Clear Calibration			
Press <Select> to clear the calibration.			
Escape	Select	Δ	∇

Sample ID

This option allows the user to assign an identification number/name. Two Sample ID options are available: ID Increment and Edit Sample ID.

ID Increment


Two choices are available for the sample ID:

None - the sample ID will be fixed and it can be set alphanumeric (see Edit Sample ID).

Automatic - the sample ID will automatically increment by one for each new log lot.

To set the ID Increment mode:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**.
- Use **△** or **▽** to select the Sample ID option.
- Press **Select** and use **△** or **▽** to highlight the ID Increment option.
- Press **None** / **Automatic** as desired.
- Press **Escape** to return to previous menu.

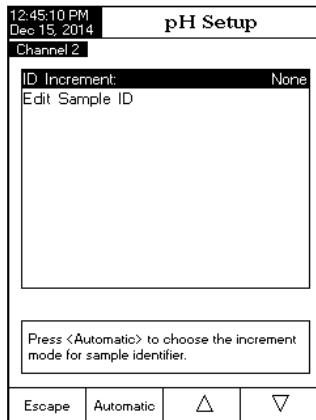

Edit Sample ID

This option allows the user to edit the sample ID.

Note: The ID Increment mode must be set to None, to use this feature.

To edit the Sample ID:


- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**.
- Use **△** or **▽** to select the Sample ID option.
- Press **Select** and use **△** or **▽** to highlight the Edit Sample ID option.
- Press **Select** to confirm your selection.
- If the selected increment is None, the Text Editor menu will be displayed on the LCD, allowing you to enter the desired sample number/name by accepting the highlighted character which is added in the text bar, using **Select**. The **▶** and **▽** keys help the user to select the desired character.



It is also possible to delete the last character; position the cursor on the Backspace character and press

Select

- Press **Escape** to return to Sample ID options. If the Saving Confirmation is enabled, press **Yes** to accept the modified option, **No** to escape without saving or **Cancel** to return to the editing mode. Otherwise, the modified options are saved automatically.
- If the selected increment mode is Automatic, the desired sample ID value can be set using **△** or **▽**.
- Press **Accept** to save the current value or press **Escape** to cancel operation.

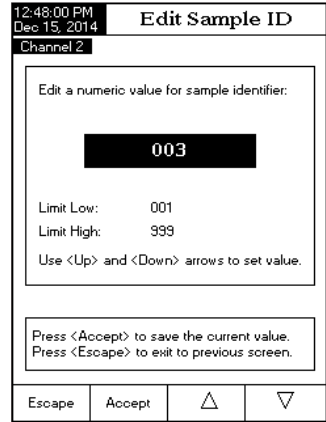

Stability Criteria

This option allows the user to select the signal stability criterion for the measured parameter (pH, mV, ISE):


Fast - this setting will give faster results with less accuracy.

Medium - this setting will give medium speed results with medium accuracy.

Accurate - this setting will give slower results with high accuracy.

To set the Stability Criteria:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **△** or **▽** to select the Stability Criteria option.
- Press **Select** and use **△** or **▽** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

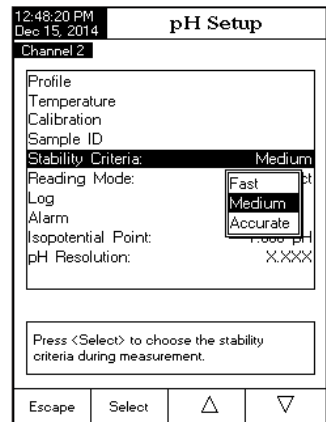


Reading Mode

This option allows the user to select between Direct and Direct/AutoHold pH reading mode. If choosing the AutoHold option, the current reading can be frozen on the LCD when **Auto Hold** is pressed and the stability criterion is reached.

To set the Reading Mode:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Reading Mode option.
- Press **Direct** / **AutoHold** to select Direct / AutoHold option as desired.
- Press **Escape** to cancel operation.

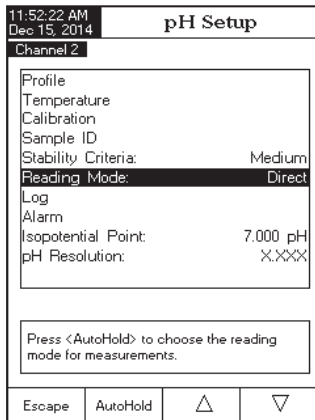

Log

Note: See section **Logging** for available types of logging.

This option allows the user to edit the logging settings: Logging Type, Logging Data Configuration, Sampling Period and New Lot.

Logging Type

Three logging types are available: Automatic, Manual and Auto Hold.


Automatic - the measurement data is logged automatically at predefined time intervals;

Manual - a snapshot of the displayed measurement data is logged with time stamp when the user manually depresses Log;

Auto Hold - this is configured along with the Direct/AutoHold Reading Mode to take a snapshot of stable measurement data. Press **Start Log** to initiate a logging session. Press **Auto Hold** to initiate an Auto Hold event. The log occurs automatically once measurement stability is reached. This type log removes subjective data, as it only captures stable measurements.

To set the Logging Type:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Log option.
- Press **Select** and use **Δ** or **∇** to highlight the Logging Type option.
- Press **Select** and use **Δ** or **∇** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

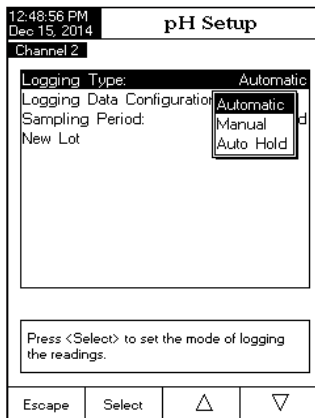


Logging Data Configuration

This option allows the user to select which additional parameters will accompany a log report: Date/Time, Calibration Data, Sample ID, Instrument ID, Operator ID, Company Name, Additional Info 1 and Additional Info 2.

To set the Logging Data Configuration:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **▽** to select the Log option.
- Press **Select** and use **Δ** or **▽** to highlight the Logging Data Configuration option.
- Press **Select** and use **Δ** or **▽** to highlight the desired parameter to be logged in file.
- Press **Yes** to enable the parameter or **No** to disable it.
- Press **Escape** to return to previous menu.

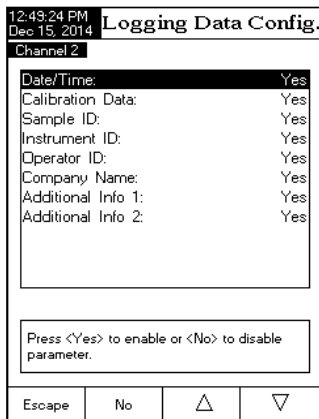


Sampling Period

This option allows the user to select the desired sampling period for automatic logging type.

To set the Sampling Period:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **▽** to select the Log option.
- Press **Select** and use **Δ** or **▽** to highlight the Sampling Period option.
- Press **Select** and use **Δ** or **▽** to select the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

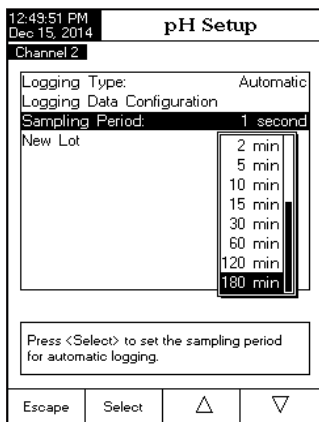


New Lot

This option is used to create a new lot when manual logging is used.

Note: If New Lot option is accessed and the Logging Type is Automatic, a warning message appears on the LCD informing the user that a new lot can be created only if the Logging Type is set as Manual.

To generate a New Lot:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Log option.
- Press **Select** and use **Δ** or **∇** to highlight the New Lot option.
- Press **Select** to generate a new manual lot. A pop-up menu will be displayed asking for confirmation.
- Press **Yes** to confirm or press **No** to escape without saving and return to the Log options.

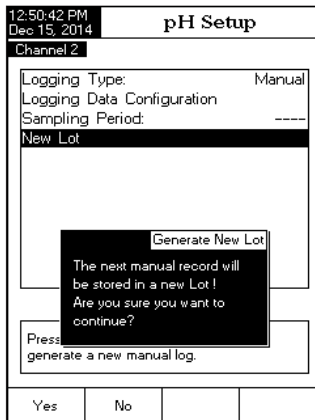

Alarm

This option allows the user to select the alarm settings: Alarm State and Alarm Limits. If the Alarm option is enabled, a continuous double beep will be heard, along with the "Alarm" indicator blinking on the LCD, each time the set limits in Measure mode are exceeded.

Note: Alarm Beeper must be set On for audible beep to be heard.

See: System Setup > Beeper > Alarm.

Alarm State

Three settings are available for the Alarm State option:


Disabled - the alarm will be disabled.

Inside Limits - the alarm state will trigger when the measured value is inside the set limits.

Outside Limits - the alarm state will trigger when the measured value is outside the set limits.

To set the Alarm State:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Alarm option.
- Press **Select** and use **Δ** or **∇** to highlight the Alarm State option.
- Press **Select** and use **Δ** or **∇** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

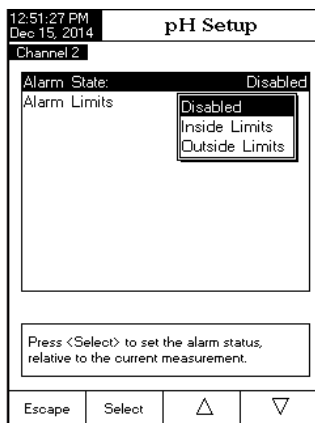


Alarm Limits

This option allows the user to set the alarm limits for the measured value.

Note: The Alarm High value can not be lower than the Alarm Low value.

To set the Alarm Limits:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Alarm option.
- Press **Select** and use **Δ** or **∇** to highlight the Alarm Limits option.
- Press **Select** and use **Next** / **Previous** to select next/previous entry to be edited.
- Press **Edit** and use **Δ** or **∇** to set the desired value, then press **Accept** to save the modified value.
- Press **Escape** to return to the Alarm options. The modified option is saved automatically.

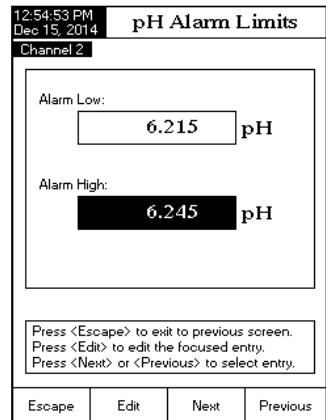

Isopotential Point


This option allows the user to edit the isopotential point of the electrode used for pH measurements. The isopotential point is the mV reading for an electrode at which temperature has no effect on the measurement. The ideal electrode has an isopotential point of 0.0 mV and 7.00 pH, while an actual electrode typically deviates slightly from the ideal values.

If the actual isopotential pH for an electrode is known, it can be set by accessing this option.

Note: If the isopotential point has been modified, recalibration must be performed.

To set the Isopotential Point:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**
- Use **Δ** or **∇** to select the Isopotential Point option.
- Press **Select** and set the desired isopotential pH value using **Δ** or **∇**.
- Press **Accept** to save the current value or press **Escape** to cancel operation.

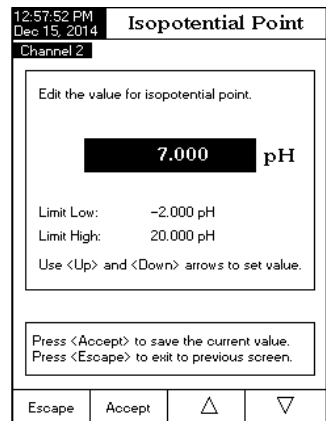


pH Resolution

Select the desired pH resolution with this option. Choose from one (X.X), two (X.XX) or three (X.XXX) digits displayed past the decimals.

To set the pH Resolution:

- Press **SETUP** while in *pH Measure* mode.
- Press **pH Setup**.
- Use **Δ** or **∇** to select the pH Resolution option.
- Press **Select** and use **Δ** or **∇** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

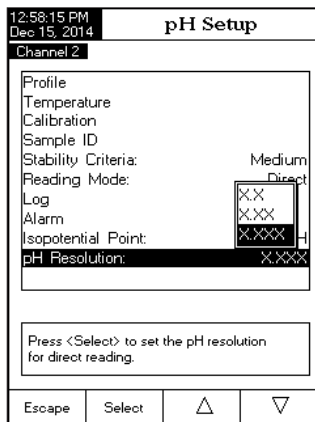

mV SETUP

The mV Setup menu allows the user to set the parameters associated with mV and Relative mV measurements. These parameters can be set specifically for each channel (HI 5222 only). The settings will be applied only to the active channel.

Accessing mV Setup

- Press **MODE** while in *Measure* mode and then **mV** or **Rel mV** to select mV / Rel mV range for the desired channel.
- Press **SETUP** and then **mV Setup** to access mV Setup menu.

To access a mV Setup option:

- Use **△** or **▽** to highlight the desired option.
- Press **Select** to access the selected option.

The following is a detailed description of the mV Setup option screens.

12:58:37 PM Dec 15, 2014	mV Setup		
Channel 2			
Profile			
Temperature			
Sample ID			
Stability Criteria:		Medium	
Reading Mode:		Direct	
Log			
Alarm			
Press <Select> to access the profiles manager.			
Escape	Select	△	▽

Profile

See pH Setup section.

Temperature

ORP measurements are not temperature compensated, although ORP values can change with temperature (e.g. reference electrode potential changes, sample equilibrium changes). It is important to report ORP values together with the reference electrode used and the temperature of measurement.

This option permits selection of the temperature source (HI 5222 only) and measurement units.

12:59:01 PM Dec 15, 2014	mV Setup		
Channel 2			
Temperature Source:		Channel 2	
Temperature Unit:		°C	
Manual Temperature:		25.0	
Press <Select> to choose the temperature source.			
Escape	Select	△	▽

Temperature Source

If using a temperature probe, sample temperature will be displayed with the “ATC” indicator displayed on the LCD. For **HI 5222** only, the ATC option can be selected from Channel 1 or Channel 2. If no temperature probe is detected, a Manually set value will be displayed (and logged) with the measurement.

Temperature Unit

Select the desired temperature unit (Celsius, Fahrenheit or Kelvin degrees) and the Meter will automatically convert to the selected unit.

Manual Temperature

If no temperature probe is connected, the desired temperature can be manually entered. The default setting is 25 °C.

Calibration (Relative mV only)

Calibration Reminder

This option allows the user to select a calibration reminder schedule if desired.

See pH Setup section > Calibration Reminder section for option access details.

Set Reminder Period

See pH Setup section > Set Reminder Period section.

Clear Calibration

This feature deletes the Relative mV calibration for the selected channel.

- Press **SETUP** while in Rel mV mode.
- Press **mV Setup** then use **Δ** or **▽** to access Calibration option.
- Press **Select** and use **Δ** or **▽** to highlight Clear Calibration option.
- Press **Select** to clear calibration. A pop-up menu will be displayed asking for confirmation (when a calibration is available).
- Press **Yes** to confirm or press **No** to escape without saving and return to the Calibration options.

Sample ID - See pH Setup section.

Stability Criteria - See pH Setup section.

Reading Mode - See pH Setup section.

Log - See Logging section, or pH Setup section.

Alarm - See pH Setup section.

ISE SETUP (HI 5222 only)

The ISE Setup menu allows the user to set the parameters regarding ISE measurement and calibration. These parameters can be set specifically for each channel. The settings will be applied only to the active channel.

Accessing ISE Setup

- Press **MODE** while in *Measure* mode and then **ISE** to select ISE range for the desired channel.
- Press **SETUP** and then **ISE Setup** to access ISE Setup menu.

To access an ISE Setup option:

- Use **Δ** or **▽** to highlight the desired option.
- Press **Select** to access the selected option.

The following is a detailed description of the ISE Setup option screens.

12:59:30 PM Dec 15, 2014		ISE Setup	
Channel 2			
Profile			
Reading Mode:	Direct		
Temperature			
Calibration			
Electrode Type:	Fluoride		
Concentration Unit:	ppm		
Sample ID			
Stability Criteria:	Medium		
Log			
Alarm			
ISE Significant Digits:	XXX		
Press <Select> to access the profiles manager.			
Escape	Select	Δ	▽

Profile - See pH Setup section.

Reading Mode

This option allows the user to select the desired reading mode: Direct, Direct/AutoHold, Known Addition, Known Subtraction, Analyte Addition and Analyte Subtraction. Four of these Reading Modes are collectively known as Incremental Methods (see ISE Theory section for details). Direct measurements and Direct/AutoHold measurements are also available.

Direct

Direct measurements are analogous to taking pH measurements. The ISE is calibrated in Ion standards and sample measurements are made directly. The ISE's manual should be consulted for tips and practices of making Direct measurements. The Ion concentration can be read directly from the instrument.

Direct/AutoHold

Direct/AutoHold measurements are made similar to Direct measurements. The advantage of using AutoHold is a measurement that has not reached equilibrium will not be used. Only after the chosen stability criteria has been met will the meter go into the AutoHold mode. Using AutoHold removes the subjective nature of stability.

Known Addition

In the Known Addition method, a sample is measured with an ISE before and after the addition of a known volume of a standard. The mV difference is then used to calculate the concentration of the Ion in the original sample.

Known Subtraction

In the Known Subtraction method, a sample is measured with an ISE before and after the addition of a known volume of a reactant standard. The reactant standard reacts with the measured Ion in the sample, reducing its concentration. The mV difference is then used to calculate the concentration of the Ion in the original sample. The stoichiometric ratio between Reactant Standard and Ion in the sample must be known.

Analyte Addition

Analyte Addition is similar to the Known Addition method, with the difference being that an aliquot of sample is added to a known volume of standard. Both solutions contain the same measured Ion. The standard is measured with an ISE before and after the addition of a known volume of a sample. The Ion concentration is then calculated using the difference in mV potential. The sample should increase the concentration of the Ion being measured.

Analyte Subtraction

In the Analyte Subtraction method, an aliquot of sample is added to a reactant standard of known concentration and volume. The sample partially reacts with the measured Ion. The stoichiometric ratio between standard and sample must be known. The Ion concentration is then calculated using the difference in mV potential.

To set the Reading Mode:

- Press **SETUP** while in *ISE Measure* mode.
- Press **ISE Setup**.
- Use **Δ** or **∇** to select the Reading Mode option.
- Press **Select** and use **Δ** or **∇** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

01:00:45 PM Dec 15, 2014		ISE Setup	
Channel 2			
Profile:	JANE		
Reading Mode:	Direct		
Temperature	Direct		
Calibration	Direct/Auto-Hold		
Electrode Type:	Known Addition		
Concentration Ur	Known Subtraction		
Sample ID	Analyte Addition		
Stability Criteria:	Analyte Subtraction		
Log			
Alarm			
ISE Significant Digits:	XXX		
Press <Select> to choose the reading mode for ISE concentration measurements.			
Escape	Select	Δ	∇

Temperature

This option permits the user to configure all parameters related to ISE temperature measurements.

Temperature Source

The options are: Manual, Channel 1 or Channel 2. If no temperature probe is detected, a manually set value will be displayed (and logged) with the measurement. If a temperature probe is connected to either channel, it may be selected. The temperature measurement will be displayed and logged with the measurement and may be used for temperature compensation calculation if Temperature Compensation is enabled.

Temperature Unit

Select the desired temperature unit (Celsius, Fahrenheit or Kelvin degrees) and the meter will automatically convert to the selected unit.

Manual Temperature

If no temperature probe is connected, the desired temperature can be set manually. The default setting is 25 °C. If the measured temperature is different, the value can be manually adjusted to obtain an accurate Ion measurement.

Temperature Compensation





ISE measurements benefit from temperature compensated corrections if:

- standards and sample temperatures differ from each other
- the Isopotential Point of the ISE is known.

If sample and standards are made at the same temperature, leave this option disabled.



Isopotential Point

If the Temperature Compensation is enabled, the isopotential point of the ISE must be added in this parameter. Verify the *Electrode Type* and *Concentration Unit* are configured for the desired application. The Isopotential point will use the selected concentration unit. Use

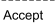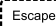 and  to edit the isopotential point value and press  to save the value or press  to cancel operation.

Notes:

- A warning message will appear on the LCD informing the user to perform a new calibration.
- A minimum of two Ion standards is required for the ISE calibration.

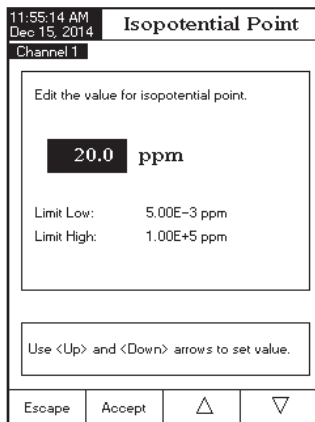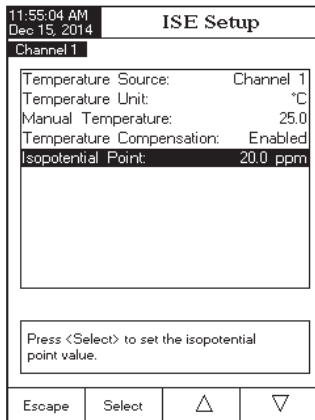

Calibration

This option allows the user to view and configure all ISE parameters related to ISE calibration.

Manual Entry


Two different standard groups can be used for calibration of ISE:

All Standards - During calibration the user can select the desired standards from a large list containing all the predefined standard values and the custom standards.

Group Standards - the user can pre-select a group of standards from the existent group of standards to be used during sensor calibration.

To set the Manual Entry:

- Press **SETUP** while in *ISE Measure* mode.
- Press **ISE Setup**.
- Use **Δ** or **▽** to select the Calibration option.
- Press **Select** and use **Δ** or **▽** to highlight the Manual Entry option.
- Press **All** or **Group** to select the desired option.

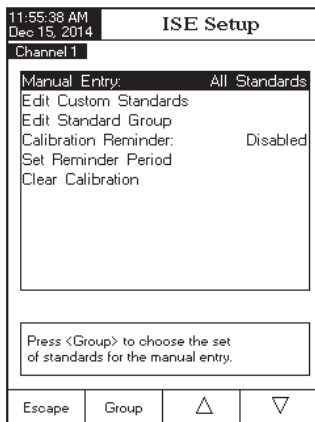


Edit Custom Standards

Use *Edit Custom Standards* function to add additional ISE standard values. Up to five custom standard values can be added. Set *Electrode Type* and *Concentration Unit* prior to adding these standards.

To edit/set the Custom Standards:

- Press **SETUP** while in *ISE Measure* mode.
- Press **ISE Setup**.
- Use **Δ** or **▽** to select the Calibration option.
- Press **Select** and use **Δ** or **▽** to highlight the Edit Custom Standards option.
- If you want to disable the custom standard, press **Invalidate Standard**. A pop-up menu will be displayed asking for confirmation. Press **Yes** to confirm (the custom standard value will turn to "----") or press **No** to cancel the operation.
- Use **Next Standard** key to select the next custom standard to be set.
- Press **Escape** to return to Edit Custom Standard options.

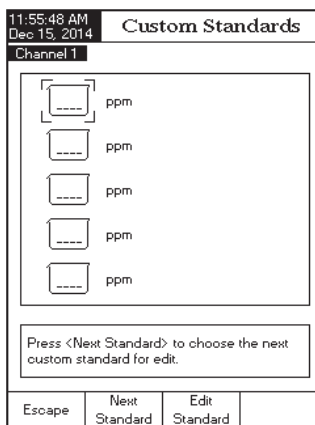


Edit Standard Group

If a *Group Standard* was selected in the parameter Manual Entry, this parameter is used to create your group of standards. If the Standard Group already contains five ISE standards, at least one ISE standard has to be removed in order to add another standard.

To edit/set the Standard Group:

- Press **SETUP** while in *ISE Measure* mode.
- Press **ISE Setup**
- Use **Δ** or **∇** to select the Calibration option.
- Press **Select** and use **Δ** or **∇** to highlight the Edit Standard Group option.
- Press **Select** and use **▶** and **∇** to choose the ISE standard to be included in the standard group.
- Press **Add** / **Remove** to add/remove the selected ISE standard to/from the standard group.
- Press **Escape** to return to Calibration options and to save the changes.

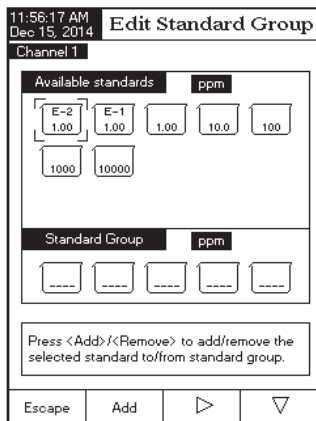

Calibration Reminder - See Calibration option from pH Setup section.

Set Reminder Period - See Calibration option from pH Setup section.

Clear Calibration - See Calibration option from pH Setup section.

Electrode Type

This option allows the user to select the desired Ion Selective Electrode used for measurements from a list: Ammonia, Bromide, Cadmium, Calcium, Carbon Dioxide, Chloride, Cupric, Cyanide, Fluoride, Iodide, Lead, Nitrate, Potassium, Silver, Sodium, Sulfate, Sulfide and five custom ISE. For the standard ISE it is possible to view the Ion constants (Name, Molar Weight and Electric Charge/Slope), while for the custom ISE all these constants can be manually set.

To set the Electrode Type:




- Press **SETUP** while in *ISE Measure* mode.
- Press **ISE Setup**
- Use **Δ** or **∇** to select the Electrode Type option.
- Press **Select** and use **Δ** or **∇** to select the desired standard ISE or a custom one from the list.

For standard ISE:

- Press **View** to visualize the Ion constants and then press **Escape** at any time to exit Ion Constants view mode.
- Press **Select** to confirm your selection and return to *ISE Setup* options.


For custom ISE:

- Press **View** to edit the Ion constants for the selected custom ISE. Use **Δ** or **∇** to select the desired constant and press **Select** to enter edit mode or **Escape** to cancel operation.
- For the Ion name the Text Editor menu will be displayed on the LCD. Enter the desired information by accepting the highlighted character which is added in the text bar, using **Select**. The **▶** and **∇** keys help the user to select the desired character. It is also possible to delete the last character by positioning the cursor on the Backspace character and pressing **Select**. Press **Escape** to return to the Ion Constants menu. If the Saving Confirmation is enabled, press **Yes** to accept the modified option, **No** to escape without saving or **Cancel** to return to the editing mode. Otherwise, the modified option is saved automatically.
- To set the appropriate Ion molar weight (in g/mol units) use **Δ** or **∇** to set the desired value and then press **Accept** to save the current value or press **Escape** to cancel operation.

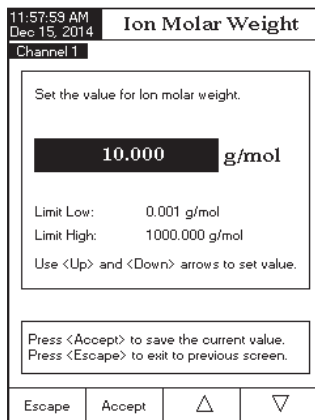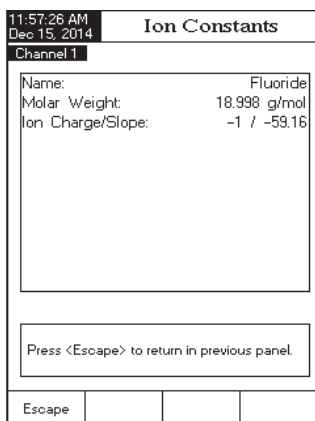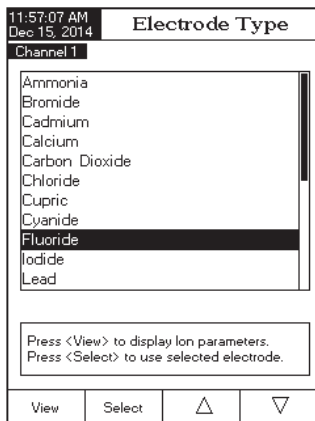

- To select the appropriate Ion Charge/Slope use Δ or ∇ and then press **Select**. If the Ion electric charge is None, its slope can be manually set by pressing **Edit**. A pop-up menu will be displayed on the LCD, in which the slope value can be set using Δ or ∇ . Press **Accept** to save the modified value or press **Escape** to return to the previous menu.

Note: If an ISE calibration was performed and a different Ion Selective Electrode is selected (standard or custom), a warning message appears on the LCD informing the user to perform a new calibration or to select the previous ISE in order to perform accurate measurements.

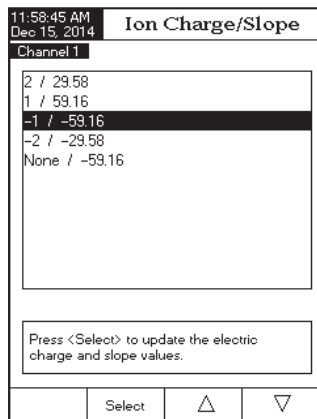


Concentration Unit

Select the desired concentration unit for the measured Ion or chemical compound. The available concentration units are: ppt, g/L, ppm, mg/L, $\mu\text{g/mL}$, ppb, $\mu\text{g/L}$, mg/mL, M, mol/L, mmol/L, %w/v and User (custom unit).

To set the Concentration Unit:

- Press **SETUP** while in *ISE Measure* mode.
- Press **ISE Setup**.
- Use Δ or ∇ to select the Concentration Unit option.
- Press **Select** and use Δ or ∇ to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

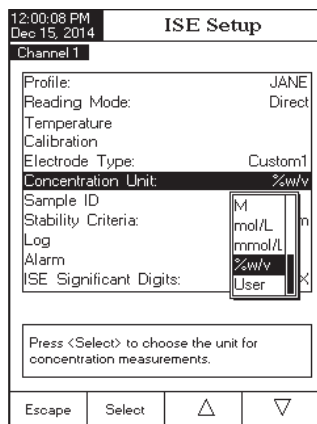

Sample ID - See pH Setup section.

Stability Criteria - See pH Setup section.

Log - See pH Setup section.

Note: The Logging Data Configuration option includes also the Ion Constants parameter. If you want it to appear in the log reports, it must be enabled.

Alarm - See pH Setup section.


Note: The Alarm Limits (Low and High) are set in the selected concentration unit of the measured Ion.

ISE Significant Digits

Accessing this option, the number of ISE significant digits can be set, with one (X), two (XX) or three (XXX) significant digits.

To set the ISE Significant Digits:

- Press **SETUP** while in *ISE Measure* mode.
- Press **ISE Setup**.
- Use **△** or **▽** to select the ISE Significant Digits option.
- Press **Select** and use **△** or **▽** to highlight the desired option.
- Press **Select** to confirm your selection or press **Escape** to cancel operation.

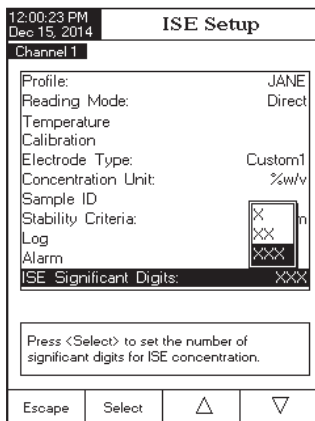

pH CALIBRATION

Calibrate the instrument often, especially if high accuracy is required.

The instrument should be recalibrated:

- Whenever the pH electrode is replaced.
- At least once a week.
- After testing aggressive chemicals.
- When “**Electrode Cond. Unknown**”, “**pH x Default Calibration**” or “**pH x Calibration Expired**” message appears on the LCD, in the Reminder messages area. The “**x**” represents channel “**1**” or channel “**2**”.

PREPARATION

Pour small quantities of the buffer solutions into clean beakers. If possible, use plastic beakers to minimize any EMC interferences.

For accurate calibration and to minimize cross-contamination, use two beakers for each buffer solution. One for rinsing the electrode and one for calibration.

If you are measuring in the acidic range, use pH 7.01 or 6.86 as first buffer and pH 4.01, 3.00 or 1.68 as second buffer. If you are measuring in the alkaline range, use pH 7.01 or 6.86 as first buffer and pH 10.01/9.18 or 12.45 as second buffer.

For extended range measurements (acidic and alkaline), perform a five points calibration by selecting five of the available buffers.

CALIBRATION PROCEDURE

There are 8 standard pH buffers that are temperature-compensated during pH calibrations: 1.68, 3.00, 4.01, 6.86, 7.01, 9.18, 10.01 and 12.45. The standard pH buffers are temperature compensated during calibration. Custom buffers require the user to use the actual buffer value at the temperature of use.

A minimum of a two point calibration using the standard pH buffers is required to determine the pH electrode condition. The buffers should bracket the sample measurement pH.

An extended pH measurement range will require calibration at multiple points. The meter is able of calibration with 5 pH buffers. For improved measurement accuracy, perform a multiple buffer calibration bracketing and including the pH range the sample measurements.

The buffer group that will be available during calibration was set in pH setup → Calibration Buffer Entry Type. The following example demonstrates pH electrode calibration if Manual selection was selected. In this case all of the 8 standard buffers will be available for calibration.

pH Calibration screen description


The screenshot shows the pH Calibration screen with the following details:

- Time: 03:02:58 PM, Date: Dec 15, 2014
- Channel 1
- Stable
- Measured buffer value: 9.965 pH
- Calibration buffer: -173.0 mV (Hanna 10.010)
- Temperature: 22.1°C (ATC1)
- Calibrated Buffers: Hanna 4.010, Hanna 7.010
- Last Calibration: Dec 15, 2014 03:02 PM
- Calibration messages: Press <Accept> to update calibration.
- Navigation buttons: Escape, Accept, Next Buffer, Previous Buffer

- Press **CAL**. If the instrument was calibrated before and calibration was not cleared, the old calibration can be cleared by pressing **Clear Cal**. After 10 seconds, **Clear Cal** will no longer be available.

Note: It is very important to clear calibration history when a new electrode is used because most errors and warning messages that appear during calibration depend on calibration history.

- Immerse the pH electrode and the temperature probe approximately 4 cm (1.5") into a buffer solution of your choice (pH 1.68, 3.00, 4.01, 6.86, 7.01, 9.18, 10.01, 12.45 or a custom buffer) and stir gently. The temperature probe should be close to the pH electrode.
- Select the pH calibration buffer used with **Next Buffer** or **Previous Buffer**. The "Please wait..." message will appear on the LCD until the reading is stable or the buffer is validated.
- If the pH buffer is validated, **Accept** will appear on the LCD. Press **Accept** to update calibration. The calibration buffer will be added to the Calibrated Buffers section.
- Immerse the pH electrode and the temperature probe into the next buffer solution and follow the above procedure or press **Escape** to exit calibration.

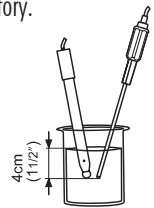


- Notes:**
- The new added calibration point will replace an old one if the difference between them is ± 0.2 pH.
 - If the existing stored calibration is full (five calibration points), a pop-up menu will be displayed on the LCD in which you can select with **<Left>** or **<Right>** keys the buffer you want to replace with the current buffer. Press **Remove** to delete the selected buffer and then press **Accept** to update calibration with the new buffer.

The screenshot shows the pH Calibration screen with the following details:

- Time: 10:50:59 AM, Date: Dec 15, 2014
- Channel 1
- Stable
- Measured buffer value: 9.175 pH
- Calibration buffer: -126.9 mV (Hanna 3.177)
- Temperature: 24.5°C (ATC1)
- Remove Buffer menu: Hanna 3.175, Hanna 3.177, Hanna 3.178, Hanna 3.179, Hanna 3.180
- Instructions: Use <Left> and <Right> arrows to select which old buffer you want to be removed. Press <Accept> to update calibration.
- Navigation buttons: Escape, Remove, <Left>, <Right>

- If using manual temperature, after selecting the standard buffer, press **SETUP**. A pop-up menu will be displayed on the LCD in which the temperature value can be adjusted using Δ or ∇ . Press **Accept** to save the new temperature value.
- If using **Custom buffers**, press **SETUP** after buffer has been accepted to change actual buffer conditions. A pop-up menu will be displayed on the LCD in which the custom buffer and the temperature value (MTC) can be adjusted by pressing **Edit** and then Δ or ∇ keys. Press **Accept** to save the modified value and then **Next** / **Previous** to select next/previous value to be adjusted.
- If the **Automatic buffer** entry type has been selected for the calibration procedure, the instrument will automatically select the closest buffer to the measured pH value from the edit buffer group (see pH Setup for details).
- If the **Semiautomatic buffer** entry type has been selected for the calibration procedure, the instrument will display only the closest buffers to the measured pH value from all the available buffers and the user must select with **Next Buffer** or **Previous Buffer** the buffer being used.

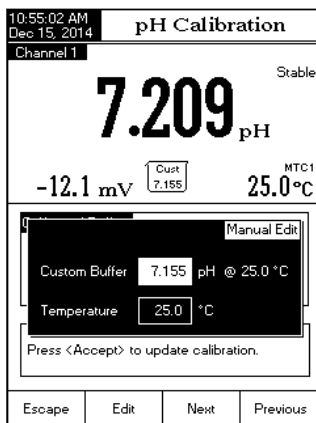

CALIBRATION MESSAGES

- **Move sensor to next buffer or check buffer:** this message appears when the difference between the pH reading and the value of the selected calibration buffer is significant. If this message is displayed, check if you have selected the appropriate calibration buffer.
- **Wrong buffer temperature:** this message appears if the buffer temperature is out of the defined buffer temperature range.
- **Clean the electrode or check the buffer. Press **Accept** to update calibration:** this message alerts the user that some dirt or deposits could be on the electrode. Refer to the electrode Cleaning Procedure.
- **Slope too low. Please check the buffer / Slope too high. Please check the buffer:** these messages appear if the current slope is under 80 % or over 110 % of default slope. Recalibrate the instrument using fresh buffers.
- **Slope too low. Press **Clear Cal** to clear old calibration / Slope too high. Press **Clear Cal** to clear old calibration:** verify the correct buffer has been selected and poured.

- **Unrecognized buffer.** Please check the buffer or the buffer list (for Semiautomatic and Automatic buffer entry type): this message appears if the current buffer value is not close to any of the buffers from the buffer list/group. Check if the current buffer is present in the buffer list or the appropriate buffer group was selected.
- **The current buffer was already calibrated:** change the buffer or press Escape to exit calibration mode.

pH MEASUREMENT


Verify the pH electrode and instrument has been calibrated before making pH measurements.

DIRECT MEASUREMENT

To measure the pH of a sample using the Direct reading mode:

- Press MODE and then pH to select *pH Measure* mode (select channel first for HI 5222).
- Select the Direct reading mode (see pH Setup for details).
- Place the electrode tip and the temperature probe approximately 4 cm (1.5") into the sample to be tested. Allow time for the electrode to stabilize. "Stable" will appear.
- The measured pH value will be displayed on the LCD, together with a short GLP information and display preferences.


Note: If the reading is out of range, "----" will be displayed on the LCD.

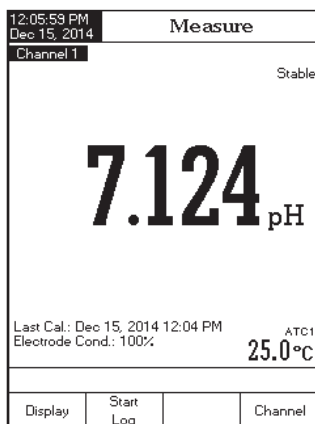

DIRECT/AUTOHOLD MEASUREMENT

To measure pH of a sample using the Direct/AutoHold reading mode:

- Press MODE and then pH to select *pH Measure* mode (for the selected channel - HI 5222 only).
- Select the Direct/AutoHold reading mode (see pH Setup for details).
- Place the electrode tip and the temperature probe approximately 4 cm (1.5") into the sample to be tested.
- The measured pH value will be displayed on the LCD. Press Auto Hold and the "AutoHold" indicator will start blinking on the LCD until the stability criterion is reached. The pH value will be frozen on the LCD, along with "AutoHold" indicator.

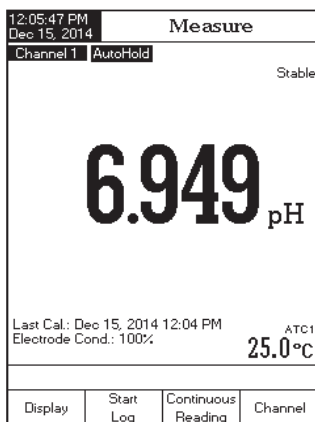

- To return to normal *Measure* mode press 





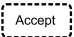
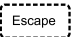
Note: If the reading is out of range, “----” will be displayed on the LCD.


Outside Cal Range warns the user if the current reading is out of the calibrated area. The calibrated area is that part of the pH range in which the calibration point ensures an accurate reading. If the reading is taken out of the calibration area, the “**Outside Cal Range**” message will start blinking on the LCD. The calibrated area is calculated in accordance with the pH resolution used during the measurement. To avoid triggering this message, the buffer values have to be well-distributed in the desired measurement range.

If measurements are taken successively in different samples, it is recommended to rinse the electrode thoroughly with deionized water or tap water and then with some of the next sample before immersing it into the next sample solution.

The pH reading is affected by temperature. In order to measure the pH accurately, the temperature effect must be compensated. To use the **Automatic Temperature Compensation (ATC)** feature, connect and place the **HI 7662-W** temperature probe into the sample as close as possible to the electrode and wait for a few seconds.

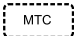If the temperature of the sample is known, **Manual Temperature Compensation (MTC)** can be used by disconnecting the temperature probe.

- Notes:**
- For single channel pH measurements (HI 5221), “MTC” or “ATC” indicators will be displayed on the LCD, while for dual channel pH measurements (HI 5222) “MTC1”/“MTC2” or “ATC1”/“ATC2” indicators will be displayed (Channel dependent).
 - For mV/Rel mV measurements, “NoProbe” or “TEMP” indicators will be displayed on the LCD (HI 5221 only), respectively, “NoProbe 1”/“NoProbe 2” or “TEMP1”/“TEMP2” indicators (HI 5222 only) for mV/Rel mV/ISE measurements, depending on the temperature probe status and channel selected.
 - When in MTC mode, the temperature can be modified by pressing  for *pH Measure* mode and  for *mV/Rel mV Measure* mode (HI 5221 only) if the Reading Mode option is Direct (HI 5222 only). The temperature value can be adjusted with  or  from -20.0 °C to 120.0 °C. Press  to save the new temperature value or press  to return to *Measure* mode.
 - When in ATC mode for pH, or TEMP for mV/Rel mV, “----” will be displayed on the LCD if the measured temperature is under or over temperature range (-20.0 °C to 120.0 °C).

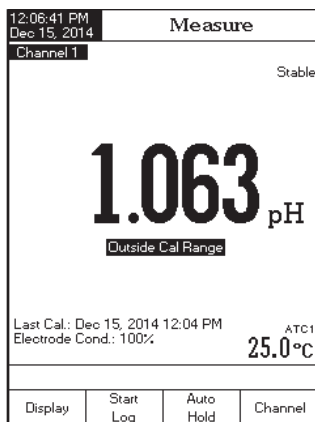

mV & Relative mV MEASUREMENTS

mV/ORP MEASUREMENTS

Oxidation-reduction potential (ORP) measurements provide the quantification of the oxidizing or reducing power of the tested sample.

To correctly perform a redox measurement, the surface of the ORP electrode must be clean and smooth.

DIRECT MEASUREMENT

To measure the mV of a sample using the Direct reading mode:

- Press **MODE** and then **mV** to enter *mV Measure* mode (Select channel first for - HI 5222 only).
- Select the Direct reading mode (see mV Setup for details).
- Place the tip of the ORP electrode 4 cm (1.5") into the sample to be tested and allow a few seconds for the reading to stabilize.
- The instrument will display the measured mV value on the LCD.

Note: If the reading is out of range, "----" will be displayed on the LCD.




DIRECT/AUTOHOLD MEASUREMENT

To measure mV of a sample using the Direct/AutoHold reading mode:

- Press **MODE** and then **mV** to select *mV Measure* mode (Select channel first for - HI 5222 only).
- Select the Direct/AutoHold reading mode (see mV Setup for details).
- Place the tip of the ORP electrode approximately 4 cm (1.5") into the sample to be tested.
- The measured mV value will be displayed on the LCD. Press **Auto Hold** and the "AutoHold" indicator will start blinking on the LCD until the stability criterion is reached. The mV value will be frozen on the LCD, along with "AutoHold" indicator.
- To return to normal *Measure* mode press **Continuous Reading**.

Note: If the reading is out of range, "----" will be displayed on the LCD.

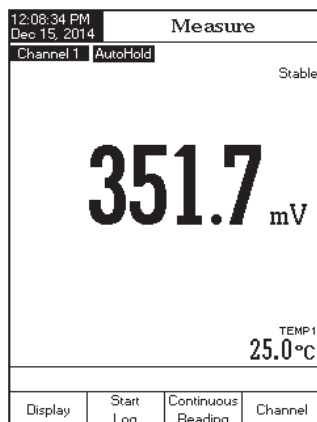


Relative mV MEASUREMENTS

To measure the Relative mV of a sample:

- Press **MODE** then **Rel mV** (select channel first for HI 5222).
- Verify if a current calibration has been made.
- If required, conduct the single point Rel mV calibration. Verify the tip of the electrode is immersed into the known solution or ORP standard.
- Press **CAL**. Use **△** and **▽** arrow keys set the standard value. Press **Accept** to store the calibration.
- Press **MODE** then **Rel mV** (select channel first for HI 5222).
- Place calibrated sensor tip into the sample to be analyzed. The instrument will display the measured Relative mV value on the LCD, together with a short GLP information about the last calibration or Offset: 0.0 mV no Rel mV calibration was performed.

Notes: If the ORP sensor is not in solution or the measured mV potential is out of range, “----” will be displayed on the LCD.

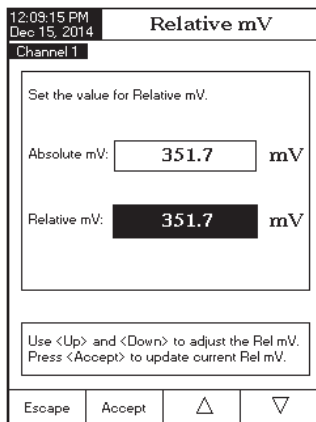

ISE CALIBRATION (HI 5222 only)

For greater accuracy, it is recommended to calibrate the ISE sensors frequently. The instrument should be recalibrated when “ISE x Calibration Expired” (the “x” represents channel “1” or channel “2”) message appears on the LCD, in the Reminder messages area.

Due to electrode conditioning time, the electrode must be kept immersed a few seconds to stabilize. The user will be guided step by step during calibration with easy-to-follow messages on the display. This will make the calibration a simple and error-free procedure.

PREPARATION

Pour small quantities of the standard solutions into clean beakers. If possible, use plastic beakers to minimize any EMC interferences.

For accurate calibration and to minimize cross-contamination, use two beakers for each standard solution. One for rinsing the electrode and one for calibration.

Note: To read concentration (not activity) ISA must be added to the standards and samples. No corrections are needed due to dilutions.

CALIBRATION PROCEDURE

The ISE calibration and measurement can be performed with or without temperature compensation. If the temperature compensation option is enabled, the isopotential point of the electrode must be set in ISE Setup in order to perform correct concentration measurements.

Before calibrating, make sure that the appropriate Electrode Type has been selected in ISE Setup according to the measured Ion/compound.

ISE Calibration screen description

The screenshot shows the ISE Calibration screen with the following elements and annotations:

- Measured concentration:** 10100 mg/L (Stable)
- Calibration Standard:** 371.2 mV (10000) and 22.3°C (ATC1)
- Calibrated Standards:** E-2, 1.00, 10.0, 100, 1000
- Calibration messages:** Press <Accept> to update calibration.
- Navigation buttons:** Escape, Accept, Next Standard, Previous Standard

Additional information from the screen: 03:26:38 PM, Dec 15, 2014, Channel 1, ISE: Sodium, Last Calibration: Dec 15, 2014 03:25 PM


The group of calibration standards is set under ISE Setup → Calibration. Select standards that are in the measurement range of the samples.

To calibrate the instrument:

- Press **CAL**. If the instrument was calibrated before and calibration was not cleared, the old calibration can be cleared by pressing **Clear Cal**. After 10 seconds, **Clear Cal** will no longer be available.

Note: It is very important to clear calibration history when a new electrode is used because most errors and warning messages that appear during calibration depend on calibration history.

- Add ISA to both standard solutions and samples.
- Immerse the Ion Selective Electrode and the temperature probe approximately 2 cm (1") into the less concentrated standard solution and stir gently.
- Select the appropriate standard solution concentration with **Next Standard** or **Previous Standard**. For *All Standards* manual entry mode, the standard concentration can be selected from a list containing all the predefined and custom standards. For *Group Standard* manual entry mode the standard concentration can be selected from the predefined group of standards. Press **Accept** to calibrate the electrode in the standard.

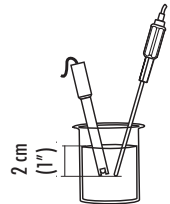


Note: To adjust standard value: Press **SETUP**. A pop-up menu will be displayed on the LCD in which the concentration value can be adjusted using **Δ** or **∇**. Press **Accept** to save the new concentration value.


- The “Please wait...” message will appear on the LCD for 10 seconds. Remove ISE from first standard, rinse tip and immerse the Ion selective electrode and the temperature probe into the next standard solution and follow the above procedure or press **Escape** to exit calibration.

Notes: • The new added calibration point will replace an old one if the difference between them is less than 20 % of the standard solution.

- If the existing stored calibration is full (five calibration points), a pop-up menu will be displayed on the LCD in which you can select with **<** or **>** the standard solution you want to replace with the current one. Press **Remove** to delete the selected calibrated point and then press **Accept** to update calibration with the new standard solution.

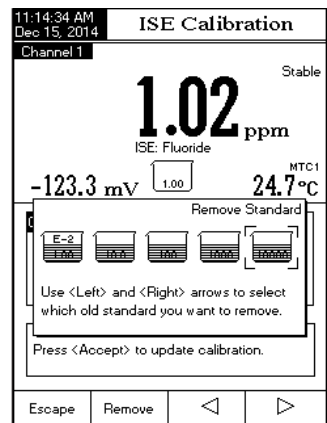

- If the isopotential point of the electrode is unknown, the ISE calibration and measurements can be performed without temperature compensation (see ISE Setup, Temperature option for details).
- When in MTC mode, after selecting a standard press **SETUP**, a pop-up menu will be displayed on the LCD in which the concentration and the temperature value can be adjusted by pressing **Edit** and then **Δ** or **∇** keys. Press **Accept** to save the modified value and then **Next** / **Previous** to select next/previous value to be adjusted. MTC value will have no effect on measurement but will be included on log data.

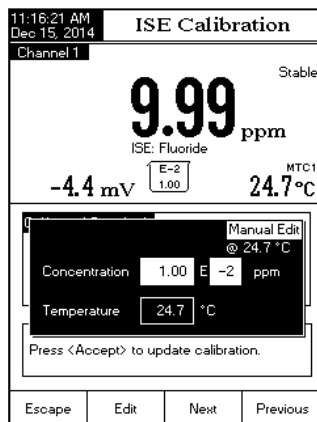

CALIBRATION MESSAGES

- **Wrong standard solution. Check the standard solution:** this message appears when the difference between the reading and the value of the selected standard solution concentration is significant. If this message is displayed, check if you have selected the appropriate calibration standard.
- **Standard too close. Check the standard or clear calibration:** this message appears when the difference between current ISE standard and the already calibrated standard is too low.
- **Slope too low/high. Check the standard solution.** Recalibrate using fresh standards.
- **Difference between standards temperature is too high.** Press **Accept** to update calibration or clear old calibration.

ISE MEASUREMENT (HI 5222 only)

Make sure the instrument and ISE sensor have been calibrated before making ISE measurements.

When using one of the incremental methods for measurement, at least a two point ISE calibration must be performed to establish the electrode slope.


For accurate measurements, add the appropriate ISA (Ionic Strength Adjuster) to both samples and standards. Consult ISE manual for sensor preparation details.

DIRECT MEASUREMENT

To measure the concentration of a sample using the Direct reading mode:

- Press **MODE** and then **ISE** to select *ISE Measure* mode for the selected channel.
- Select the Direct reading mode (see ISE Setup for details).
- Add ISA to the sample solution.
- Submerge the Ion Selective Electrode tip and the temperature probe approximately 2 cm (1") into the sample. Allow time for the electrode to stabilize.
- The measured concentration value will be displayed on the LCD in the selected units.


Note: If the reading is out of range, "----" will be displayed on the LCD.

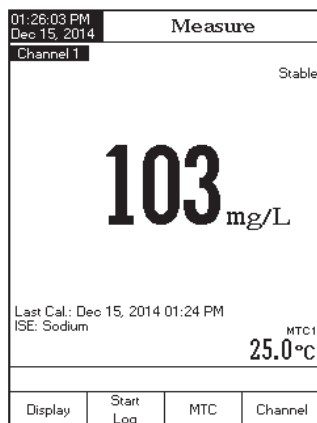

DIRECT/AUTOHOLD MEASUREMENT

To measure the concentration of a sample using the Direct/AutoHold reading mode:

- Press **MODE** and then **ISE** to select *ISE Measure* mode for the selected channel.
- Select the Direct/AutoHold reading mode (see ISE Setup for details).
- Add ISA to the sample solution.
- Dip the Ion Selective Electrode tip and the temperature probe approximately 2 cm (1") into the sample to be tested.
- The measured concentration value will be displayed on the LCD. Press **Auto Hold** the "AutoHold" indicator will blink on the LCD until the stability criterion is reached. The concentration value will be frozen on the LCD, along with "AutoHold" indicator.

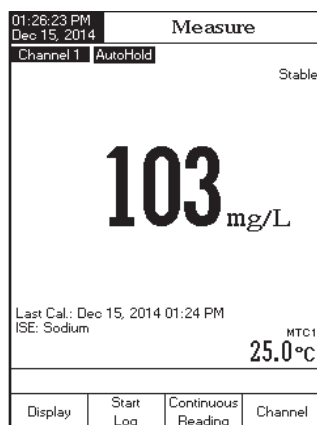

- To return to normal Measure mode press Continuous Reading.

Note: If the reading is out of range, “----” will be displayed on the LCD.

KNOWN ADDITION

To measure the concentration of a sample using the Known Addition incremental method:

- Press MODE and then ISE to select ISE Measure mode for the selected channel.
- Select the Known Addition method (see ISE Setup for details).
- Prior to starting a KA procedure, the ISE sensor must be calibrated with a minimum of two standards containing ISA. The slope of the electrode will be used in all calculations involved in KA.
- If following an established procedure: Press KA. Edit the method variables and follow the procedure below.
- Press Edit to set the method parameters. Press Next / Previous to select next/previous parameter to edit, then press Edit and use Δ or ▽ to set the desired parameter value. Press Accept to save the modified value and then press Escape to exit method parameters edit menu.
- If developing a procedure: Before attempting Known addition analysis it is important to determine what sample volume, standard concentration and standard volume will produce the best results. As a general rule, the addition of standard should change the mV value of the sample by 15 - 20 mV. For a positively charged Ion (i.e. Sodium, Potassium, Calcium), the standard addition should increase the mV. For a negatively charged Ion (i.e. Sulfide, Fluoride, Chloride), the standard addition should decrease the mV. Start with a small trial. For example: Measure 50 mL of sample, add a magnetic stir bar, place on a stirrer, add ISA (consult ISE manual) and place ISE electrode tip into the sample. Put instrument in mV mode and record the observed mV. Using a micropipet, add a volume of the highest ISE standard available (i.e. 0.1M or 1000 ppm). Start by adding 500 μ L at a time (for example) Watch the change in mV. When you have observed approximately a 15 mV change from the original sample. Calculate the total volume added. Adjust sample and standard volumes proportionally to standard volumes that can be measured with accuracy. Use volumetric pipettes for standard, ISA and sample addition.
- Press KA: Edit the procedure variables to the volumes determined in the prior step.

01:27:38 PM Dec 15, 2014		Known Addition	
Channel 1		244.6	Stable TEMP1 25.0 °C
<div style="border: 1px solid black; padding: 5px;"> <p style="text-align: center;">First Step First Reading</p> <p style="text-align: right;">Manual Edit</p> <p>Sample Vol. 100.000 mL</p> <p>ISA Vol. 2.000 mL</p> <p>Std. Vol. 10.000 mL</p> <p>Std. Conc. 100 mg/L</p> <p style="text-align: center;">then press <Continue>.</p> </div>			
Escape	Edit	Next	Previous

01:27:55 PM Dec 15, 2014		Known Addition	
Channel 1		244.5	Stable TEMP1 25.0 °C
<div style="border: 1px solid black; padding: 5px;"> <p style="text-align: center;">First Step First Reading Second Step Second Reading</p> </div>			
Sample Volume:		100.000 mL	
ISA Buffer Vol. :		2.000 mL	
Reagent Volume:		10.000 mL	
Reagent Conc.:		100 mg/L	
<div style="border: 1px solid black; padding: 5px;"> <p>Add 2.000 mL ISA to 100.000 mL of Sample, and immerse the electrode, then press <Continue>.</p> </div>			
Escape	Continue	Edit	

Procedure:

- Press **KA** to enter Known Addition mode.
- Volumetrically add sample to a clean beaker. Add Stir bar and place on a magnetic stir plate. Stir sample. The method will prompt user to add ISA. Place ISE sensor tip into the solution and a mV value will show on the display.
- Press **Continue** to take the first mV reading.
- When the reading is stable, press **Read** to store the first mV reading. The second step of the method will be displayed on the LCD in which the user is notified to add the Volume of Standard to the sample.
- Press **Continue** to take the second mV reading.
- When the reading is stable, press **Read** to store the second mV reading. The ISE measurement results will be displayed on the LCD.
- Press **Save** to log the current results into a ISE Method Report. Press **Direct Measure** to return to ISE Measure mode. Press **Start KA** to measure additional samples. Rinse ISE sample between samples.
- Press **Edit** to modify the method parameters.

Note: Press **Escape** at any time to stop the measurement and return to *ISE Measure* mode.

01:28:39 PM Dec 15, 2014		ISE Results	
Channel 1		102 mg/L	
Sample ID:			
Calculated Slope:		98.2 %	
Reading 1:		244.5 mV	
Reading 2:		244.5 mV	
Sample Volume:		100.000 mL	
Reagent Volume:		10.000 mL	
ISA Volume:		2.000 mL	
Reagent Conc.:		100 mg/L	
Press <Direct Measure> to return in main measurement panel.			
Press <Save> to log the current results.			
Direct Measure	Save	Edit	Start KA

KNOWN SUBTRACTION

To measure the concentration of a sample using the Known Subtraction method:

- Press **MODE** and then **ISE** to select *ISE Measure* mode for the selected channel.
- Select the Known Subtraction method (see ISE Setup → Reading Mode).
- Prior to starting a KS procedure, the ISE sensor must be calibrated with a minimum of two standards containing ISA. The slope of the electrode will be used in all calculations involved in KS.
- If following an established procedure: Press **KS** then edit the method variables and follow the procedure below.
- Press **Edit** to set the method parameters. Press **Next** / **Previous** to select next/previous parameter to edit, then press **Edit** and use **Δ** or **▽** to set the desired parameter value. Press **Accept** to save the modified value and then press **Escape** to exit method parameters edit menu.
- If developing a procedure: Before attempting Known Subtraction analysis it is important to determine what sample volume, standard reactant concentration and standard volume will produce the best results and the way the Reagent will react with the measured Ion on a molar basis (Stoichiometric factor). As a general rule, the addition of standard should change the mV value of the sample by 15-20 mV.

For a positively charged Ion (i.e. Calcium), the reactant addition should decrease the mV. For a negatively charged Ion (i.e. Sulfide, Fluoride, Chloride), the reactant addition should increase the mV. Start with a small trial. For example: Measure 50 mL of sample, add a magnetic stir bar, place on a stirrer, add ISA (consult ISE manual) and place ISE electrode tip in to the sample. Put instrument in mV mode and record the observed mV. Using a micropipet, add a volume of the Reactant standard. Start by adding 500 μL at a time (for example). Watch the change in mV. When you have observed approximately a 15 mV change from the original sample, calculate the total volume added. Adjust sample and standard volumes proportionally to standard volumes that can be measured with accuracy. Use volumetric pipettes for standard, ISA and reagent addition.

- Press **KS** then edit the procedure variables to the volumes determined in the prior step.

Procedure:

- Press **KS** to enter Known Subtraction mode.
- Volumetrically add sample to a clean beaker. Add stir bar and place on a magnetic stir plate. Stir sample. The method will prompt user to add ISA. Place ISE sensor tip into the solution and a mV value will show on the display.
- Press **Continue** to take the first mV reading.
- When the reading is stable, press **Read** to store the first mV reading. The second step of the method will be displayed on the LCD in which the user is notified to add the Volume of Reagent to the sample.
- Press **Continue** to take the second mV reading.
- When the reading is stable, press **Read** to store the second mV reading. The ISE measurement results will be displayed on the LCD.
- Press **Save** to log the current results into an ISE Method Report. Press **Direct Measure** to return to *ISE Measure* mode. Press **Start KS** to start another measurement. Rinse ISE sensor between samples.
- Press **Edit** to modify parameters.

Note: Press **Escape** at any time to stop the measurement and return to *ISE Measure* mode.

ANALYTE ADDITION

To measure the concentration of a sample using Analyte Addition method:

- Press **MODE** and then **ISE** to select *ISE Measure* mode for the selected channel.
- Select the Analyte Addition method (see ISE Setup \rightarrow Reading Mode).
- Prior to starting an AA procedure, the ISE sensor must be calibrated with a minimum of two standards containing ISA. The slope of the electrode will be used in all calculations involved in AA.
- If following an established procedure: Press **AA** then edit the method variables and follow the procedure below.

- Press **Edit** to set the method parameters. Press **Next** / **Previous** to select next/previous parameter to edit, then press **Edit** and use **Δ** or **∇** to set the desired parameter value. Press **Accept** to save the modified value and then press **Escape** to exit method parameters edit menu.
- If developing a procedure: Before attempting Analyte Addition analysis, it is important to determine which standard volume, concentration and sample size will produce the best results. As a general rule, the standard must be less concentrated than the sample so the addition of sample will increase the total Ion content of the beaker and change the mV value by at least 10 mV. For a positively charged Ion (i.e. Sodium), the AA increases the mV. For a negatively charged Ion (i.e. Sulfide, Fluoride, Chloride), the AA should decrease the mV. Start with a small trial. For example: Measure 50 mL of standard, add a magnetic stir bar and place on a stirrer, add ISA (consult ISE manual) and place ISE electrode tip into the sample. Put instrument in mV mode and record the observed mV. Using a micropipet, add a volume of the sample. Start by adding 500 μL at a time (for example). Watch the change in mV. When you have observed approximately a 10 mV change from the original standard, calculate the total volume added. Adjust sample and standard volumes proportionally to standard volumes that can be measured with accuracy. Use volumetric pipettes for standard, ISA and sample addition.
- Press **AA** then edit the procedure variables to the volumes determined in the prior step.

Procedure:

- Press **AA** to enter Analyte Addition mode.
- Volumetrically add standard to a clean beaker. Add Stir bar and place on a magnetic stir plate. Stir standard. The method will prompt user to add ISA. Place ISE sensor tip into the solution and a mV value will show on the display.
- Press **Continue** to take the first mV reading.
- When the reading is stable, press **Read** to store the first mV reading. The second step of the method will be displayed on the LCD, in which the user is notified to add the Sample Volume to the standard solution. The method parameters are also displayed on the LCD.
- Press **Continue** to take the second mV reading.
- When the reading is stable, press **Read** to store the second mV reading. The ISE measurement results will be displayed on the LCD.
- Press **Save** to log the current results into an ISE Method Report. Press **Direct Measure** to return to *ISE Measure* mode.
- Press **Start AA** to start another measurement. Rinse ISE sensor between samples.
- Press **Edit** to modify the method parameters.

Note: Press **Escape** at any time to stop the measurement and return to *ISE Measure* mode.

ANALYTE SUBTRACTION

To measure the concentration of a sample using Analyte Subtraction method:

- Press **MODE** and then **ISE** to select *ISE Measure* mode for the selected channel.
- Select the Analyte Subtraction method (see ISE Setup → Reading Mode).
- Prior to starting an AS procedure, the ISE sensor must be calibrated with a minimum of two standards containing ISA. The slope of the electrode will be used in all calculations involved in AS.
- If following an established procedure: Press **AS** then edit the method variables and follow the procedure below.
- Press **Edit** to set the method parameters. Press **Next** / **Previous** to select next/previous parameter to edit, then press **Edit** and use **Δ** or **▽** to set the desired parameter value. Press **Accept** to save the modified value and then press **Escape** to exit method parameters edit menu.
- If developing a procedure: Before attempting Analyte Subtraction analysis, it is important to determine which sample volume, Reactant volume and concentration, will produce the best results and the way the Reagent will react with the measured Ion on a molar basis (Stoichiometric factor). As a general rule, the reactant should contain the measured Ion so the sample addition will react with the Ion and reduce the measured concentration of the sample. The change of the mV value, before and after the sample addition, should be at least 10 mV. Start with a small trial. For example: Measure 50 mL of reactant, add a magnetic stir bar and place on a stirrer, add ISA (consult ISE manual) and place ISE electrode tip into the sample. Put instrument in mV mode and record the observed mV. Using a micropipet, add a volume of the sample. Start by adding 500 μL at a time (for example) Watch the change in mV. When you have observed approximately a 10 mV change from the original value, calculate the total volume added. Adjust sample and standard volumes proportionally to standard volumes that can be measured with accuracy. Use volumetric pipettes for standard, ISA and sample addition.
- Press **AS** then edit the procedure variables to the volumes determined in the prior step.

Procedure:


- Press **AS** to enter Analyte Subtraction mode.
- Volumetrically add Reactant to a clean beaker. Add Stir bar and place on a magnetic stir plate. Stir standard. The method will prompt user to add ISA. Place ISE sensor tip into the solution and a mV value will show on the display.
- Press **Continue** to take the first mV reading.
- When the reading is stable, press **Read** to store the first mV reading. The second step of the method will be displayed on the LCD in which the user is notified to add the Sample Volume to the standard solution.
- Press **Continue** to take the second mV reading.

- When the reading is stable, press **Read** to store the second mV reading. The ISE measurement results will be displayed on the LCD.
- Press **Save** to log the current results into an ISE Method Report. Press **Direct Measure** to return to *ISE Measure* mode. Press **Start AS** to start another measurement. Rinse sensor between samples.
- Press **Edit** to modify the method parameters.

Note: Press **Escape** at any time to stop the measurement and return to *ISE Measure* mode.

TEMPERATURE CALIBRATION

The user temperature calibration menu can be accessed during meter startup by simultaneously pressing three keys as shown in the drawing below. Press the keys after the short beep is heard at the meter power on. Keep all three keys pressed until Temp. Calibration menu appears.

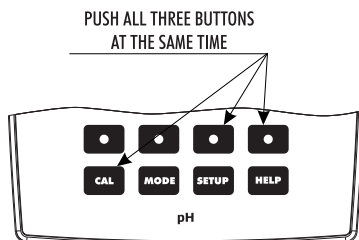


Note: The user temperature calibration is performed at three points: around 0 °C, 50 °C and 100 °C.

To perform the user temperature calibration:

- Select the desired temperature channel by pressing **Channel**.
- Press **Start User Calib** to start the temperature calibration. Adjust the temperature preset value using **Δ** or **▽** when necessary.
- Insert the temperature probe into the beaker with water at 0 °C.
- Wait for measurement to stabilize and then press **Accept** to confirm the calibration point.
- Repeat the previous steps for 50 °C and 100 °C.
- Save the calibration.
- Press **Escape** to return to measure mode.

Note: Press **Clear User Calib** if you want to clear the temperature user calibration.

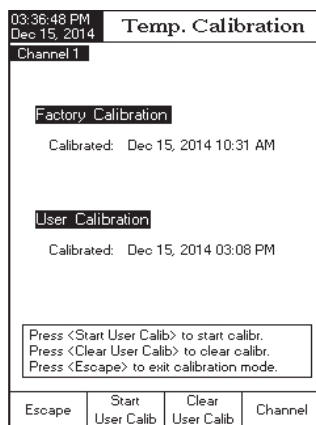



LOGGING

There are 5 ways the Reading Mode and Log may be configured together. The table below shows the combinations and indicates where the completed log will be stored.



Reading Mode	Log	log Recall
Direct	Automatic (1)	Automatic Log
	Manual (2)	Manual Log
	Auto Hold (NA)	Not Applicable
Direct/AutoHold	Automatic (3)	Automatic Log
	Manual (4)	Manual Log
	Auto Hold (5)	Manual Log


1) Direct Reading Mode and Automatic Log:

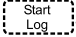
Real time continuous measurements are on display with continuous logging to meter memory. Press .

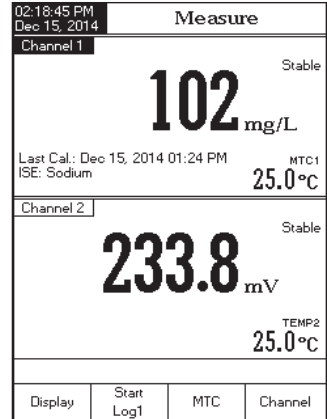

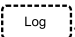



2) Direct Reading Mode and Manual Log:


Real time continuous measurements are on display and snapshots of measurement data are stored in the Manual log when the user presses . Subsequent snapshots will be added to the same Manual Log every time the  is depressed unless **New Lot** is selected under Log options.

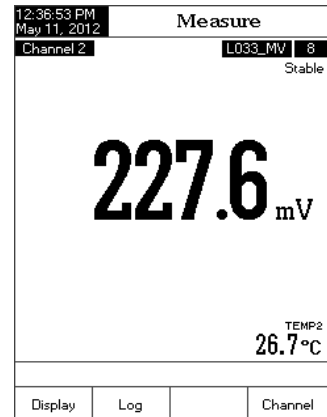Note: When the  is pressed the lot ID along with the current record number will appear for short time on the selected channel window on the top/left corner (e.g. L033_MV 8 - this means lot ID L033_mV and record number 8).



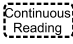

3) Direct/AutoHold Reading Mode and Automatic Log

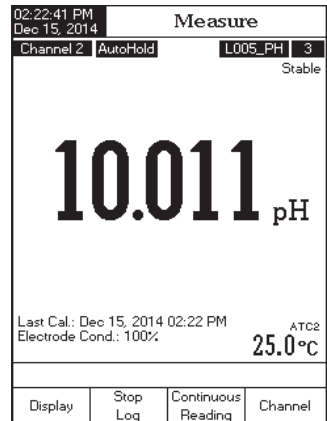



Press  and then  keys must be pressed on front display to initiate this function. Real time continuous measurements are on display with “AutoHold” flashing and real time continuous logging into meter memory, until the meter reaches the stability criteria to go into Auto Hold mode. The stored sample logs will be marked with an “H” to indicate the Auto Hold mode. The virtual key  returns operation to real time continuous measurements and  stops the logging session.



4) Direct/AutoHold Reading Mode and Manual Log

Press  in order to add one new record in the log report. The manual log is working even if it is in Auto Hold or Continuous reading mode. Press  to initiate the Auto Hold event. “AutoHold” will flash until the stability criteria is reached and then the screen freezes in Auto Hold mode, the data is marked with an “H”.


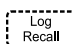
5) Direct/AutoHold Reading Mode and Auto Hold Log

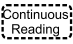
Press  and then  keys initiate and automate the capture of stable data which is stored in the Recall Manual Log file. During the process, “AutoHold” will flash until the stability criteria is reached and then the screen freezes in Auto Hold mode, the data is logged and marked with an “H”. The virtual key  returns operation to Real time continuous measurement. Press  again to log a second stable data point. The lot ID along with the record index will appear for short time on the top/left corner on the selected channel window, every time a record will be added to the lot.

LOG RECALL



This feature allows the user to view all stored data. If no data was logged, the “No records were found” message will be displayed on the LCD in the Log Recall screen. Otherwise, the instrument will display all the memorized lots in accordance with the selected option: Automatic Log, Manual Log or ISE Method Report (HI 5222 only).

To view the memorized data:

- Press  while in *Measure* mode.
- Press . The “Choose Log Report Type” message will be displayed in the Reminder messages area.

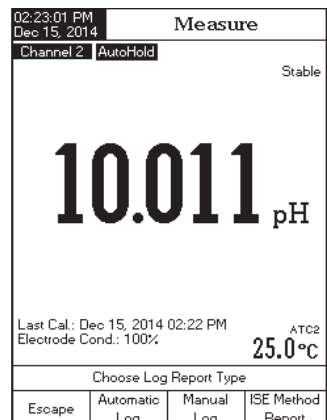


- Press **Automatic Log**, **Manual Log** or **ISE Method Report** to select the desired Log Report type. All logged lots for the selected Log Report type will be displayed on the LCD.
- To filter the displayed lots, press **MODE** and then the desired parameter **pH**, **mV** or **ISE** - HI 5222 only. Only the selected measurement parameter lots will be displayed on the LCD.
- Select the desired lot with **▲** or **▼** and press **View** to display the logged data from the highlighted lot. The "Please wait..." message will be displayed on the LCD for one second. The selected Logging Data Configuration options will be displayed on the LCD, together with GLP information (last calibration date and calibrated buffers/standards) if a calibration has been performed on the selected mode and the logged values (measured value, mV value, temperature value, temperature compensation mode and the logging time).

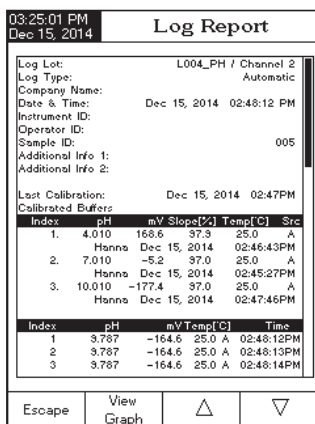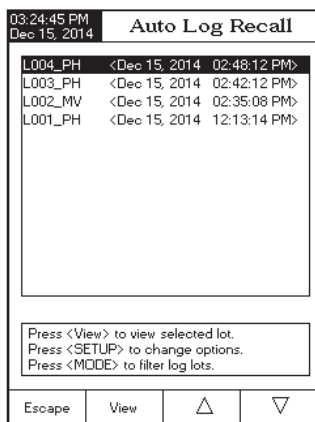

Note: For automatic logging only, it is possible to view the plotted graph.

- Press **View Graph** to display the graph.
- By pressing **Shift Axes** it is possible to move the graph along the X or Y axis with the arrow keys.
- If pressing **SETUP** while the graph is displayed, the zoom menu for the X and Y axes will be accessed. Press **Zoom Time** or **Zoom pH** / **Zoom mV** / **Zoom Rel.mV** / **Zoom Ion** to switch between the active zooming axes and then zoom in or out on the selected axis by pressing the appropriate virtual key.
- Press **Escape** to return to the previous menu at any time.

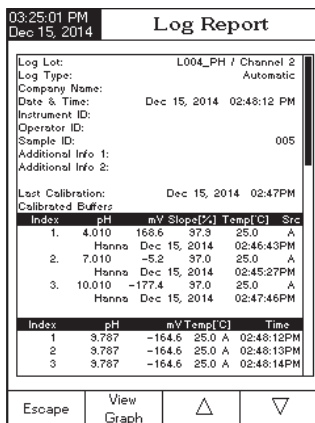

To delete lots:

- Press **SETUP** while in *Log Recall* mode.
- Press **Delete** or **Delete All** to access delete or delete all mode. Otherwise, press **View** to return to *Log Recall* view mode.
- After selecting one of the deleting modes, use **Δ** or **∇** to select one lot and then press **Delete** or **Delete All** to delete the selected lot or all lots. The “Please wait...” message will be displayed on the LCD until the selected lot or all lots are deleted.
- Press **SETUP** and then press **View** to exit deleting mode and return to *Log Recall* view mode.
- Press **Escape** to exit *Log Recall* mode and return to *Measure* mode.

Note: Logged lots should also be deleted whenever “Please Delete Old Log Files” or “Low Data Logging Space” message appears on the LCD, in the Reminder messages area.

03:26:12 PM Dec 15, 2014		Auto Log Recall	
L004_PH	<Dec 15, 2014 02:48:12 PM>		
L003_PH	<Dec 15, 2014 02:42:12 PM>		
L002_MV	<Dec 15, 2014 02:35:08 PM>		
L001_PH	<Dec 15, 2014 12:13:14 PM>		
Press <View> to select view mode. Press <Delete> for delete mode. Press <Delete All> for delete all mode.			
	View	Delete	Delete All

PC INTERFACE

Data transmission from the instrument to the PC can be done with the **HI 92000** Windows® compatible software (optional). **HI 92000** also offers graphing and on-line help features.

Data logged on the **HI 5221** and **HI 5222** meters can be exported to the most popular spreadsheet programs for further analysis.

HI 5221 and **HI 5222** instruments have an USB interface available.

Use a standard USB cable to connect your instrument to the PC.

Make sure that the instrument and the **HI 92000** software have the same baud rate and the appropriate communication port.

The PC software may also be used for real time logging.

ADDITIONAL INFORMATION

ISE THEORY

An Ion Selective Electrode (ISE) is an electrochemical sensor that changes voltage with the activity or concentration of ions in solutions. The change in voltage is a logarithmic relationship with concentration, and is expressed by the Nernst equation:

$$E = E^{\circ} + S \log(a)$$

where: E - the measured voltage;

E° - standard voltage and other standard system voltages;

a - the activity of the Ion being measured;

$$S = \frac{2.303RT}{nF}$$

S - the Nernst slope factor and is derived from thermodynamic principles:

R - the universal gas constant (8.314 J/Kmol);

T - the temperature in degrees Kelvin;

F - the Faraday's constant (96,485 C/mol);

n - the Ion charge.

The slope may be positive or negative depending upon the Ion charge (n).

SPECIES	SLOPE (mV/decade)
Monovalent cation	+59.16
Monovalent anion	-59.16
Divalent cation	+29.58
Divalent anion	-29.58

Activity and concentration are related by an "activity coefficient", expressed as:

$$a = \gamma \cdot C$$

where: a - the activity of the Ion being measured;

γ - the activity coefficient;

C - the concentration of the Ion being measured.

In very dilute solutions γ approaches 1 so activity and concentration are the same.

Actual samples that are more concentrated have much smaller activity coefficients ($\gamma < 1$). The addition of an inert background salt to standards and samples stabilizes the activity coefficient so that concentration measurements may be made directly. Some of Hanna's Ionic Strength Adjuster formulations also may optimize pH and complex interferences, in addition to standardizing the ionic strength.

The Nernst equation can be rewritten:

$$E = E^o + S \log(C)$$

ION SELECTIVE ANALYSIS METHODS

Direct Analysis

This method is a simple procedure for measuring multiple samples. It should only be used in the linear working regions of the sensor. A direct reading instrument such as the **HI 5222** determines concentration of the unknown by a direct reading after calibrating the instrument with the standards. The instrument is calibrated as described in "ISE CALIBRATION & MEASUREMENTS" section, with two or more freshly made standards that are in the measurement range of the unknowns. Ionic strength adjustment is made to samples and standards. Unknowns are measured directly by the instrument.

At lower concentrations, in non-linear regions of the electrode response, multiple calibration points will extend measurements to a practical detection limit. Calibrations must be performed more frequently in these cases.

Incremental Methods

Incremental methods are useful for the measurement of samples whose constituents are variable or concentrated. Incremental techniques can reduce errors from such variables as temperature, viscosity, or pH extremes and will provide indirect analysis of ions for which there is no ISE sensor for a direct measurement. There are four commonly used different incremental methods for sample measurement. They are Known Addition, Known Subtraction, Analyte Addition and Analyte Subtraction. **HI 5222** allows the analyst to use these techniques as a simple routine procedure, thus eliminating calculations or tables. The method once set up can be used for repetitive measurements on multiple samples.

Known Addition and Known Subtraction

With Known addition, *standard is added* to a sample being measured. The standard and sample contain the same ion. mV are taken before and after the standard addition. From the change in mV, the sample concentration is determined.

$$C_{SAMP} = \frac{C_{SAMP} \cdot V_{STD}}{(V_{SAMP} + V_{STD} + V_{ISA}) \cdot 10^{\frac{\Delta E}{S}} - (V_{SAMP} + V_{ISA})} \cdot \frac{(V_{SAMP} + V_{ISA})}{V_{SAMP}}$$

With Known subtraction, *a known standard is added to an ionic sample* being measured. The standard reacts with the measured ion in the sample in a known manner, thus removing measured ions from the solution. From the change in mV, the concentration of the sample is determined.


$$C_{SAMP} = \frac{C_{STD} \cdot V_{STD} \cdot f}{(V_{SAMP} + V_{ISA}) - (V_{SAMP} + V_{STD} + V_{ISA}) \cdot 10^{\frac{\Delta E}{S}}} \cdot \frac{(V_{SAMP} + V_{ISA})}{V_{SAMP}}$$

where: C_{SAMP} - the sample concentration;
 C_{STD} - the standard concentration;
 V_{SAMP} - the sample volume;
 V_{STD} - the standard volume;
 V_{ISA} - ISA volume

ΔE - the difference of potential from the electrode;
 S - the electrode slope, determined in a previous calibration;
 f - the stoichiometric ratio between sample and standard;

Example 1


You have sulfide samples and you are adding Ag^+ . The reaction is:

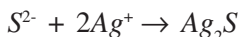

One mole sulfide sample reacts with 2 moles silver standard ($f = 1/2$).

Example 2

You have sulfide samples and you are adding Pb^{2+} . The reaction is:

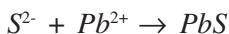

One mole sulfide sample reacts with 1 mole lead standard ($f = 1$).

Analyte Addition and Analyte Subtraction

Analyte Addition and Subtraction are variations of the previous two methods.

With Analyte Addition, sample (analyte) is added to an Ion standard being measured. The standard and sample contain the same Ion. mV are taken before and after the sample addition. From the mV the analyte concentration is determined.

$$C_{SAMP} = \frac{C_{STD} \cdot V_{STD}}{(V_{STD} + V_{ISA})} \cdot \frac{(V_{STD} + V_{SAMP} + V_{ISA}) \cdot 10^{\frac{\Delta E}{S}} - (V_{STD} + V_{ISA})}{V_{SAMP}}$$

With Analyte Subtraction, sample (analyte) is added to an Ion standard being measured. The analyte reacts with the measured Ion in a known manner thus removing measured ions from the solution. From the change in mV the concentration of the analyte is determined.

$$C_{SAMP} = f \cdot \left\{ \frac{(V_{STD} + V_{ISA})}{V_{SAMP}} - \left[1 + \frac{(V_{STD} + V_{ISA})}{V_{SAMP}} \right] \cdot 10^{\frac{\Delta E}{S}} \right\} \cdot \left(\frac{C_{STD} \cdot V_{STD}}{V_{STD} + V_{ISA}} \right)$$

where: C_{SAMP} - the sample concentration;
 C_{STD} - the standard concentration;
 V_{SAMP} - the sample volume;
 V_{STD} - the standard volume;
 V_{ISA} - ISA volume;

ΔE - the difference of potential from the electrode;
 S - the electrode slope, determined in a previous calibration;
 f - the stoichiometric ratio between sample and standard;

pH BUFFER TEMPERATURE DEPENDENCE


Temperature has an effect on pH. The calibration buffer solutions are affected by temperature changes to a lower degree than normal solutions.

During calibration the instrument will automatically calibrate to the pH value corresponding to the measured or set temperature.

TEMP			pH BUFFERS							
°C	°K	°F	1.679	3.000	4.010	6.862	7.010	9.177	10.010	12.454
0	273	32	1.670	3.072	4.007	6.982	7.130	9.459	10.316	13.379
5	278	41	1.670	3.051	4.002	6.949	7.098	9.391	10.245	13.178
10	283	50	1.671	3.033	4.000	6.921	7.070	9.328	10.180	12.985
15	288	59	1.673	3.019	4.001	6.897	7.046	9.273	10.118	12.799
20	293	68	1.675	3.008	4.004	6.878	7.027	9.222	10.062	12.621
25	298	77	1.679	3.000	4.010	6.862	7.010	9.177	10.010	12.450
30	303	86	1.683	2.995	4.017	6.851	6.998	9.137	9.962	12.286
35	308	95	1.688	2.991	4.026	6.842	6.989	9.108	9.919	12.128
40	313	104	1.693	2.990	4.037	6.837	6.983	9.069	9.881	11.978
45	318	113	1.700	2.990	4.049	6.834	6.979	9.040	9.847	11.834
50	323	122	1.707	2.991	4.062	6.834	6.978	9.014	9.817	11.697
55	328	131	1.715	2.993	4.076	6.836	6.979	8.990	9.793	11.566
60	333	140	1.724	2.995	4.091	6.839	6.982	8.969	9.773	11.442
65	338	149	1.734	2.998	4.107	6.844	6.987	8.948	9.757	11.323
70	343	158	1.744	3.000	4.123	6.850	6.993	8.929	9.746	11.211
75	348	167	1.755	3.002	4.139	6.857	7.001	8.910	9.740	11.104
80	353	176	1.767	3.003	4.156	6.865	7.010	8.891	9.738	11.003
85	358	185	1.780	3.002	4.172	6.873	7.019	8.871	9.740	10.908
90	363	194	1.793	3.000	4.187	6.880	7.029	8.851	9.748	10.819
95	368	203	1.807	2.996	4.202	6.888	7.040	8.829	9.759	10.734

During calibration the instrument will display the pH buffer value at 25 °C.

ELECTRODE CONDITIONING AND MAINTENANCE

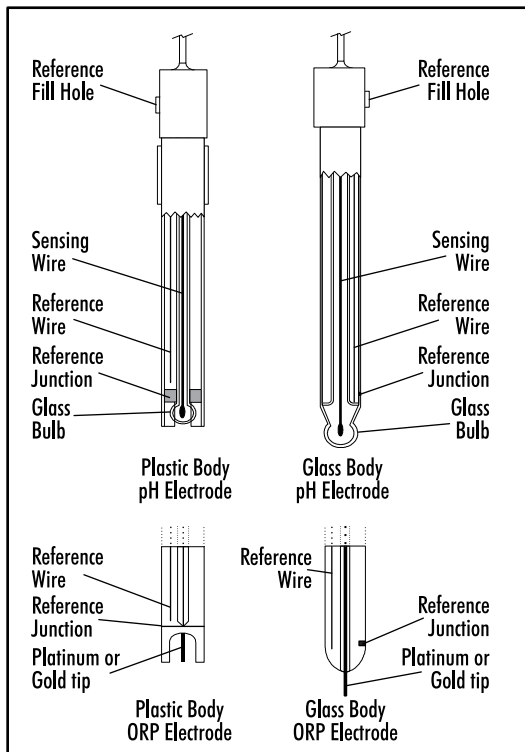

PREPARATION PROCEDURE

Remove the protective cap off the pH electrode.

SALT DEPOSITS MAY BE PRESENT. This is normal with electrodes. They will disappear when rinsed with water. During transport, tiny bubbles of air may form inside the glass bulb affecting proper functioning of the electrode. These bubbles can be removed by “shaking down” the electrode as you would do with a glass thermometer.

If the bulb and/or junction is dry, soak the electrode in **HI 70300** or **HI 80300** Storage Solution for at least one hour.

For refillable electrodes:

If the filling solution (electrolyte) is more than 2” cm (1”) below the fill hole, add **HI 7082** or **HI 8082** 3.5M KCl Electrolyte Solution for double junction or **HI 7071** or **HI 8071** 3.5M KCl+AgCl Electrolyte Solution for single junction electrodes.

Unscrew the fill hole screw during measurements. This will allow electrolyte to flow out of the junction.

For AmpHel® electrodes:

If the electrode does not respond to pH changes, the battery may have run down and the electrode or battery (if replaceable) should be replaced.

MEASURE

Rinse the pH electrode tip with distilled water. Immerse the tip bottom 4 cm (1.5") in the sample and stir gently for a few seconds.

For a faster response and to avoid cross-contamination of the samples, rinse the electrode tip with a few drops of the solution to be tested, before taking measurements.

STORAGE PROCEDURE

To minimize clogging and ensure a quick response time, the glass bulb and the junction of pH electrode should be kept moist and not allowed to dry out.

Replace the solution in the protective cap with a few drops of **HI 70300** or **HI 80300** Storage Solution or, in its absence, Filling Solution (**HI 7071** or **HI 8071** for single junction and **HI 7082** or **HI 8082** for double junction electrodes). Follow the Preparation Procedure before taking measurements.

Note: NEVER STORE THE ELECTRODE IN DISTILLED OR DEIONIZED WATER.

PERIODIC MAINTENANCE

Inspect the electrode and the cable. The cable used for connection to the instrument must be intact and there must be no points of broken insulation on the cable or cracks on the electrode stem or bulb. Connectors must be perfectly clean and dry. If any scratches or cracks are present, replace the electrode. Rinse off any salt deposits with water.

pH PROBE MAINTENANCE

For refillable electrodes:

Refill the reference chamber with fresh electrolyte (**HI 7071** or **HI 8071** for single junction or **HI 7082** or **HI 8082** for double junction electrodes). Allow the electrode to stand upright for 1 hour. Follow the Storage Procedure above.

pH CLEANING PROCEDURE

- *General* Soak in Hanna **HI 7061** or **HI 8061** General Cleaning Solution for approximately 30 minutes.
- *Protein* Soak in Hanna **HI 7073** Protein Cleaning Solution for 15 minutes.
- *Inorganic* Soak in Hanna **HI 7074** Inorganic Cleaning Solution for 15 minutes. This is good at cleaning a black ceramic junction.
- *Oil/grease* Rinse with Hanna **HI 7077** or **HI 8077** Oil and Fat Cleaning Solution.


IMPORTANT: After performing any of the cleaning procedures, rinse the electrode thoroughly with distilled water, refill the reference chamber with fresh electrolyte (not necessary for gel-filled electrodes) and soak the electrode in **HI 70300** or **HI 80300** Storage Solution for at least 1 hour before taking measurements.

TROUBLESHOOTING GUIDE

SYMPTOMS	PROBLEM	SOLUTION
Slow response/excessive drift.	Dirty pH electrode. Dirty reference junction.	Soak the electrode tip in HI 7061 solution for 30 minutes and then clean the electrode. Soak in HI 7074 .
Readings fluctuate up and down (noise).	Clogged/dirty junction. Low electrolyte level (refillable electrodes only).	Clean the electrode. Refill with fresh solution (for refillable electrodes only).
The LCD displays "----" during measurements (pH, mV, mV Rel or ISE).	Out of range in the appropriate scale.	Check to see if the pH bulb and reference junction are in solution. Make sure the sample is in the specified range. Check the electrolyte level and the general state of the pH/ORP or ISE electrode.
Out of range in the mV scale.	Dry reference junction.	Soak in HI 70300 Storage solution for at least one hour.
The instrument does not work with the temperature probe.	Out of order temperature probe.	Replace the probe.
The meter fails to calibrate or gives faulty readings.	Broken or out of order electrode.	Replace the electrode.
Explicit warnings are displayed during calibration.	Dirty/broken electrode, contaminated buffers.	Follow displayed instructions.
The electrode condition is not displayed after calibration.	Only one-point calibration has been performed.	Perform at least a two-point calibration.
The instrument does not start up.	Internal or software error.	Restart the instrument using the power switch or unplug adapter. If the error persists, contact your vendor.

TEMPERATURE CORRELATION FOR pH SENSITIVE GLASS

The resistance of glass electrodes partially depends on the temperature. The lower the temperature, the higher the resistance. It takes more time for the reading to stabilize if the resistance is higher.

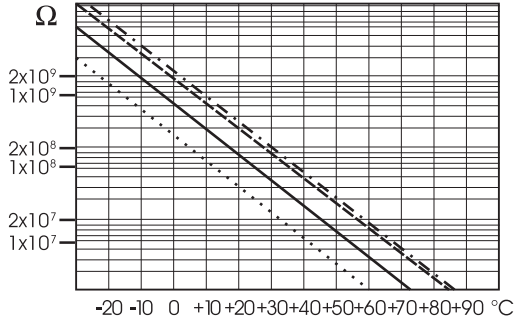

Since the resistance of the pH electrode is in the range of 50 – 200 Mohm, the current across the membrane is in the pico Ampere range. Large currents can disturb the calibration of the electrode for many hours.

The pH electrode's life also depends on the temperature. If constantly used at high temperatures, the electrode life is drastically reduced.

Typical Electrode Life

Ambient Temperature	1 – 3 years
90 °C (194 °F)	Less than 4 months
120 °C (248 °F)	Less than 1 month

Alkaline Error

High concentrations of sodium ions interfere with readings in alkaline solutions. The pH at which the interference starts to be significant depends upon the composition of the glass. This interference is called alkaline error and causes the pH to be underestimated.

ACCESSORIES

pH BUFFER SOLUTIONS

HI 6016	pH 1.679 Buffer Solution, 500 mL bottle
HI 6003	pH 3.000 Buffer Solution, 500 mL bottle
HI 8004L	pH 4.01 Buffer Solution in FDA approved bottle, 500 mL
HI 6004	pH 4.010 Buffer Solution, 500 mL bottle
HI 8006L	pH 6.86 Buffer Solution in FDA approved bottle, 500 mL
HI 6068	pH 6.862 Buffer Solution, 500 mL bottle
HI 8007L	pH 7.01 Buffer Solution in FDA approved bottle, 500 mL
HI 6007	pH 7.010 Buffer Solution, 500 mL bottle
HI 8009L	pH 9.18 Buffer Solution in FDA approved bottle, 500 mL
HI 8010L	pH 10.01 Buffer Solution in FDA approved bottle, 500 mL
HI 6010	pH 10.010 Buffer Solution, 500 mL bottle
HI 6124	pH 12.450 Buffer Solution, 500 mL bottle

ELECTRODE STORAGE SOLUTIONS

HI 70300L	Storage Solution, 500 mL bottle
HI 80300L	Storage Solution in FDA approved bottle, 500 mL

ELECTRODE CLEANING SOLUTIONS

HI 70000P	Electrode Rinse Sachets, 20 mL, 25 pcs
HI 7061L	General Purpose Solution, 500 mL bottle
HI 7073L	Protein Cleaning Solution, 500 mL bottle
HI 7074L	Inorganic Cleaning Solution, 500 mL bottle
HI 7077L	Oil & Fat Cleaning Solution, 500 mL bottle
HI 8061L	General Purpose Solution in FDA approved bottle, 500 mL
HI 8073L	Protein Cleaning Solution in FDA approved bottle, 500 mL
HI 8077L	Oil & Fat Cleaning Solution in FDA approved bottle, 500 mL

ELECTRODE REFILL ELECTROLYTE SOLUTIONS

HI 7071	3.5M KCl + AgCl Electrolyte, 4x30 mL, for single junction electrodes
HI 7072	1M KNO ₃ Electrolyte, 4x30 mL
HI 7082	3.5M KCl Electrolyte, 4x30 mL, for double junction electrodes
HI 8071	3.5M KCl + AgCl Electrolyte in FDA approved bottle, 4x30 mL, for single junction electrodes
HI 8072	1M KNO ₃ Electrolyte in FDA approved bottle, 4x30 mL
HI 8082	3.5M KCl Electrolyte in FDA approved bottle, 4x30 mL, for double junction electrodes
HI 8093	1M KCl + AgCl Electrolyte in FDA approved bottle, 4x30 mL

ORP SOLUTIONS

HI 7020L Test Solution 200-275 mV, 500 mL bottle

HI 7021L Test Solution 240 mV, 500 mL bottle


HI 7022L Test Solution 470 mV, 500 mL bottle

HI 7091L Reducing Pretreatment Solution, 500 mL

HI 7092L Oxidizing Pretreatment Solution, 500 mL

pH ELECTRODES


All electrodes part numbers ending in B are supplied with a BNC connector and 1 m (3.3') cable, as shown below:

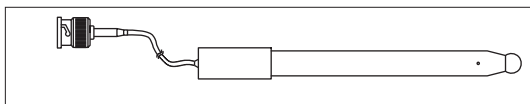

HI 1043B

Glass body, double junction, refillable, combination pH electrode.


Use: strong acid/alkali.

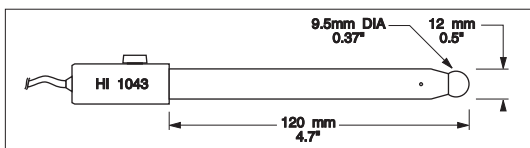

HI 1053B

Glass body, triple ceramic, conical shape, refillable, combination pH electrode.


Use: emulsions.

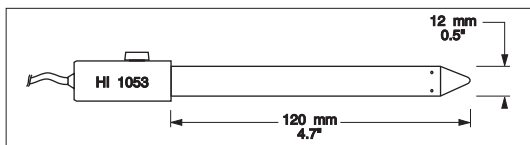

HI 1083B

Glass body, micro, Viscolene, non refillable, combination pH electrode.


Use: biotechnology, micro titration.

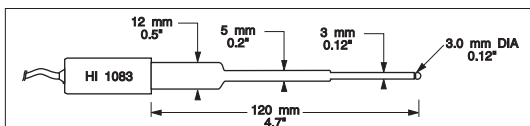

HI 1131B

Glass body, refillable, double junction, combination pH electrode.


Use: general purpose.

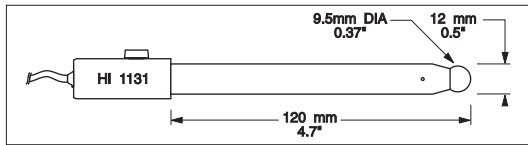

HI 1330B

Glass body, semimicro, single junction, refillable, combination pH electrode.


Use: laboratory, vials.

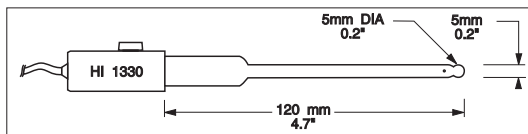

HI 1331B

Glass body, semimicro, single junction, refillable, combination pH electrode.


Use: flasks.

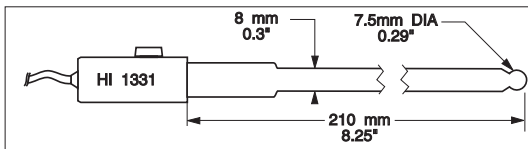

HI 1230B

Plastic body (PEI), double junction, gel filled, combination pH electrode.


Use: general, field.

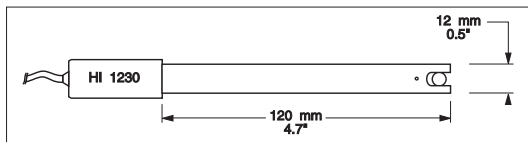

HI 2031B

Glass body, semimicro, conical, refillable, combination pH electrode.


Use: semisolid products.

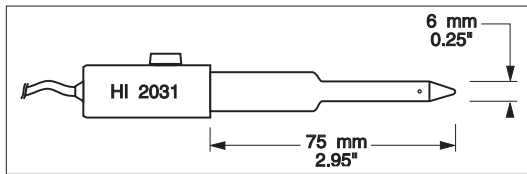

HI 1332B

Plastic body (PEI), double junction, refillable, combination pH electrode.


Use: general purpose.

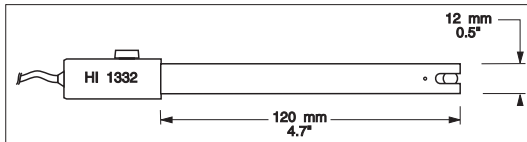

HI 1413B

Glass body, single junction, flat tip, Viscolene, non refillable, combination pH electrode.


Use: surface measurement.

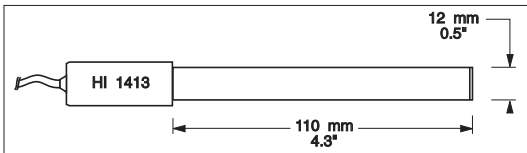

FC 100B

Plastic body (PVDF), double junction, refillable, combination pH electrode.


Use: general purpose for food industry.

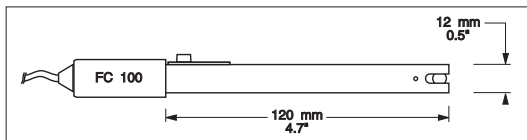

FC 200B

Plastic body (PVDF), open junction, conical, Viscolene, non refillable, combination pH electrode.


Use: meat & cheese.

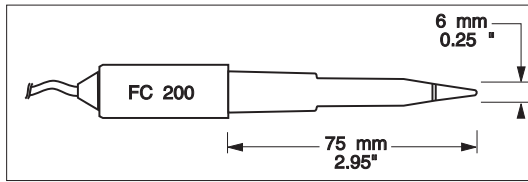

FC 210B

Glass body, double junction, conical, Viscolene, non refillable, combination pH electrode.


Use: milk, yogurt.

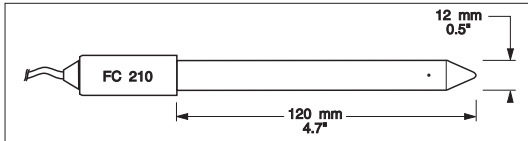

FC 220B

Glass body, triple ceramic, single junction, refillable, combination pH electrode.


Use: food processing.

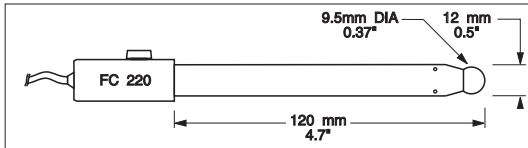

FC 911B

Plastic body (PVDF), double junction, refillable with built-in amplifier, combination pH electrode.

Use: very high humidity.

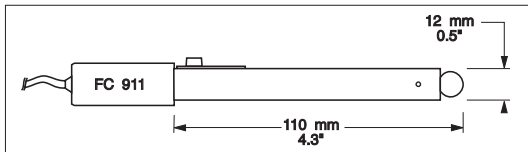


ORP ELECTRODES

HI 3131B

Glass body, refillable, combination platinum ORP electrode.


Use: titration.

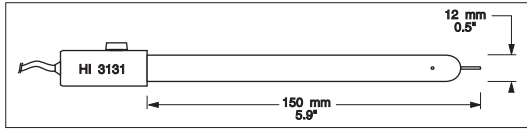

HI 3230B

Plastic body (PEI), gel filled, combination platinum ORP electrode.


Use: general purpose.

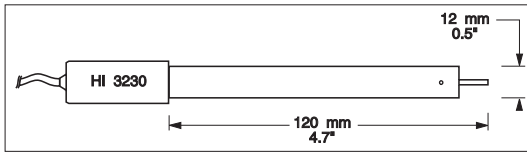

HI 4430B

Plastic body (PEI), gel filled, combination gold ORP electrode.

Use: general purpose.

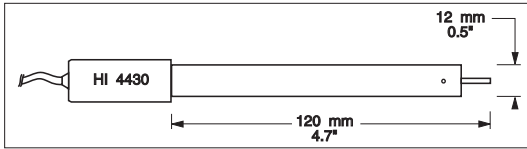


Consult the Hanna General Catalog for more electrodes with screw-type or BNC connectors.

EXTENSION CABLE FOR SCREW-TYPE ELECTRODES (SCREW TO BNC ADAPTER)

HI 7855/1 Extension cable 1 m (3.3') long

HI 7855/3 Extension cable 3 m (9.9') long

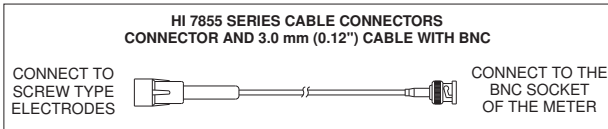

OTHER ACCESSORIES

HI 710005/8	Voltage adapter from 115 Vac / 12 Vdc 800 mA (USA plug)
HI 710006/8	Voltage adapter from 230 Vac / 12 Vdc 800 mA (European plug)
HI 76404W	Electrode holder
HI 8427	pH and ORP electrode simulator with 1 m (3.3') coaxial cable ending in female BNC connectors
HI 931001	pH and ORP electrode simulator with LCD and 1 m (3.3') coaxial cable ending in female BNC connectors
HI 7662-W	Temperature probe with 1 m (3.3') cable
HI 92000	Windows® compatible software
HI 920013	USB cable

RECOMMENDATIONS FOR USERS

Before using these products, make sure they are entirely suitable for the environment in which they are used. Operation of these instruments in residential areas could cause unacceptable interferences to radio and TV equipment, requiring the operator to follow all necessary steps to correct interferences.

The glass bulb at the end of the pH electrode is sensitive to electrostatic discharges. Avoid touching this glass bulb at all times.

Any variation introduced by the user to the supplied equipment may degrade the instruments' EMC performance.

To avoid electrical shock, do not use these instruments when voltages at the measurement surface exceed 24 Vac or 60 Vdc.

To avoid damage or burns, do not perform any measurement in microwave ovens.

Hanna Instruments reserves the right to modify the design, construction or appearance of its products without advance notice.



Hanna Instruments Inc.
Highland Industrial Park
584 Park East Drive
Woonsocket, RI 02895 USA

Technical Support for Customers

Tel. (800) 426 6287
Fax (401) 765 7575
E-mail tech@hannainst.com
www.hannainst.com

Local Sales and Customer Service Office

