


# Dino-Lite Edge 5216 Series Premier 5116 Series Quick Start

## System Configuration

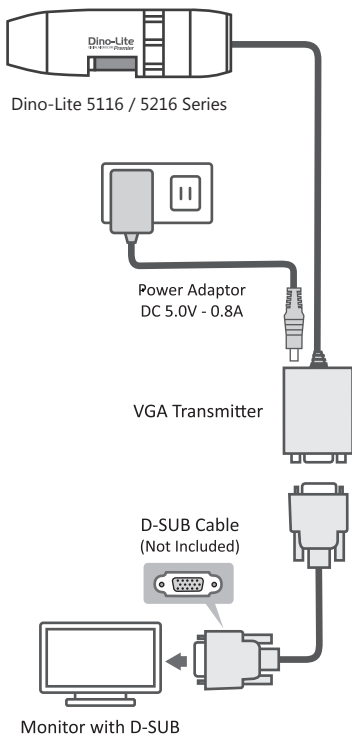

## How to Use

### NOTICE

- The higher temperature of the top metal part is normal. Please avoid touching it when in use.
- It is recommended to use Dino-Lite in a well-ventilated environment.
- Be sure to use the supplied power adaptor connecting to Dino-Lite for preventing any damage to the device.

- Use the adjustment knob (see Fig. 1) to obtain a focus or magnification with the Dino-Lite.
- Tap the MicroTouch (see Fig. 1) to freeze /unfreeze frame.
- Tap and hold the MicroTouch (see Fig. 1) for 2 seconds to switch the LEDs.
- Slide the Scroll Lock (see Fig. 1) lever down, if want to fix the position of the adjustment knob.

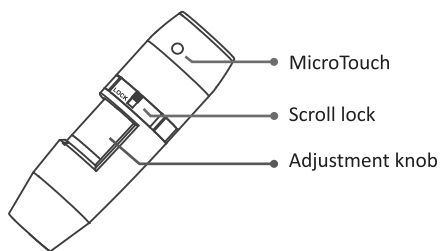


Fig.1: Location of Dino-Lite Controls