WCT

Kältefalle

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Cold Trap Bath

Operating Manual

gilt für / written for
WCT -40/ -80

witeg Labortechnik GmbH

(English manual at page 12)

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Contents

1 Introduction	13
2 Safety Instructions	13
2.1 Description of symbols	13
2.2 Instructions.....	13
3 Product description	15
3.1 Key features.....	15
3.2 General survey	15
3.3 Operating surface	16
3.4 Term description	16
4 Packing contents	16
5 Positioning the device and commissioning	17
6 Job description	17
6.1 Menu structure	17
6.2 Set turn-off delay.....	18
6.3 Set turn-on delay.....	18
6.4 Set temperature.....	18
6.5 Set Offset	18
6.6 Key lock.....	19
6.7 Storage function.....	19
7 Maintenance and cleaning	19
8 Troubleshooting Guide	20
9 Technical data	21
10 Certificate	22

1 Introduction

Thank you for purchasing our **WCT Cold Trap Barh**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to low temperatures**

Instructions to avoid injuries caused by low temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the usage.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.

- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Warn all attendees if you leave the device unattended.
- Do not clean the scorching or frozen device.
- Pay attention and only work with gloves, otherwise you may freeze on the device's metal parts.
- Do not put thermal-sensitive items at the device.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface to avoid vibrations.
- If you set the device in a site with vibrations, make sure that the device can not slip away.
- The device could get damaged if the vibrations are too strong.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superior to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overrefrigerator, short circuits or fires.
- Make sure the device stands upright within the last 24 hours.
- Do not tilt the device.

- Do not use this device in area with potentially explosive atmosphere.

- Reparations and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty.

3 Product description

3.1 Key features

The **Wisdom WCT Cold Trap Bath** is designed for chemical, biological, medical and pharmaceutical applications. The device was developed for simplest usage and a very long life.


Key features are as follows:

Simplest operating by MODE button and Jog-Shuttle knob.


- Fast commissioning.
- Long life by simply and solid construction.
- Temperature and delay time showed on illuminated LCD.
- Storage functions for set time and temperature values.
- Turn-off an -on delay from one minute up to over 99 hours in one minute steps adjustable.
- Signal sound after end of turn-off delay.
- Easy to justify.
- Key lock against inadvertent changing of set speed and timer values.
- A lot of laboratory devices are mountable very quickly.

3.2 General survey

The functionality of the described components is the same at each type of the **WCT**-series. Only the position may differ from type to type.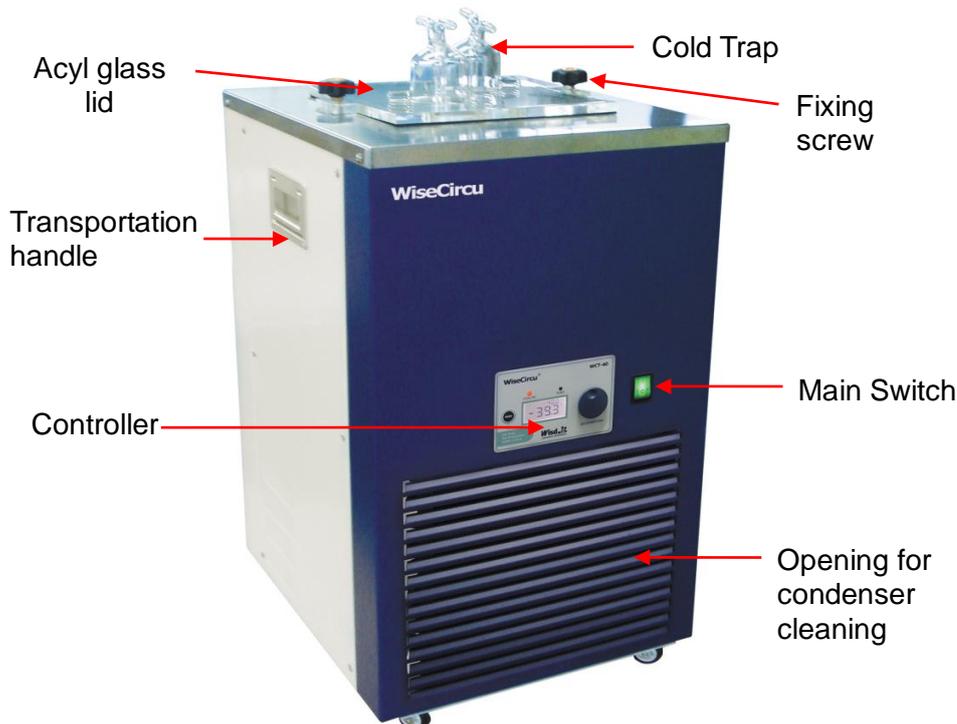

3.3 Operating surface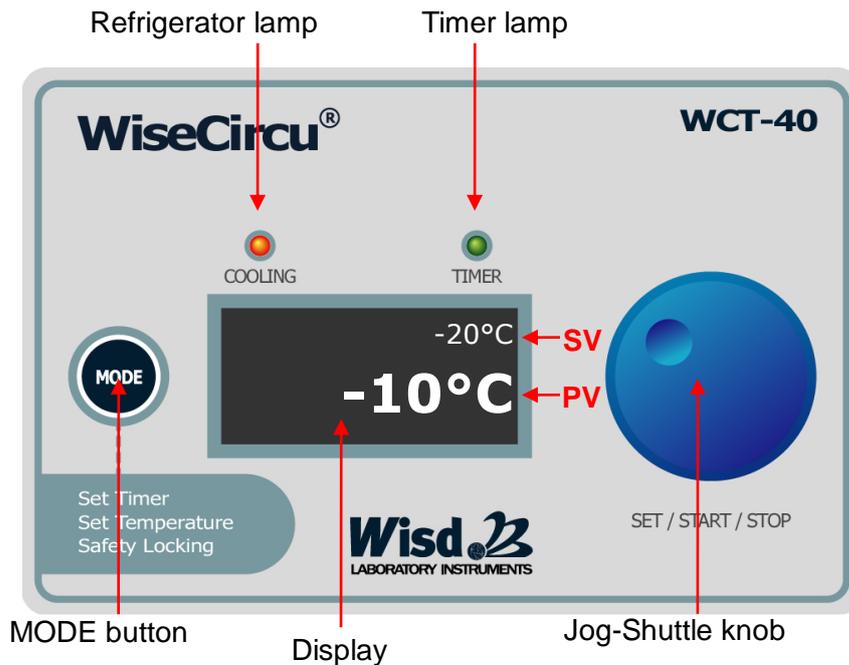

3.4 Term description

Main switch

- On / off switch.

Timer lamp

- Is blinking in menu items for timer modulation
- Shines continuous in the key lock menu item when timer runs.

Refrigerator lamp

- Is blinking in menu items for temperature modulation
- Shines continuous in the key lock menu item when refrigerator runs.

Jog-Shuttle knob

- Change values by turning.
- Push the knob for start and stop or probate set values.

MAIN button

- For changing between the different menu items.

Key combination

- Push and hold the MODE button first and in addition the Jog-Shuttle knob to reach the sub menu items, then release both buttons.

4 Packing contents

quantity	denomination
1	WCT Cold Trap Bath
1	Power supply cord
1	Stainless steel lid
1	Acrylic glass lid
2	Fixing screw
2	Glass made cold trap
1	Operating manual


5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package and perhaps soiling carefully.
3. Do not tilt the device. The cooling agent must not disperse in the cooling unit, otherwise the device could get damaged during the operation. Make sure that the device stands upright for the last 12 hours.
4. If there are any unexpected transport damages, please inform your forwarder and your customer service.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Fill the operation liquid in. The cooling coil must be covered completely.
7. Make sure that you follow all safety instructions.
8. Connect the device with the power outlet.
9. Switch the device on and start the process (look at job description).

6 Job description

6.1 Menu structure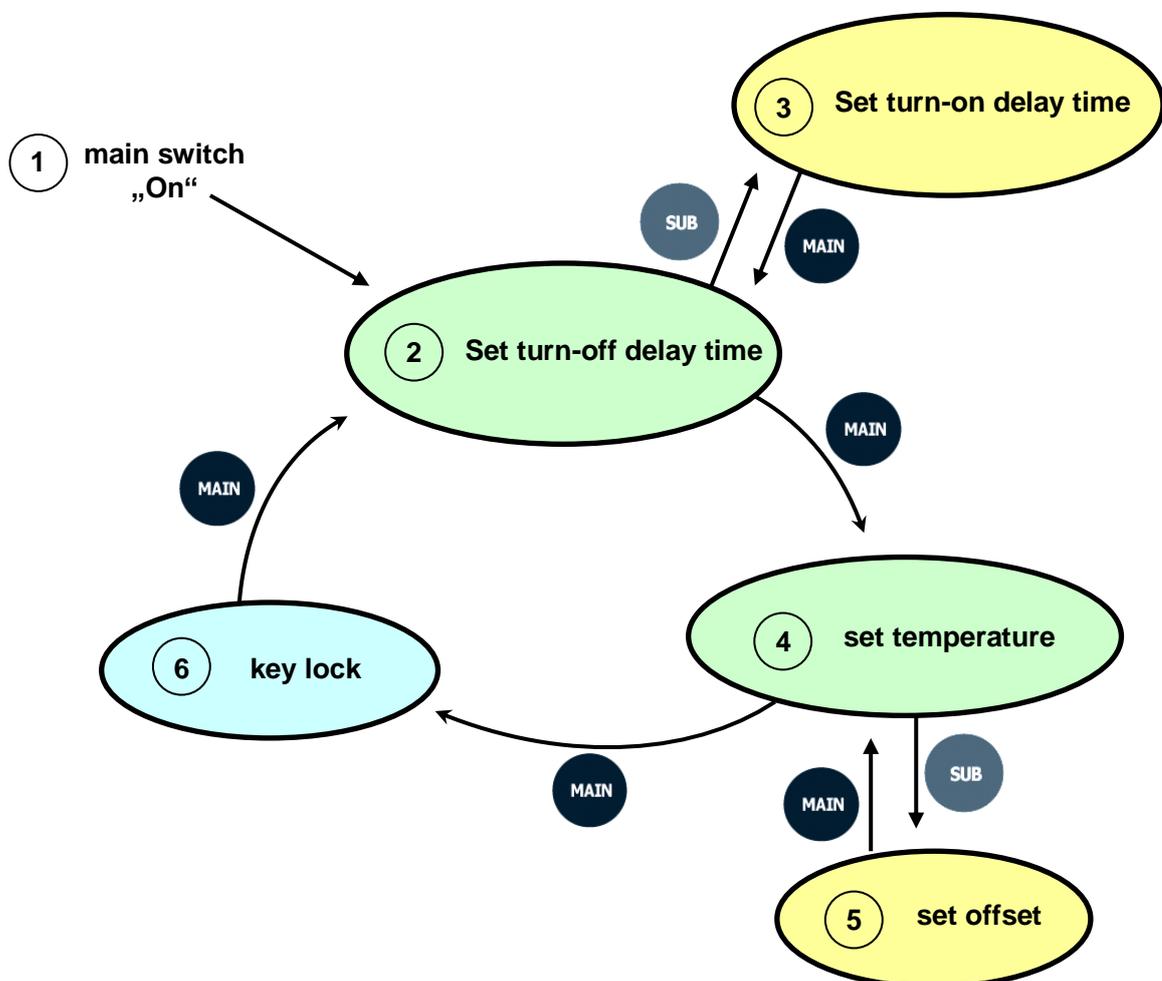

6.2 Set turn-off delay

The turn-off delay timer serves as automatically stop function. It turns off the refrigerator after the timeframe you set is elapsed.

- First menu item after switching the device on is the menu “set turn-off delay”. The timer lamp is blinking.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- “00:00” stands for endurance run.
- Turn the Jog-Shuttle to choose your desired value and probate it by pushing the Jog-Shuttle. The delay time starts to run by probating the value.
- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.

6.3 Set turn-on delay

The turn-on delay timer serves as automatically start function. It turns on the refrigerator after elapsing of the timeframe you set.

- To reach this sub menu item you have use the key combination (look at 3.4) in the menu “set turn-off delay”. **LCD shows “d-tr”**.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Shuttle to choose your desired value and probate it by pushing the Jog-Shuttle. The delay time starts to run by probating the value.
- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.
- Turn the Shuttle knob clockwise to increase the time and counter-clockwise to reduce it.
- The turn-off delay time will not elapse before the turn-on delay time is executed.

6.4 Set temperature

- To reach this menu item you have to push the MAIN button in the menu “set turn-off delay”
- The temperature is adjustable in 0,1 °C-steps.
- Turn the Jog-Shuttle to choose your desired value and probate it by pushing the Jog-Shuttle. The delay time starts to run by probating the value.
- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.
- Turn the Shuttle knob clockwise to increase the temperature and counter-clockwise to reduce it.
- The refrigerator lamp shines, if the heater is activated.

6.5 Set Offset

If the temperature measured by you differs from the temperature the device is showing (PV) you have the possibility to compensate the divergence by the Offset. So you can balance differences of about $\pm 10^{\circ}\text{C}$. The Offset is calculated as follows:

Temperature measured by you – showed Temperature (PV) = Offset

- To reach this menu you have to push the SUB button in the menu item “set temperature”.
- You can recognize this menu at the PV showed in the SV-line.
- If you turn the Jog-Shuttle knob.
- Set the calculated value by turning the Jog-Shuttle. The Offset will be added to the PV immediately.
- Save and confirm the Offset by pushing the MAIN button, you will also reach the previous menu.

6.6 Key lock

- To avoid accidental changing of the set values, push the MAIN button in the menu item "set temperature".
- In this menu item can you see the actually temperature in the SV-line and the remaining time of turn-off delay in the PV-line. The function of the Jog-Shuttle knob is deactivated.
- Press the MAIN button once more to leave the key lock. Then you are in the menu item "set turn-off delay".

6.7 Storage function

The **WCT Cold Trap Bath** includes a storage function for temperature and time values. That is practical if you want to work with the same parameters often. Then you didn't need to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Any stored parameters are erasable. Turn the value with the Jug-Shuttle till "0" and then hold the knob pressed for two seconds.
- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Shuttle knob for storing.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown.
Unit Not Refrigerating	<ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check if the set temperature (SV) confirmed with the Jog-Shuttle knob (by turn the Jog-Shuttle knob cannot change the set temperature – see point 6.4) ✓ Check that the set temperature (SV) is lower than 35 °C.
Refrigerating but it doesn't fall to the SV temperature	<ul style="list-style-type: none"> ✓ Problem with the cooling circuit.
ERROR 1	<ul style="list-style-type: none"> ✓ Temperature Sensor or main controller Error.

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

9 Technical data

		WCT-40	WCT-80
Dimension (w×d×h)	Bath (cm)	220×220×210	220×220×210
	Extern (cm)	377×453×705	503×553×740
Capacity		10 Lit.	
Temp.	Range	Down to -40 °C	Down to -80 °C
	Accuracy	PT100	
Controller		Digital fuzzy controller with Jog-Shuttle knob (turn + push)	
Display		Illuminated LCD	
Timer		99 h 59 min, turn-on & -off delay + endurance operation	
Material	Intern	Stainless steel	
	Extern	Powder coated steel	
	Insulation	Polystyrene (20 mm)	
Pump		14 lit / min (intern / extern)	
Refrigerator		900 W	1.800 W
Refrigerant		404 A	
Others		<ul style="list-style-type: none"> - Storage function (temp., timer) - Key lock (Jog-Shuttle input disabled) - Easy justification - Alarm (for error status and timer-end) 	
Power supply		AC 230 V, 1N~, 50/60 Hz	

10 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	WCT Cold Trap Bath
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de